

## PEARL Procedures

Igor procedures for the analysis of PEARL data

Version rev-distro-1.4.0-1-g0a436db-dirty

Thu Sep 21 2017 12:35:04



# Contents

<b>1</b>	<b>Introduction</b>	<b>1</b>
1.1	Introduction . . . . .	1
1.2	Installation . . . . .	1
1.3	License Information . . . . .	1
<b>2</b>	<b>Projections</b>	<b>3</b>
<b>3</b>	<b>Todo List</b>	<b>5</b>
<b>4</b>	<b>Module Index</b>	<b>7</b>
4.1	Packages . . . . .	7
<b>5</b>	<b>Namespace Index</b>	<b>9</b>
5.1	Namespace List . . . . .	9
<b>6</b>	<b>Data Structure Index</b>	<b>11</b>
6.1	Data Structures . . . . .	11
<b>7</b>	<b>File Index</b>	<b>13</b>
7.1	File List . . . . .	13
<b>8</b>	<b>Module Documentation</b>	<b>15</b>
8.1	ARPES package . . . . .	15
8.1.1	Detailed Description . . . . .	15
<b>9</b>	<b>Namespace Documentation</b>	<b>17</b>
9.1	PearlAnglescanProcess Namespace Reference . . . . .	17
9.1.1	Detailed Description . . . . .	17
9.2	PearlAreaDisplay Namespace Reference . . . . .	17
9.2.1	Detailed Description . . . . .	17
9.2.2	Miscellaneous . . . . .	17
9.3	PearlAreaImport Namespace Reference . . . . .	18
9.3.1	Detailed Description . . . . .	18
9.4	PearlAreaProfiles Namespace Reference . . . . .	18

9.4.1 Detailed Description . . . . .	18
9.5 PearlArpes Namespace Reference . . . . .	18
9.5.1 Detailed Description . . . . .	18
9.6 PearlDataExplorer Namespace Reference . . . . .	18
9.6.1 Detailed Description . . . . .	19
9.7 PearlElog Namespace Reference . . . . .	19
9.7.1 Detailed Description . . . . .	19
9.8 PearlMatrixImport Namespace Reference . . . . .	19
9.8.1 Detailed Description . . . . .	19
9.9 PearlPShellImport Namespace Reference . . . . .	19
9.9.1 Detailed Description . . . . .	19
9.10 PearlScientaPreprocess Namespace Reference . . . . .	19
9.10.1 Detailed Description . . . . .	19
<b>10 Data Structure Documentation . . . . .</b>	<b>21</b>
10.1 DoniachSunjicStruct Struct Reference . . . . .	21
10.1.1 Detailed Description . . . . .	21
10.1.2 Field Documentation . . . . .	21
10.1.2.1 broadening . . . . .	21
10.1.2.2 convolution . . . . .	21
10.1.2.3 model . . . . .	21
10.1.2.4 oversampling . . . . .	21
10.1.2.5 precision . . . . .	22
10.1.2.6 pw . . . . .	22
10.1.2.7 xdw . . . . .	22
10.1.2.8 xw . . . . .	22
10.1.2.9 yw . . . . .	22
10.2 errorCode Struct Reference . . . . .	22
10.2.1 Detailed Description . . . . .	22
10.2.2 Field Documentation . . . . .	23
10.2.2.1 ALREADY_FILE_OPEN . . . . .	23
10.2.2.2 EMPTY_RESULTFILE . . . . .	23
10.2.2.3 FILE_NOT_READABLE . . . . .	23
10.2.2.4 INTERNAL_ERROR_CONVERTING_DATA . . . . .	23
10.2.2.5 INVALID_RANGE . . . . .	23
10.2.2.6 NO_FILE_OPEN . . . . .	23
10.2.2.7 NO_NEW_BRICKLETS . . . . .	23
10.2.2.8 SUCCESS . . . . .	23
10.2.2.9 UNKNOWN_ERROR . . . . .	23
10.2.2.10 WAVE_EXIST . . . . .	23

10.2.2.11 WRONG_PARAMETER . . . . .	23
<b>11 File Documentation</b>	<b>25</b>
11.1 fermi-edge-analysis.ipf File Reference . . . . .	25
11.1.1 Detailed Description . . . . .	25
11.1.2 Function Documentation . . . . .	26
11.1.2.1 analyse_curved_edge(wave data) . . . . .	26
11.1.2.2 analyser_energy_resolution(variable epass, variable slit) . . . . .	26
11.1.2.3 FermiFuncLinDOS2D_corr(variable w, threadsafe x, wave y) . . . . .	26
11.1.2.4 integrate_curved_edge(wave data, wave data_sig) . . . . .	26
11.1.2.5 record_results(variable index) . . . . .	26
11.1.2.6 show_shift(wave data) . . . . .	27
11.1.2.7 slit_correction(wave data, wave data_out, variable epass) . . . . .	27
11.1.2.8 slit_shift(variable ypix, variable epass) . . . . .	27
11.1.3 Variable Documentation . . . . .	27
11.1.3.1 hemi_radius_mm . . . . .	27
11.1.3.2 mcp_radius_epass . . . . .	27
11.1.3.3 mcp_radius_mm . . . . .	27
11.1.3.4 mcp_radius_pix . . . . .	27
11.2 mainpage.dox File Reference . . . . .	27
11.3 pearl-anglescan-process.ipf File Reference . . . . .	27
11.3.1 Detailed Description . . . . .	30
11.3.2 Function Documentation . . . . .	31
11.3.2.1 calc_graph_azi(variable x, variable y, variable projection=defaultValue, variable zeroAngle=defaultValue) . . . . .	31
11.3.2.2 calc_graph_polar(variable x, variable y, variable projection=defaultValue) . . . . .	32
11.3.2.3 calc_graph_radius(variable polar, variable projection=defaultValue) . . . . .	32
11.3.2.4 calc_nth(variable Theta_st, variable Theta_in, variable th, variable Phi_ran, variable Phi_ref, string Holomode) . . . . .	32
11.3.2.5 calc_phi_step(variable Theta_in, variable th, variable Theta_st, variable Phi_ran, variable Phi_ref, string Holomode) . . . . .	33
11.3.2.6 Calc_The_step(variable th, variable Theta_st, string Holomode) . . . . .	33
11.3.2.7 CalcN_Theta(string HoloMode, variable Theta_in, variable Theta_ran, variable Theta_st) . . . . .	33
11.3.2.8 check_contrast(wave values, variable pccmin, variable pccmax, variable *vmin, variable *vmax) . . . . .	33
11.3.2.9 clear_hemi_grid(string nickname) . . . . .	33
11.3.2.10 convert_angles_ttpa2polar(wave theta, wave tilt, wave phi, wave analyser, wave polar, wave azi) . . . . .	33
11.3.2.11 convert_angles_ttpd2polar(wave theta, wave tilt, wave phi, wave data, wave polar, wave azi) . . . . .	34
11.3.2.12 crop_strip(wave strip, variable xlo, variable xhi) . . . . .	34

11.3.2.13 display_hemi_scan(string nickname, variable projection=defaultValue, variable graphtype=defaultValue, variable do_ticks=defaultValue, variable do_↵ grids=defaultValue, string graphname=defaultValue) . . . . .	34
11.3.2.14 display_polar_graph(string graphname, variable angle_offset=defaultValue, variable do_ticks=defaultValue) . . . . .	35
11.3.2.15 display_scanlines(string nickname, variable alpha_lo, variable alpha_hi, wave m_theta, wave m_tilt, wave m_phi, variable folding=defaultValue, variable projection=defaultValue) . . . . .	36
11.3.2.16 draw_diffraction_cone(string graphname, string groupname, variable theta_axis, variable theta_inner, variable phi) . . . . .	37
11.3.2.17 draw_hemi_axes(string graphname, variable do_grids=defaultValue) . . . . .	37
11.3.2.18 duplicate_hemi_scan(string source_nickname, dfref dest_folder, string dest_↵ nickname, variable xpdplot=defaultValue) . . . . .	38
11.3.2.19 find_hemi_data(string nickname, string *prefix, string *intwave) . . . . .	38
11.3.2.20 get_hemi_nickname(wave w) . . . . .	38
11.3.2.21 get_hemi_prefix(wave w) . . . . .	39
11.3.2.22 hemi_add_anglescan(string nickname, wave values, wave polar, wave azi, wave weights=defaultValue) . . . . .	39
11.3.2.23 hemi_add_aziscan(string nickname, wave values, variable polar, wave azi, wave weights=defaultValue) . . . . .	39
11.3.2.24 hemi_azi_cut(string nickname, variable pol) . . . . .	39
11.3.2.25 hemi_polar_cut(string nickname, variable azim) . . . . .	40
11.3.2.26 import_tpi_scan(string nickname, wave theta, wave phi, wave intensity, variable folding=defaultValue, variable npolar=defaultValue, variable nograph=default_↵ Value, variable xpdplot=defaultValue) . . . . .	40
11.3.2.27 interpolate_hemi_scan(string nickname) . . . . .	41
11.3.2.28 line_average(wave source, wave dest) . . . . .	41
11.3.2.29 load_hemi_scan(string nickname, string pathname, string filename) . . . . .	41
11.3.2.30 make_hemi_grid(variable npol, string nickname, variable xpdplot=defaultValue) .	41
11.3.2.31 normalize_strip_2d(wave strip, wave theta, variable theta_offset=defaultValue, variable smooth_method=defaultValue, variable smooth_factor=defaultValue, variable check=defaultValue) . . . . .	42
11.3.2.32 normalize_strip_phi(wave strip, wave theta, wave phi, variable theta_↵ offset=defaultValue, variable theta_range=defaultValue, variable check=default_↵ Value) . . . . .	42
11.3.2.33 normalize_strip_theta(wave strip, wave theta, variable theta_offset=defaultValue, variable smooth_method=defaultValue, variable smooth_factor=defaultValue, variable check=defaultValue) . . . . .	43
11.3.2.34 normalize_strip_x(wave strip, variable smooth_method=defaultValue, variable smooth_factor=defaultValue, variable check=defaultValue) . . . . .	44
11.3.2.35 pizza_service(wave data, string nickname, variable theta_offset, variable tilt_offset, variable phi_offset, variable npolar=defaultValue, variable nograph=defaultValue, variable folding=defaultValue, variable xpdplot=defaultValue) . . . . .	44
11.3.2.36 pizza_service_2(wave data, string nickname, wave m_theta, wave m_tilt, wave m_phi, variable npolar=defaultValue, variable nograph=defaultValue, variable folding=defaultValue, variable xpdplot=defaultValue) . . . . .	45
11.3.2.37 polar_graph_hook(WMWinHookStruct *s) . . . . .	46

11.3.2.38 quick_pizza_image(wave data, string nickname, variable theta_offset, variable tilt_offset, variable phi_offset, variable npolar=defaultValue, variable nograph=defaultValue, variable folding=defaultValue)	46
11.3.2.39 rotate_hemi_scan(string nickname, variable angle)	47
11.3.2.40 save_hemi_scan(string nickname, string pathname, string filename)	47
11.3.2.41 set_contrast(variable pccmin, variable pccmax, string graphname=defaultValue, string colortable=defaultValue)	47
11.3.2.42 set_polar_graph_cursor(string nickname, string cursorname, variable polar_angle, variable azimuth_angle, string graphname=defaultValue)	47
11.3.2.43 show_analyser_line(variable theta, variable tilt, variable phi, variable theta_offset, variable tilt_offset, variable phi_offset, variable npolar=defaultValue, variable nograph=defaultValue, variable xpdplot=defaultValue)	47
11.3.2.44 strip_delete_frames(wave strip, variable qlo, variable qhi, wave theta, wave tilt, wave phi)	48
11.3.2.45 trim_hemi_scan(string nickname, variable theta_max)	48
11.3.2.46 update_polar_info(string graphname)	49
11.3.3 Variable Documentation	49
11.3.3.1 kProjArea	49
11.3.3.2 kProjDist	49
11.3.3.3 kProjGnom	49
11.3.3.4 kProjOrtho	49
11.3.3.5 kProjScaleArea	49
11.3.3.6 kProjScaleDist	49
11.3.3.7 kProjScaleGnom	50
11.3.3.8 kProjScaleOrtho	50
11.3.3.9 kProjScaleStereo	50
11.3.3.10 kProjStereo	50
11.4 pearl-anglescan-tracker.ipf File Reference	50
11.4.1 Function Documentation	52
11.4.1.1 add_image_data(wave image, variable theta, variable tilt, variable phi)	52
11.4.1.2 AfterCompiledHook()	52
11.4.1.3 ast_add_image(wave image, variable theta, variable tilt, variable phi)	52
11.4.1.4 ast_callback_data(variable chan)	52
11.4.1.5 ast_callback_detector(variable chan)	53
11.4.1.6 ast_callback_manip(variable chan)	53
11.4.1.7 ast_close()	53
11.4.1.8 ast_export(dfref folder, string nickname, variable xpdplot=defaultValue)	53
11.4.1.9 ast_import(string nickname)	53
11.4.1.10 ast_prepare(variable theta_offset=defaultValue, variable tilt_offset=defaultValue, variable phi_offset=defaultValue)	53
11.4.1.11 ast_set_processing(string reduction_func, string reduction_params)	54
11.4.1.12 ast_setup()	54

11.4.1.13	ast_update_detector(variable theta, variable tilt, variable phi, variable range)	54
11.4.1.14	ast_window_hook(WMWinHookStruct *s)	54
11.4.1.15	bp_capture(WMButtonAction *ba)	55
11.4.1.16	edit_offsets()	55
11.4.1.17	edit_reduction_params()	55
11.4.1.18	epics_connect()	55
11.4.1.19	epics_disconnect()	55
11.4.1.20	epics_disconnect_chid(string chid_var_name)	55
11.4.1.21	export_tracker_data()	55
11.4.1.22	extend_data(variable num_slices)	55
11.4.1.23	IgorQuitHook(string app)	56
11.4.1.24	import_tracker_data()	56
11.4.1.25	init_package()	56
11.4.1.26	load_prefs()	56
11.4.1.27	load_tracker_data()	56
11.4.1.28	pmp_data(WMPopupAction *pa)	56
11.4.1.29	pmp_data_mouseup(WMPopupAction *pa)	56
11.4.1.30	pmp_parameters(WMPopupAction *pa)	56
11.4.1.31	pmp_parameters_mouseup(WMPopupAction *pa)	56
11.4.1.32	process_image_data()	56
11.4.1.33	save_prefs()	57
11.4.1.34	save_tracker_data()	57
11.4.1.35	setup_data()	57
11.4.1.36	setup_detector()	57
11.4.1.37	setup_graph()	57
11.4.1.38	toggle_capture()	57
11.4.1.39	update_capture()	57
11.4.1.40	update_data_graph()	57
11.4.1.41	update_detector(variable theta, variable tilt, variable phi, variable range)	57
11.4.1.42	update_detector_graph()	58
11.4.2	Variable Documentation	58
11.4.2.1	package_path	58
11.4.2.2	prefs_objects	58
11.4.2.3	version	58
11.5	pearl-area-display.ipf File Reference	58
11.5.1	Detailed Description	60
11.5.2	2D data	60
11.5.3	3D data	60
11.5.4	Function Documentation	60
11.5.4.1	ad_box_filter(wave image, string options)	60

11.5.4.2	<a href="#">ad_brick_slicer(wave data)</a>	61
11.5.4.3	<a href="#">ad_calc_cursor_profiles(wave image)</a>	61
11.5.4.4	<a href="#">ad_calc_histogram(wave image)</a>	61
11.5.4.5	<a href="#">ad_calc_profiles(wave image, variable pa, variable qa, variable pb, variable qb)</a>	61
11.5.4.6	<a href="#">ad_default_image_filter(wave image, string options)</a>	62
11.5.4.7	<a href="#">ad_display(wave image)</a>	62
11.5.4.8	<a href="#">ad_display_brick(wave data)</a>	62
11.5.4.9	<a href="#">ad_display_histogram(wave image)</a>	63
11.5.4.10	<a href="#">ad_display_profiles(wave image, string filter=defaultValue)</a>	63
11.5.4.11	<a href="#">ad_display_slice(wave data)</a>	63
11.5.4.12	<a href="#">ad_export_profile(wave view_image, variable dim, variable trace=defaultValue, variable show=defaultValue, variable overwrite=defaultValue)</a>	64
11.5.4.13	<a href="#">ad_gizmo_set_plane(wave brick, variable dim, variable value)</a>	64
11.5.4.14	<a href="#">ad_profiles_crosshairs(wave image, variable clear=defaultValue)</a>	65
11.5.4.15	<a href="#">ad_profiles_cursor_mode(wave image, variable mode)</a>	65
11.5.4.16	<a href="#">ad_profiles_hook(WMWinHookStruct *s)</a>	65
11.5.4.17	<a href="#">ad_profiles_set_cursor(wave image, string cursormode, variable xa, variable ya, variable pscale=defaultValue)</a>	66
11.5.4.18	<a href="#">ad_profiles_set_slice(wave brick, variable dim, variable value)</a>	66
11.5.4.19	<a href="#">ad_slicer_init_bg()</a>	66
11.5.4.20	<a href="#">ad_slicer_move_bg(WMBackgroundStruct *s)</a>	66
11.5.4.21	<a href="#">ad_slicer_start_bg(wave brick, variable dimension, string posvariable, variable delta)</a>	66
11.5.4.22	<a href="#">ad_slicer_stop_bg(string posvariable)</a>	67
11.5.4.23	<a href="#">ad_transpose_filter(wave image, string options)</a>	67
11.5.4.24	<a href="#">ad_update_profiles(wave image)</a>	67
11.5.4.25	<a href="#">bp_extract_slice(WMButtonAction *ba)</a>	67
11.5.4.26	<a href="#">bp_move_slice(WMButtonAction *ba)</a>	68
11.5.4.27	<a href="#">bp_move_slice_center(wave brick, variable dim, string posvariable)</a>	68
11.5.4.28	<a href="#">bp_reset_cursors(WMButtonAction *ba)</a>	68
11.5.4.29	<a href="#">get_source_image(wave view)</a>	68
11.5.4.30	<a href="#">get_view_folder(wave source)</a>	68
11.5.4.31	<a href="#">get_view_image(wave source)</a>	68
11.5.4.32	<a href="#">graphname_from_dfref(dfref df, string prefix)</a>	68
11.5.4.33	<a href="#">make_view_folder(wave source)</a>	69
11.5.4.34	<a href="#">pmp_export(WMPopupAction *pa)</a>	69
11.5.4.35	<a href="#">set_trace_colors(string graphname)</a>	69
11.5.4.36	<a href="#">slp_slice_position(WMSliderAction *sa)</a>	69
11.5.4.37	<a href="#">svp_slice_position(WMSetVariableAction *sva)</a>	69
11.5.4.38	<a href="#">svp_smoothing(WMSetVariableAction *sva)</a>	69
11.5.4.39	<a href="#">update_slice_info()</a>	69

11.6	pearl-area-import.ipf File Reference	69
11.6.1	Detailed Description	71
11.6.2	Function Documentation	71
11.6.2.1	ad_load_dialog(string APathName)	71
11.6.2.2	ad_suggest_foldername(string filename, variable ignoredate=defaultValue, string sourcename=defaultValue, variable unique=defaultValue)	71
11.6.2.3	adh5_default_reduction(wave source, wave dest1, wave dest2, string *param)	72
11.6.2.4	adh5_list_reduction_funcs()	72
11.6.2.5	adh5_load_complete(string ANickName, string APathName, string AFileName, variable load_data=defaultValue, variable load_attr=defaultValue)	72
11.6.2.6	adh5_load_detector(variable fileID, string detectorpath)	73
11.6.2.7	adh5_load_detector_image(variable fileID, string detectorpath, variable dim2start, variable dim2count, variable dim3start, variable dim3count)	73
11.6.2.8	adh5_load_detector_slabs(variable fileID, string detectorpath, variable progress=defaultValue)	73
11.6.2.9	adh5_load_info(string APathName, string AFileName)	74
11.6.2.10	adh5_load_preview(string ANickName, string APathName, string AFileName, variable load_data=defaultValue, variable load_attr=defaultValue)	74
11.6.2.11	adh5_load_reduced(string ANickName, string APathName, string AFileName, funcref reduction_func, string reduction_param, variable load_data=defaultValue, variable load_attr=defaultValue, variable progress=defaultValue)	74
11.6.2.12	adh5_load_reduced_detector(variable fileID, string detectorpath, funcref reduction_func, string reduction_param, variable progress=defaultValue, variable nthreads=defaultValue)	75
11.6.2.13	adh5_loadattr_all(variable fileID, string attributespath)	75
11.6.2.14	adh5_redim(wave data)	76
11.6.2.15	adh5_scale(wave data, string source=defaultValue)	76
11.6.2.16	adh5_scale_scan(wave data)	76
11.6.2.17	adh5_scale_scienta(wave data)	76
11.6.2.18	adh5_setup_profile(wave image, wave profile, variable dim)	76
11.6.2.19	adh5_test_reduction_func(wave source, wave dest1, wave dest2, funcref reduction_func, string reduction_param)	77
11.6.2.20	BeforeFileOpenHook(variable refNum, string fileName, string path, string type, string creator, variable kind)	77
11.6.2.21	GetAttrDataFolderDFR(wave data)	77
11.6.2.22	read_attribute_info(string datawavename, string source, variable *idest)	77
11.6.2.23	reduce_slab_image(wave slabdata, wave image, wave profile1, wave profile2, funcref reduction_func, string reduction_param)	78
11.6.2.24	reduce_slab_worker(funcref reduction_func)	78
11.7	pearl-area-profiles.ipf File Reference	78
11.7.1	Detailed Description	79
11.7.2	Function Documentation	79
11.7.2.1	ad_collect_multiscan_y(wave dataset, wave positions, wave destwave, variable noavg=defaultValue)	79

11.7.2.2	<code>ad_extract_rod(wave dataset, variable x1, variable x2, variable y1, variable y2, variable z1, variable z2, string destname, variable noavg=defaultValue, variable sdev=defaultValue, variable pscale=defaultValue)</code>	79
11.7.2.3	<code>ad_extract_rod_x(wave dataset, variable q1, variable q2, variable r1, variable r2, string destname, variable noavg=defaultValue, variable sdev=defaultValue)</code>	80
11.7.2.4	<code>ad_extract_rod_y(wave dataset, variable p1, variable p2, variable r1, variable r2, string destname, variable noavg=defaultValue, variable sdev=defaultValue)</code>	80
11.7.2.5	<code>ad_extract_rod_z(wave dataset, variable p1, variable p2, variable q1, variable q2, string destname, variable noavg=defaultValue, variable sdev=defaultValue)</code>	80
11.7.2.6	<code>ad_extract_slab(wave dataset, variable x1, variable x2, variable y1, variable y2, variable z1, variable z2, string destname, variable noavg=defaultValue, variable pscale=defaultValue)</code>	80
11.7.2.7	<code>ad_extract_slab_x(wave dataset, variable p1, variable p2, string destname, variable noavg=defaultValue)</code>	81
11.7.2.8	<code>ad_extract_slab_y(wave dataset, variable q1, variable q2, string destname, variable noavg=defaultValue)</code>	81
11.7.2.9	<code>ad_extract_slab_z(wave dataset, variable r1, variable r2, string destname, variable noavg=defaultValue)</code>	81
11.7.2.10	<code>ad_profile_x(wave dataset, variable q1, variable q2, string destname, variable noavg=defaultValue)</code>	81
11.7.2.11	<code>ad_profile_x_w(wave dataset, variable q1, variable q2, wave destwave, variable noavg=defaultValue)</code>	81
11.7.2.12	<code>ad_profile_y(wave dataset, variable p1, variable p2, string destname, variable noavg=defaultValue)</code>	81
11.7.2.13	<code>ad_profile_y_w(wave dataset, variable p1, variable p2, wave destwave, variable noavg=defaultValue)</code>	81
11.7.2.14	<code>calc_y_profile_mins(wave image)</code>	81
11.8	pearl-arpes.ipf File Reference	81
11.8.1	Detailed Description	82
11.8.2	Function Documentation	82
11.8.2.1	<code>AfterCompiledHook()</code>	82
11.8.2.2	<code>UnloadPearlArpesPackage()</code>	83
11.9	pearl-data-explorer.ipf File Reference	83
11.9.1	Detailed Description	84
11.9.2	Function Documentation	85
11.9.2.1	<code>attributes_notebook(wave attr_names, wave attr_values, string title)</code>	85
11.9.2.2	<code>bp_attr_notebook(WMButtonAction *ba)</code>	85
11.9.2.3	<code>bp_browse_filepath(WMButtonAction *ba)</code>	85
11.9.2.4	<code>bp_dataset_display(WMButtonAction *ba)</code>	85
11.9.2.5	<code>bp_dataset_folder(WMButtonAction *ba)</code>	85
11.9.2.6	<code>bp_dataset_next(WMButtonAction *ba)</code>	85
11.9.2.7	<code>bp_dataset_prev(WMButtonAction *ba)</code>	85
11.9.2.8	<code>bp_file_next(WMButtonAction *ba)</code>	85
11.9.2.9	<code>bp_file_prev(WMButtonAction *ba)</code>	85

11.9.2.10 bp_load_files(WMButtonAction *ba) . . . . .	85
11.9.2.11 bp_load_files_opt(WMButtonAction *ba) . . . . .	85
11.9.2.12 bp_load_prefs(WMButtonAction *ba) . . . . .	85
11.9.2.13 bp_save_prefs(WMButtonAction *ba) . . . . .	86
11.9.2.14 bp_update_datasets(WMButtonAction *ba) . . . . .	86
11.9.2.15 bp_update_filelist(WMButtonAction *ba) . . . . .	86
11.9.2.16 display_dataset(string datasetname) . . . . .	86
11.9.2.17 display_preview_trace(wave xtrace, wave ytrace) . . . . .	86
11.9.2.18 extract_preview_image(wave data, wave preview) . . . . .	86
11.9.2.19 init_package() . . . . .	86
11.9.2.20 itx_suggest_foldername(string filename, variable ignoredate=defaultValue, string sourcename=defaultValue, variable unique=defaultValue) . . . . .	86
11.9.2.21 lbp_datasets(WMListboxAction *lba) . . . . .	86
11.9.2.22 lbp_filelist(WMListboxAction *lba) . . . . .	86
11.9.2.23 load_file(string filename, string options=defaultValue) . . . . .	86
11.9.2.24 load_hdf_file(string filename, string options=defaultValue) . . . . .	87
11.9.2.25 load_itx_file(string filename, string options=defaultValue) . . . . .	87
11.9.2.26 load_mtrx_file(string filename, string options=defaultValue) . . . . .	87
11.9.2.27 load_prefs() . . . . .	87
11.9.2.28 load_pshell_file(string filename, string options=defaultValue) . . . . .	87
11.9.2.29 load_selected_files(string options=defaultValue) . . . . .	87
11.9.2.30 notebook_add_attributes(string notebook_name, wave attr_filter, wave attr_↔ names, wave attr_values) . . . . .	87
11.9.2.31 pearl_data_explorer() . . . . .	87
11.9.2.32 pearl_file_type(string filename) . . . . .	87
11.9.2.33 PearlDataExplorer() . . . . .	88
11.9.2.34 preview_attributes(dfref attr_folder, dfref dest_folder=defaultValue, wave attr_↔ filter=defaultValue, variable include_datawaves=defaultValue, variable include_↔ infowaves=defaultValue) . . . . .	88
11.9.2.35 preview_datafolder() . . . . .	88
11.9.2.36 preview_dataset(string datasetname) . . . . .	88
11.9.2.37 preview_file(string filename) . . . . .	88
11.9.2.38 preview_hdf_file(string filename) . . . . .	88
11.9.2.39 preview_itx_file(string filename) . . . . .	88
11.9.2.40 preview_mtrx_file(string filename) . . . . .	88
11.9.2.41 preview_pshell_file(string filename) . . . . .	89
11.9.2.42 preview_setscale_x(wave data, wave preview) . . . . .	89
11.9.2.43 prompt_default_process(string *param) . . . . .	89
11.9.2.44 prompt_func_params(string func_name, string *func_param) . . . . .	89
11.9.2.45 prompt_hdf_options(string *options) . . . . .	89
11.9.2.46 save_prefs() . . . . .	90

11.9.2.47	<code>show_preview_graph(wave data, wave xdata=defaultValue)</code>	90
11.9.2.48	<code>test_attributes_notebook()</code>	90
11.9.2.49	<code>update_datasets()</code>	90
11.9.2.50	<code>update_filelist()</code>	90
11.9.3	Variable Documentation	90
11.9.3.1	<code>ks_filematch_adh5</code>	90
11.9.3.2	<code>ks_filematch_itx</code>	90
11.9.3.3	<code>ks_filematch_mtrx</code>	90
11.9.3.4	<code>ks_filematch_pshell</code>	90
11.9.3.5	<code>package_name</code>	90
11.9.3.6	<code>package_path</code>	90
11.10	pearl-elog.ipf File Reference	91
11.10.1	Detailed Description	93
11.10.2	Function Documentation	94
11.10.2.1	<code>AfterFileOpenHook(variable refNum, string file, string pathName, string type, string creator, variable kind)</code>	94
11.10.2.2	<code>bp_attach(WMButtonAction *ba)</code>	94
11.10.2.3	<code>bp_attach_allnone(WMButtonAction *ba)</code>	94
11.10.2.4	<code>bp_attach_top(WMButtonAction *ba)</code>	94
11.10.2.5	<code>bp_attach_updown(WMButtonAction *ba)</code>	94
11.10.2.6	<code>bp_clear(WMButtonAction *ba)</code>	95
11.10.2.7	<code>bp_login(WMButtonAction *ba)</code>	95
11.10.2.8	<code>bp_logout(WMButtonAction *ba)</code>	95
11.10.2.9	<code>bp_save_graphs(WMButtonAction *ba)</code>	95
11.10.2.10	<code>bp_submit(WMButtonAction *ba)</code>	95
11.10.2.11	<code>cleanup_temp_files()</code>	95
11.10.2.12	<code>create_cmd_file(string cmd)</code>	95
11.10.2.13	<code>create_graph_file(string graphname, variable fileindex)</code>	95
11.10.2.14	<code>create_message_file(string message)</code>	96
11.10.2.15	<code>elog_add_attachment(string logbook, variable id, string graphs)</code>	96
11.10.2.16	<code>elog_config(string elog_path=defaultValue, string hostname=defaultValue, variable port=defaultValue, string subdir=defaultValue)</code>	97
11.10.2.17	<code>elog_create_entry(string logbook, string attributes, string message, variable encoding=defaultValue, string graphs=defaultValue, variable replyto=defaultValue)</code>	97
11.10.2.18	<code>elog_create_logbook(string name, string template=defaultValue)</code>	97
11.10.2.19	<code>elog_init_pearl_templates()</code>	98
11.10.2.20	<code>elog_login(string logbook, string username, string password)</code>	98
11.10.2.21	<code>elog_logout(string logbook)</code>	98
11.10.2.22	<code>elog_panel_hook(WMWinHookStruct *s)</code>	98
11.10.2.23	<code>elog_prompt_logbook()</code>	98
11.10.2.24	<code>elog_prompt_login(string logbook)</code>	99

11.10.2.25	elog_validate_attributes(string logbook, string attributes)	99
11.10.2.26	format_url(string logbook)	99
11.10.2.27	get_default_panel_name()	99
11.10.2.28	get_elog_df(string name, variable category)	99
11.10.2.29	get_log_path()	100
11.10.2.30	get_panel_attributes(string windowname)	100
11.10.2.31	get_panel_graphs(string windowname)	100
11.10.2.32	get_panel_message(string windowname)	100
11.10.2.33	get_timestamp(string sep)	100
11.10.2.34	igorBeforeNewHook(string igorApplicationNameStr)	101
11.10.2.35	igorQuitHook(string igorApplicationNameStr)	101
11.10.2.36	init_package(variable clean=defaultValue)	101
11.10.2.37	init_volatile_vars()	101
11.10.2.38	list_logbooks(variable templates=defaultValue)	101
11.10.2.39	load_prefs()	102
11.10.2.40	move_attach_item(string logbook, variable item, variable distance)	102
11.10.2.41	parse_result()	102
11.10.2.42	pearl_elog(string logbook)	102
11.10.2.43	PearlElogPanel(string logbook)	102
11.10.2.44	prepare_command_line(string logbook)	102
11.10.2.45	prepare_graph_attachments(string graphs)	103
11.10.2.46	save_prefs()	103
11.10.2.47	set_panel_attributes(string windowname, string attributes, variable clear=defaultValue)	103
11.10.2.48	set_panel_graphs(string windowname, string graphs)	103
11.10.2.49	set_panel_message(string windowname, string message)	104
11.10.2.50	update_attach_items(string logbook)	104
11.10.3	Variable Documentation	104
11.10.3.1	elog_parse_id	104
11.10.3.2	elog_success_msg	104
11.10.3.3	kAttachColName	104
11.10.3.4	kAttachColSel	104
11.10.3.5	kAttachColTitle	104
11.10.3.6	kdfPersistent	104
11.10.3.7	kdfRoot	105
11.10.3.8	kdfTemplates	105
11.10.3.9	kdfVolatile	105
11.10.3.10	package_name	105
11.10.3.11	package_path	105
11.11	pearl-fitfuncs.ipf File Reference	105

11.11.1 Function Documentation	106
11.11.1.1 Au4f(wave w, variable x)	106
11.11.1.2 Au4f_2p2(wave w, variable x)	106
11.11.1.3 Au4f_2p3(wave w, variable x)	106
11.11.1.4 Calc_DoniachSunjicBroad(wave pw, wave yw)	106
11.11.1.5 DoniachSunjicBroad(wave pw, wave yw, wave xw)	106
11.11.1.6 DoniachSunjicBroadS(DoniachSunjicStruct *s)	106
11.11.1.7 ds1_bg(wave w, variable x)	106
11.11.1.8 ds2_bg(wave w, variable x)	106
11.11.1.9 ds4_bg(wave w, variable x)	106
11.11.1.10 ds6_bg(wave w, variable x)	106
11.11.1.11 FermiGaussConv(wave pw, wave yw, wave xw)	106
11.11.1.12 Fit_DoniachSunjicBroad(wave pw, wave yw, wave xw, wave ww)	107
11.11.1.13 ShowComponents_Au4f_2p2(wave coef_wave, wave fit_wave)	107
11.11.1.14 ShowComponents_Au4f_2p3(wave coef_wave, wave fit_wave)	107
11.11.2 Variable Documentation	107
11.11.2.1 version	107
11.12 pearl-gui-tools.ipf File Reference	107
11.12.1 Function Documentation	107
11.12.1.1 display_progress_panel(string title, string message, variable progress_max)	107
11.12.1.2 kill_progress_panel()	107
11.12.1.3 update_progress_panel(variable progress, string message=defaultValue, variable progress_max=defaultValue)	107
11.13 pearl-matrix-import.ipf File Reference	108
11.13.1 Detailed Description	109
11.13.2 Function Documentation	109
11.13.2.1 AfterFileOpenHook(variable refNum, string file, string pathName, string type, string creator, variable kind)	109
11.13.2.2 BeforeFileOpenHook(variable refNum, string fileName, string path, string type, string creator, variable kind)	109
11.13.2.3 check_package_folder()	110
11.13.2.4 init_package()	110
11.13.2.5 initStruct(errorCode *errorCode)	110
11.13.2.6 matrix_format_elog_message(wave metadata)	110
11.13.2.7 matrix_preview_2d(wave data, wave metadata)	110
11.13.2.8 mtrx_create_folder(string fileName, dfref df_base=defaultValue)	110
11.13.2.9 mtrx_file_brickletID(string resultFile, variable runCycle, variable scanCycle, string channel)	111
11.13.2.10 mtrx_get_cycle_folder(dfref df_base=defaultValue, variable runCycle=defaultValue, variable scanCycle=defaultValue)	111
11.13.2.11 mtrx_load_all()	111

11.13.2.12	<code>mtx_load_file(string pathName, string fileName, string traces=defaultValue)</code>	111
11.13.2.13	<code>mtx_load_info(string APathName, string AFileName)</code>	112
11.13.2.14	<code>mtx_load_preview(string destName, string pathName, string fileName, string traces=defaultValue)</code>	112
11.13.2.15	<code>mtx_open_file(string pathName, string fileNameOrPath)</code>	113
11.13.2.16	<code>mtx_parse_filename(string fileName, string *resultFile, variable *runCycle, variable *scanCycle, string *channel)</code>	114
11.13.2.17	<code>mtx_scale_dataset(wave data)</code>	114
11.13.2.18	<code>mtx_split_filename(string fileName, string *prefix, string *datepart, string *timepart)</code>	114
11.13.2.19	<code>preview_matrix_file(string filename)</code>	114
11.13.2.20	<code>subtract_line_bg(wave img)</code>	115
11.13.3	Variable Documentation	115
11.13.3.1	<code>ks_filematch_mtx</code>	115
11.13.3.2	<code>package_name</code>	115
11.13.3.3	<code>package_path</code>	115
11.14	pearl-menu.ipf File Reference	115
11.14.1	Function Documentation	115
11.14.1.1	<code>Display2dProfiles()</code>	115
11.14.1.2	<code>Display3dSlicer()</code>	115
11.14.1.3	<code>DisplayGizmoSlicer()</code>	116
11.14.1.4	<code>LoadPearlArpes()</code>	116
11.14.1.5	<code>LoadPearlOptics()</code>	116
11.14.1.6	<code>LoadPearlPreparation()</code>	116
11.14.1.7	<code>PearlAnglescanTracker(string epicsname, string wbRGB)</code>	116
11.14.1.8	<code>PearlLiveDisplay(string epicsname, string nickname, string wbRGB)</code>	116
11.14.1.9	<code>PearlMenuEnableFunc(string funcname)</code>	116
11.15	pearl-otf-import.ipf File Reference	116
11.15.1	Function Documentation	116
11.15.1.1	<code>gather_batch(string foldermatch, string ywavematch, string xwavematch, string destfolder)</code>	116
11.15.1.2	<code>otf_gather_batch(string ywavematch, string xwavematch, string destfolder)</code>	117
11.15.1.3	<code>otf_gather_iterator(dfref df, string *sdata)</code>	117
11.15.1.4	<code>otf_interp(variable e1, variable e2, variable npts, variable smo)</code>	117
11.15.1.5	<code>otf_load_itx(string pathname, string filename)</code>	117
11.15.1.6	<code>otf_load_itx_all(string pathname)</code>	117
11.15.1.7	<code>otf_load_itx_match(string pathname, string matchstr)</code>	117
11.15.1.8	<code>otf_rename_folders(string pattern, variable unique_index=defaultValue, string new_suffix=defaultValue, string match_str=defaultValue)</code>	117
11.15.1.9	<code>otf_rename_folders_iterator(dfref df, string *sdata)</code>	117
11.15.1.10	<code>otf_smo_int(wave win, wave wout, wave wpe, variable smo)</code>	117

11.16pearl-polar-coordinates.ipf File Reference . . . . .	117
11.16.1 Function Documentation . . . . .	118
11.16.1.1 cart2polar(variable xx, variable yy, variable zz, variable *radius, variable *theta, variable *phi) . . . . .	118
11.16.1.2 cart2polar_wave(wave in, wave out) . . . . .	118
11.16.1.3 polar2cart(variable radius, variable theta, variable phi, variable *xx, variable *yy, variable *zz) . . . . .	118
11.16.1.4 polar2cart_wave(wave in, wave out) . . . . .	118
11.16.1.5 polar_distance(variable polar1, variable azim1, variable polar2, variable azim2) . . . . .	118
11.17pearl-pshell-import.ipf File Reference . . . . .	118
11.17.1 Detailed Description . . . . .	120
11.17.2 Function Documentation . . . . .	121
11.17.2.1 find_attr_folder(dfref dataDF) . . . . .	121
11.17.2.2 find_scale_wave(string name, dfref dataDF, dfref scanDF, dfref attrDF) . . . . .	121
11.17.2.3 find_scan_folder(dfref dataDF) . . . . .	121
11.17.2.4 ps_detect_scale(wave ax, wave lo, wave hi, wave un) . . . . .	121
11.17.2.5 ps_scale_dataset(wave data) . . . . .	122
11.17.2.6 ps_scale_dataset_2(wave data, wave ax, wave lo, wave hi, wave un) . . . . .	122
11.17.2.7 ps_scale_datasets() . . . . .	123
11.17.2.8 ps_set_dimlabels(wave data) . . . . .	123
11.17.2.9 ps_set_dimlabels2(wave data, string name) . . . . .	124
11.17.2.10ps5_close_file(variable fileID) . . . . .	124
11.17.2.11ps5_list_scan_datasets(variable fileID, string scanpath, variable include_↵ regions=defaultValue) . . . . .	124
11.17.2.12ps5_list_scan_regions(variable fileID, string scanpath) . . . . .	125
11.17.2.13ps5_list_scans(variable fileID) . . . . .	125
11.17.2.14ps5_load_complete(string ANickName, string APathName, string AFileName, variable load_data=defaultValue, variable load_attr=defaultValue) . . . . .	125
11.17.2.15ps5_load_dataset(variable fileID, string scanpath, string datasetname, variable set_scale=defaultValue) . . . . .	126
11.17.2.16ps5_load_dataset_meta(variable fileID, string datapath, string datasetname, wave datawave) . . . . .	127
11.17.2.17ps5_load_dataset_reduced(variable fileID, string scanpath, string datasetname, funcref reduction_func, string reduction_param, variable progress=defaultValue, variable nthreads=defaultValue) . . . . .	127
11.17.2.18ps5_load_dataset_slab(variable fileID, string datapath, string datasetname, vari- able dim2start, variable dim2count, variable dim3start, variable dim3count) . . . . .	128
11.17.2.19ps5_load_dataset_slabs(variable fileID, string datapath, string datasetname, variable progress=defaultValue) . . . . .	129
11.17.2.20ps5_load_info(string APathName, string AFileName) . . . . .	129
11.17.2.21ps5_load_preview(string APathName, string AFileName, variable load_↵ data=defaultValue, variable load_attr=defaultValue, string pref_scans=default↵ Value, string pref_datasets=defaultValue) . . . . .	129

11.17.2.22	psh5_load_reduced(string ANickName, string APathName, string AFileName, funcref reduction_func, string reduction_param, variable progress=defaultValue)	130
11.17.2.23	psh5_load_scan_attrs(variable fileID, string scanpath, variable attr_sets=defaultValue)	131
11.17.2.24	psh5_load_scan_complete(variable fileID, string scanpath, variable load_data=defaultValue, variable load_attr=defaultValue)	131
11.17.2.25	psh5_load_scan_data(variable fileID, string scanpath)	132
11.17.2.26	psh5_load_scan_info(variable fileID, string scanpath)	132
11.17.2.27	psh5_load_scan_meta(variable fileID, string scanpath)	133
11.17.2.28	psh5_load_scan_preview(variable fileID, string scanpath, variable set_scale=defaultValue, string pref_datasets=defaultValue)	133
11.17.2.29	psh5_load_scan_section(variable fileID, string scanpath, variable dim, variable set_scale=defaultValue, string pref_datasets=defaultValue)	134
11.17.2.30	psh5_open_file(string ANickName, string APathName, string AFileName)	134
11.17.2.31	reduce_slab_image(wave slabdata, wave image, wave profile1, wave profile2, funcref reduction_func, string reduction_param)	135
11.17.2.32	reduce_slab_worker(funcref reduction_func)	135
11.17.2.33	select_dataset(string file_datasets, string pref_datasets)	135
11.17.2.34	wave2list(wave wt, string sep)	135
11.17.2.35	wave2list(wave w, string format, string sep)	135
11.17.3	Variable Documentation	136
11.17.3.1	kAngleDimLabel	136
11.17.3.2	kDataDimLabel	136
11.17.3.3	kDetectorSensitivity	136
11.17.3.4	kEnergyDimLabel	136
11.17.3.5	kPreviewDatasets	136
11.17.3.6	kScanDimLabel	136
11.17.3.7	kScientaScalingDatasets	136
11.17.3.8	kTransposedDatasets	136
11.18	pearl-scienta-preprocess.ipf File Reference	137
11.18.1	Detailed Description	137
11.18.2	Function Documentation	138
11.18.2.1	capture_int_linbg_cursors()	138
11.18.2.2	csr_int_linbg_reduction(string win)	138
11.18.2.3	fit_scienta_ang_transm(wave data, wave params)	138
11.18.2.4	fit_scienta_poly_bg(wave data, wave params, variable bgterms)	138
11.18.2.5	gauss2_reduction(wave source, wave dest1, wave dest2, string *param)	138
11.18.2.6	int_linbg_reduction(wave source, wave dest1, wave dest2, string *param)	139
11.18.2.7	int_quadbg_reduction(wave source, wave dest1, wave dest2, string *param)	139
11.18.2.8	prompt_gauss2_reduction(string *param)	139
11.18.2.9	prompt_int_linbg_reduction(string *param)	139
11.18.2.10	prompt_int_quadbg_reduction(string *param)	139

11.18.2.11	<code>prompt_redim_linbg_reduction(string *param)</code>	139
11.18.2.12	<code>prompt_Shockley_anglefit(string *param)</code>	139
11.18.2.13	<code>redim_linbg_reduction(wave source, wave dest1, wave dest2, string *param)</code>	139
11.18.2.14	<code>scienta_ang_transm(wave w, variable x)</code>	140
11.18.2.15	<code>scienta_norm(wave w, variable x)</code>	140
11.18.2.16	<code>scienta_poly_bg(wave w, variable e, variable a)</code>	140
11.18.2.17	<code>Shockley_anglefit(wave source, wave dest1, wave dest2, string *param)</code>	140
11.18.2.18	<code>test_gauss2_reduction(wave image)</code>	141
11.18.2.19	<code>test_int_linbg(wave image)</code>	141
11.18.2.20	<code>test_int_quadbg(wave image)</code>	141
11.18.2.21	<code>test_shockley_anglefit(wave image, variable branch)</code>	141
11.19	pearl-tools.ipf File Reference	141
11.19.1	Function Documentation	141
11.19.1.1	<code>AppendToGraphIterator(wave w, string *sdata)</code>	141
11.19.1.2	<code>DefaultFolderIterator(dfref df, string *sdata)</code>	141
11.19.1.3	<code>DefaultWaveIterator(wave w, string *sdata)</code>	141
11.19.1.4	<code>IterateDataFolders(string matchStr, funcref iterator, string sdata, string progress↵ _title=defaultValue)</code>	142
11.19.1.5	<code>IterateWaves(string matchStr, funcref iterator, string sdata)</code>	142
11.19.1.6	<code>SumWavesIterator(wave w, string *sdata)</code>	142
11.20	pearl-vector-operations.ipf File Reference	142
11.20.1	Function Documentation	142
11.20.1.1	<code>create_rotation_matrix_free()</code>	142
11.20.1.2	<code>rotate2d_x(variable xx, variable yy, variable angle)</code>	142
11.20.1.3	<code>rotate2d_y(variable xx, variable yy, variable angle)</code>	142
11.20.1.4	<code>rotate_x_wave(wave inout, variable angle)</code>	142
11.20.1.5	<code>rotate_y_wave(wave inout, variable angle)</code>	142
11.20.1.6	<code>rotate_z_wave(wave inout, variable angle)</code>	143
11.20.1.7	<code>set_rotation_x(wave matrix, variable angle)</code>	143
11.20.1.8	<code>set_rotation_y(wave matrix, variable angle)</code>	143
11.20.1.9	<code>set_rotation_z(wave matrix, variable angle)</code>	143
	<b>Index</b>	<b>145</b>



# Chapter 1

## Introduction

### 1.1 Introduction

PEARL Procedures is a suite of Igor Pro procedures developed for data acquisition and data processing at the PEARL beamline at the Swiss Light Source.

### 1.2 Installation

PEARL Procedures should be installed according to the regular Igor Pro guidelines. Please read the Igor help About Igor Pro User Files for details.

- Make a `pearl-procs` directory in your private or shared User Procedures folder, and copy the PEARL Procedures distribution there.
- Create shortcuts of the `pearl-arpes.ipf` and `pearl-menu.ipf` files, and move them to the Igor Procedures folder next to your User Procedures folder.
- Find the `HDF5.XOP` extension in the Igor Pro Folder under More Extensions/File Loaders, create a shortcut, and move the shortcut to the Igor Extensions folder next to your User Procedures folder.
- Find the `HDF5 Help.ihf` next to `HDF5.XOP`, create a shortcut, and move the shortcut to the Igor Help Files folder next to your User Procedures folder.

### 1.3 License Information

An open distribution of PEARL Procedures is available under the [Apache License, Version 2.0](https://git.psi.ch/pearl-public/igor-procs) at <https://git.psi.ch/pearl-public/igor-procs>. Users of PEARL Procedures are requested to coordinate and share the development of the code with the original author. Please read and respect the respective license agreements.

#### Author

Matthias Muntwiler, [matthias.muntwiler@psi.ch](mailto:matthias.muntwiler@psi.ch)

#### Version

This documentation is compiled from version `rev-distro-1.4.0-1-g0a436db-dirty`.

**Copyright**

2009-2016 by [Paul Scherrer Institut](#)  
Licensed under the [Apache License, Version 2.0](#)

## Chapter 2

# Projections

the functions of the anglescan package support the following map projections.

for a description of the different projections, see, for example, [https://en.wikipedia.org/wiki/Map\\_projection](https://en.wikipedia.org/wiki/Map_projection)

Selector	Projection	Function	Properties
kProjDist = 0	azimuthal equidistant	$r = c * \text{theta}$	radius is proportional to polar angle.
kProjStereo = 1	stereographic	$r = c * \tan \text{theta}/2$	circles on sphere map to circles.
kProjArea = 2	azimuthal equal-area	$r = c * \sin \text{theta}/2$	preserves area measure.
kProjGnom = 3	gnomonic	$r = c * \tan \text{theta}$	great circles map to straight lines.
kProjOrtho = 4	orthographic	$r = c * \sin \text{theta}$	k-space mapping in ARPES and LEED.

the projections in this package are defined for  $0 \leq \text{theta} < 90$ .



## Chapter 3

# Todo List

Global **adh5\_scale\_scan** (wave data)

incomplete

Global **elog\_validate\_attributes** (string logbook, string attributes)

function currently not implemented, always returns 0

Global **epics\_connect** ()

the X03DA channel names are hard-coded.

Global **load\_hemi\_scan** (string nickname, string pathname, string filename)

function not implemented

Global **mtrx\_open\_file** (string pathName, string fileNameOrPath)

fix possible cache issues, add an option to override the cache.

Global **psh5\_load\_reduced** (string ANickName, string APathName, string AFileName, funcref reduction\_↔  
func, string reduction\_param, variable progress=defaultValue)

load scan positions.



## Chapter 4

# Module Index

### 4.1 Packages

Packages are collections of procedure files to be loaded together. PEARL defines the following packages:

ARPES package . . . . . [15](#)



## Chapter 5

# Namespace Index

### 5.1 Namespace List

Here is a list of all namespaces with brief descriptions:

<a href="#">PearlAnglescanProcess</a>	Processing and holographic mapping of angle scanned XPD data . . . . .	17
<a href="#">PearlAreaDisplay</a>	Instant visualization of angle scan and manipulator position . . . . .	17
<a href="#">PearlAreaImport</a>	HDF5 file import from EPICS area detectors . . . . .	18
<a href="#">PearlAreaProfiles</a>	Profile extraction for multi-dimensional datasets acquired from area detectors . . . . .	18
<a href="#">PearlArpes</a>	Data acquisition and analysis package for ARPES at PEARL . . . . .	18
<a href="#">PearlDataExplorer</a>	Preview and import panel for PEARL data . . . . .	18
<a href="#">PearlElog</a>	Interface for writing ELOG entries with Igor graphs as attachment . . . . .	19
<a href="#">PearlMatrixImport</a>	Data file import for omicron matrix (STM) files . . . . .	19
<a href="#">PearlPShellImport</a>	Import data from PShell . . . . .	19
<a href="#">PearlScientaPreprocess</a>	Preprocessing functions for Scienta detector images . . . . .	19



## Chapter 6

# Data Structure Index

### 6.1 Data Structures

Here are the data structures with brief descriptions:

<a href="#">DoniachSunjicStruct</a>	21
<a href="#">errorCode</a>	
From matrixfilereader help	22



## Chapter 7

# File Index

### 7.1 File List

Here is a list of all files with brief descriptions:

<a href="#">fermi-edge-analysis.ipf</a>	
Tools for analysing the Fermi edge measured by the Scienta EW4000 analyser . . . . .	25
<a href="#">pearl-anglescan-process.ipf</a>	
Processing and holographic mapping of angle scanned XPD data . . . . .	27
<a href="#">pearl-anglescan-tracker.ipf</a>	
<a href="#">pearl-area-display.ipf</a>	
Visualization tools for 2D and 3D data . . . . .	58
<a href="#">pearl-area-import.ipf</a>	
HDF5 file import from EPICS area detectors . . . . .	69
<a href="#">pearl-area-profiles.ipf</a>	
Profile extraction for multi-dimensional datasets acquired from area detectors . . . . .	78
<a href="#">pearl-arpes.ipf</a>	
Data acquisition and analysis package for ARPES at PEARL . . . . .	81
<a href="#">pearl-data-explorer.ipf</a>	
Preview and import panel for PEARL data . . . . .	83
<a href="#">pearl-eelog.ipf</a>	
Interface for writing ELOG entries with Igor graphs as attachment . . . . .	91
<a href="#">pearl-fitfuncs.ipf</a>	
<a href="#">pearl-gui-tools.ipf</a>	
<a href="#">pearl-matrix-import.ipf</a>	
Data file import for omicron matrix (STM) files . . . . .	108
<a href="#">pearl-menu.ipf</a>	
<a href="#">pearl-otf-import.ipf</a>	
<a href="#">pearl-polar-coordinates.ipf</a>	
<a href="#">pearl-pshell-import.ipf</a>	
Import data from PShell . . . . .	118
<a href="#">pearl-scienta-preprocess.ipf</a>	
Preprocessing functions for Scienta detector images . . . . .	137
<a href="#">pearl-tools.ipf</a>	
<a href="#">pearl-vector-operations.ipf</a>	
	142



## Chapter 8

# Module Documentation

### 8.1 ARPES package

data acquisition and analysis package for ARPES at PEARL.

#### Files

- file [fermi-edge-analysis.ipf](#)  
*tools for analysing the Fermi edge measured by the Scienta EW4000 analyser.*
- file [pearl-anglescan-process.ipf](#)  
*processing and holographic mapping of angle scanned XPD data.*
- file [pearl-area-display.ipf](#)  
*visualization tools for 2D and 3D data.*
- file [pearl-area-import.ipf](#)  
*HDF5 file import from EPICS area detectors.*
- file [pearl-area-profiles.ipf](#)  
*profile extraction for multi-dimensional datasets acquired from area detectors.*
- file [pearl-data-explorer.ipf](#)  
*preview and import panel for PEARL data*
- file [pearl-elog.ipf](#)  
*interface for writing ELOG entries with Igor graphs as attachment.*
- file [pearl-pshell-import.ipf](#)  
*import data from PShell*
- file [pearl-scienta-preprocess.ipf](#)  
*preprocessing functions for Scienta detector images.*

#### 8.1.1 Detailed Description

data acquisition and analysis package for ARPES at PEARL.

The purpose of a package is to load a bunch of dependent procedure files. The ARPES package loads the following files which are helpful when working with ARPES data (any data from the Scienta analyser, that is) from PEARL.

- [pearl-area-import.ipf](#)
- [pearl-area-display.ipf](#)
- [pearl-area-profiles.ipf](#)

- [pearl-data-explorer.ipf](#)
- [pearl-scienta-preprocess.ipf](#)
- [pearl-anglescan-process.ipf](#)
- [pearl-anglescan-tracker.ipf](#)
- [pearl-elog.ipf](#)

Most of these files require the HDF5.XOP. The following files are loaded if the EPICS.XOP is present:

- [pearl-area-live.ipf](#)
- [pearl-epics.ipf](#)
- [pearl-arpes-scans.ipf](#)

## Chapter 9

# Namespace Documentation

### 9.1 PearlAnglescanProcess Namespace Reference

processing and holographic mapping of angle scanned XPD data.

#### 9.1.1 Detailed Description

processing and holographic mapping of angle scanned XPD data.

[PearlAnglescanProcess](#) is declared in [pearl-anglescan-process.ipf](#).

### 9.2 PearlAreaDisplay Namespace Reference

instant visualization of angle scan and manipulator position.

#### 9.2.1 Detailed Description

instant visualization of angle scan and manipulator position.

- 3D graphics in a Gizmo window
- 2D graphics with profiles
- Slicer panel

data dictionary of global variables in view\_XXXX folder:

- gizmo\_graphname = name of the gizmo window
- slice\_graphname = name of the slice/profiles graph window
- slice\_panelname = name of the slicer panel

#### 9.2.2 Miscellaneous

scienta HDF5 import, display, and slicer how-to

```
/// ad_display_brick(root:x03da_scienta_20130821_01560:data)
/// ad_scale_scienta(root:x03da_scienta_20130821_01560:data)
/// ad_scale_extra(root:x03da_scienta_20130821_01560:data,root:x03da_scienta_20130821_01560:MonoEnergy,root:x03da_scienta_20130821_01560:ManipulatorPosition)
```

```
/// ad_display_brick(root:x03da_scienta_20130821_01560:data)
/// ad_brick_slicer(root:x03da_scienta_20130821_01560:data, "gizmo_x03da_scienta_20130821_01")
///
```

```
@author matthias muntwiler, matthias.muntwiler@psi.ch
```

```
@copyright 2013-15 Paul Scherrer Institut @n
Licensed under the Apache License, Version 2.0 (the "License"); @n
you may not use this file except in compliance with the License. @n
You may obtain a copy of the License at
http://www.apache.org/licenses/LICENSE-2.0
```

[PearlAreaDisplay](#) is declared in [pearl-area-display.ipf](#).

## 9.3 PearlAreaImport Namespace Reference

HDF5 file import from EPICS area detectors.

### 9.3.1 Detailed Description

HDF5 file import from EPICS area detectors.

[PearlAreaImport](#) is declared in [pearl-area-import.ipf](#).

## 9.4 PearlAreaProfiles Namespace Reference

profile extraction for multi-dimensional datasets acquired from area detectors.

### 9.4.1 Detailed Description

profile extraction for multi-dimensional datasets acquired from area detectors.

[PearlAnglescanTracker](#) is declared in [pearl-area-profiles.ipf](#).

## 9.5 PearlArpes Namespace Reference

data acquisition and analysis package for ARPES at PEARL.

### 9.5.1 Detailed Description

data acquisition and analysis package for ARPES at PEARL.

[PearlArpes](#) is declared in [pearl-arpes.ipf](#).

## 9.6 PearlDataExplorer Namespace Reference

preview and import panel for PEARL data

### 9.6.1 Detailed Description

preview and import panel for PEARL data

[PearlDataExplorer](#) is declared in [pearl-data-explorer.ipf](#).

## 9.7 PearlElog Namespace Reference

interface for writing ELOG entries with Igor graphs as attachment.

### 9.7.1 Detailed Description

interface for writing ELOG entries with Igor graphs as attachment.

[PearlElog](#) is declared in [pearl-elog.ipf](#).

## 9.8 PearlMatrixImport Namespace Reference

data file import for omicron matrix (STM) files

### 9.8.1 Detailed Description

data file import for omicron matrix (STM) files

[PearlMatrixImport](#) is declared in [pearl-matrix-import.ipf](#).

## 9.9 PearlPShellImport Namespace Reference

import data from PShell

### 9.9.1 Detailed Description

import data from PShell

[PearlPShellImport](#) is declared in [pearl-pshell-import.ipf](#).

## 9.10 PearlScientaPreprocess Namespace Reference

preprocessing functions for Scienta detector images.

### 9.10.1 Detailed Description

preprocessing functions for Scienta detector images.

[PearlScientaPreprocess](#) is declared in [pearl-scienta-preprocess.ipf](#).



## Chapter 10

# Data Structure Documentation

### 10.1 DoniachSunjicStruct Struct Reference

#### Data Fields

- wave [pw](#)
- wave [yw](#)
- wave [xw](#)
- variable [precision](#)
- variable [oversampling](#)
- wave [xdw](#)
- wave [model](#)
- wave [broadening](#)
- wave [convolution](#)

#### 10.1.1 Detailed Description

Definition at line 191 of file `pearl-fitfuncs.ipf`.

#### 10.1.2 Field Documentation

##### 10.1.2.1 wave DoniachSunjicStruct::broadening

Definition at line 207 of file `pearl-fitfuncs.ipf`.

##### 10.1.2.2 wave DoniachSunjicStruct::convolution

Definition at line 208 of file `pearl-fitfuncs.ipf`.

##### 10.1.2.3 wave DoniachSunjicStruct::model

Definition at line 206 of file `pearl-fitfuncs.ipf`.

##### 10.1.2.4 variable DoniachSunjicStruct::oversampling

Definition at line 201 of file `pearl-fitfuncs.ipf`.

#### 10.1.2.5 variable `DoniachSunjicStruct::precision`

Definition at line 200 of file `pearl-fitfuncs.ipf`.

#### 10.1.2.6 wave `DoniachSunjicStruct::pw`

Definition at line 191 of file `pearl-fitfuncs.ipf`.

#### 10.1.2.7 wave `DoniachSunjicStruct::xdw`

Definition at line 205 of file `pearl-fitfuncs.ipf`.

#### 10.1.2.8 wave `DoniachSunjicStruct::xw`

Definition at line 197 of file `pearl-fitfuncs.ipf`.

#### 10.1.2.9 wave `DoniachSunjicStruct::yw`

Definition at line 196 of file `pearl-fitfuncs.ipf`.

The documentation for this struct was generated from the following file:

- [pearl-fitfuncs.ipf](#)

## 10.2 `errorCode` Struct Reference

from `matrixfilereader` help

### Data Fields

- int32 [SUCCESS](#)
- int32 [UNKNOWN\\_ERROR](#)
- int32 [ALREADY\\_FILE\\_OPEN](#)
- int32 [EMPTY\\_RESULTFILE](#)
- int32 [FILE\\_NOT\\_READABLE](#)
- int32 [NO\\_NEW\\_BRICKLETS](#)
- int32 [WRONG\\_PARAMETER](#)
- int32 [INTERNAL\\_ERROR\\_CONVERTING\\_DATA](#)
- int32 [NO\\_FILE\\_OPEN](#)
- int32 [INVALID\\_RANGE](#)
- int32 [WAVE\\_EXIST](#)

### 10.2.1 Detailed Description

from `matrixfilereader` help

Definition at line 197 of file `pearl-matrix-import.ipf`.

### 10.2.2 Field Documentation

#### 10.2.2.1 int32 errorCode::ALREADY\_FILE\_OPEN

Definition at line 200 of file pearl-matrix-import.ipf.

#### 10.2.2.2 int32 errorCode::EMPTY\_RESULTFILE

Definition at line 201 of file pearl-matrix-import.ipf.

#### 10.2.2.3 int32 errorCode::FILE\_NOT\_READABLE

Definition at line 202 of file pearl-matrix-import.ipf.

#### 10.2.2.4 int32 errorCode::INTERNAL\_ERROR\_CONVERTING\_DATA

Definition at line 205 of file pearl-matrix-import.ipf.

#### 10.2.2.5 int32 errorCode::INVALID\_RANGE

Definition at line 207 of file pearl-matrix-import.ipf.

#### 10.2.2.6 int32 errorCode::NO\_FILE\_OPEN

Definition at line 206 of file pearl-matrix-import.ipf.

#### 10.2.2.7 int32 errorCode::NO\_NEW\_BRICKLETS

Definition at line 203 of file pearl-matrix-import.ipf.

#### 10.2.2.8 int32 errorCode::SUCCESS

Definition at line 197 of file pearl-matrix-import.ipf.

#### 10.2.2.9 int32 errorCode::UNKNOWN\_ERROR

Definition at line 199 of file pearl-matrix-import.ipf.

#### 10.2.2.10 int32 errorCode::WAVE\_EXIST

Definition at line 208 of file pearl-matrix-import.ipf.

#### 10.2.2.11 int32 errorCode::WRONG\_PARAMETER

Definition at line 204 of file pearl-matrix-import.ipf.

The documentation for this struct was generated from the following file:

- [pearl-matrix-import.ipf](#)



# Chapter 11

## File Documentation

### 11.1 fermi-edge-analysis.ipf File Reference

tools for analysing the Fermi edge measured by the Scienta EW4000 analyser.

```
#include "pearl-area-profiles"
```

#### Functions

- variable [analyse\\_curved\\_edge](#) (wave data)
- variable [record\\_results](#) (variable index)
- variable [integrate\\_curved\\_edge](#) (wave data, wave data\_sig)
- variable [slit\\_correction](#) (wave data, wave data\_out, variable epass)
- threadsafe variable [FermiFuncLinDOS2D\\_corr](#) (variable w, threadsafe x, wave y)
- threadsafe variable [slit\\_shift](#) (variable ypix, variable epass)
- variable [show\\_shift](#) (wave data)
- variable [analyser\\_energy\\_resolution](#) (variable epass, variable slit)

*calculate the energy resolution of the analyser*

#### Variables

- static const variable [mcp\\_radius\\_pix](#) = 555  
*MCP radius seen by the camera in pixels.*
- static const variable [mcp\\_radius\\_mm](#) = 20  
*physical size (radius) of the MCP in mm*
- static const variable [hemi\\_radius\\_mm](#) = 200  
*physical size (radius) of the hemisphere in mm*
- static const variable [mcp\\_radius\\_epass](#) = 0.04  
*energy range imaged on MCP relative to the pass energy*

#### 11.1.1 Detailed Description

tools for analysing the Fermi edge measured by the Scienta EW4000 analyser.

proposed procedure

- angular normalization

- fit curved fermi function
- calculate corrected energy coordinates and map to single independent variable
- fit normal fermi function

#### Author

matthias muntwiler, [matthias.muntwiler@psi.ch](mailto:matthias.muntwiler@psi.ch)  
thomas diene

#### Copyright

2013-15 Paul Scherrer Institut  
Licensed under the Apache License, Version 2.0 (the "License");  
you may not use this file except in compliance with the License.  
You may obtain a copy of the License at <http://www.apache.org/licenses/LICENSE-2.0>

### 11.1.2 Function Documentation

#### 11.1.2.1 variable `analyse_curved_edge` ( *wave data* )

Definition at line 27 of file `fermi-edge-analysis.ipf`.

#### 11.1.2.2 variable `analyser_energy_resolution` ( *variable epass*, *variable slit* )

calculate the energy resolution of the analyser

##### Parameters

<i>epass</i>	pass energy in eV
<i>slit</i>	analyser entrance slit in mm

##### Returns

energy resolution (FWHM)

Definition at line 306 of file `fermi-edge-analysis.ipf`.

#### 11.1.2.3 threadsafe variable `FermiFuncLinDOS2D_corr` ( *variable w*, *threadsafe x*, *wave y* )

Definition at line 216 of file `fermi-edge-analysis.ipf`.

#### 11.1.2.4 variable `integrate_curved_edge` ( *wave data*, *wave data\_sig* )

Definition at line 116 of file `fermi-edge-analysis.ipf`.

#### 11.1.2.5 variable `record_results` ( *variable index* )

Definition at line 98 of file `fermi-edge-analysis.ipf`.

#### 11.1.2.6 variable `show_shift` ( *wave data* )

Definition at line 284 of file `fermi-edge-analysis.ipf`.

#### 11.1.2.7 variable `slit_correction` ( *wave data*, *wave data\_out*, variable *epass* )

Definition at line 173 of file `fermi-edge-analysis.ipf`.

#### 11.1.2.8 threadsafe variable `slit_shift` ( variable *ypix*, variable *epass* )

Definition at line 267 of file `fermi-edge-analysis.ipf`.

### 11.1.3 Variable Documentation

#### 11.1.3.1 `const variable hemi_radius_mm = 200` [static]

physical size (radius) of the hemisphere in mm

Definition at line 263 of file `fermi-edge-analysis.ipf`.

#### 11.1.3.2 `const variable mcp_radius_epass = 0.04` [static]

energy range imaged on MCP relative to the pass energy

Definition at line 265 of file `fermi-edge-analysis.ipf`.

#### 11.1.3.3 `const variable mcp_radius_mm = 20` [static]

physical size (radius) of the MCP in mm

Definition at line 261 of file `fermi-edge-analysis.ipf`.

#### 11.1.3.4 `const variable mcp_radius_pix = 555` [static]

MCP radius seen by the camera in pixels.

Definition at line 259 of file `fermi-edge-analysis.ipf`.

## 11.2 mainpage.dox File Reference

### 11.3 `pearl-anglescan-process.ipf` File Reference

processing and holographic mapping of angle scanned XPD data.

```
#include "pearl-vector-operations"
#include "pearl-polar-coordinates"
#include <New Polar Graphs>
```

#### Namespaces

- [PearlAnglescanProcess](#)

*processing and holographic mapping of angle scanned XPD data.*

## Functions

- variable [strip\\_delete\\_frames](#) (wave strip, variable qlo, variable qhi, wave theta, wave tilt, wave phi)  
*delete a contiguous range of frames from a strip.*
- variable [normalize\\_strip\\_x](#) (wave strip, variable smooth\_method=defaultValue, variable smooth\_factor=defaultValue, variable check=defaultValue)  
*divide the strip by the average X distribution.*
- variable [normalize\\_strip\\_phi](#) (wave strip, wave theta, wave phi, variable theta\_offset=defaultValue, variable theta\_range=defaultValue, variable check=defaultValue)  
*divide the strip by a sine function in phi (wobble correction).*
- variable [normalize\\_strip\\_theta](#) (wave strip, wave theta, variable theta\_offset=defaultValue, variable smooth\_method=defaultValue, variable smooth\_factor=defaultValue, variable check=defaultValue)  
*divide the strip by the average polar distribution.*
- variable [normalize\\_strip\\_2d](#) (wave strip, wave theta, variable theta\_offset=defaultValue, variable smooth\_method=defaultValue, variable smooth\_factor=defaultValue, variable check=defaultValue)  
*divide the strip by a two-dimensional normalization function.*
- variable [crop\\_strip](#) (wave strip, variable xlo, variable xhi)  
*crop a strip at the sides.*
- variable [pizza\\_service](#) (wave data, string nickname, variable theta\_offset, variable tilt\_offset, variable phi\_offset, variable npolar=defaultValue, variable nograph=defaultValue, variable folding=defaultValue, variable xpdplot=defaultValue)  
*create a pizza plot from a measured (energy-integrated) data strip*
- variable [pizza\\_service\\_2](#) (wave data, string nickname, wave m\_theta, wave m\_tilt, wave m\_phi, variable npolar=defaultValue, variable nograph=defaultValue, variable folding=defaultValue, variable xpdplot=defaultValue)  
*create a pizza plot from a measured (energy-integrated) data strip*
- variable [show\\_analyser\\_line](#) (variable theta, variable tilt, variable phi, variable theta\_offset, variable tilt\_offset, variable phi\_offset, variable npolar=defaultValue, variable nograph=defaultValue, variable xpdplot=defaultValue)  
*calculate and display the line seen by the analyser for a specific emission angle*
- variable [convert\\_angles\\_ttpd2polar](#) (wave theta, wave tilt, wave phi, wave data, wave polar, wave azi)  
*convert angles from TTPA (theta-tilt-phi-analyser) scheme to polar coordinates.*
- variable [convert\\_angles\\_ttpa2polar](#) (wave theta, wave tilt, wave phi, wave analyser, wave polar, wave azi)  
*convert angles from TTPA (theta-tilt-phi-analyser) scheme to polar coordinates.*
- static variable [line\\_average](#) (wave source, wave dest)
- static variable [calc\\_nth](#) (variable Theta\_st, variable Theta\_in, variable th, variable Phi\_ran, variable Phi\_ref, string Holomode)  
*calculate the number of phis for a given theta*
- static variable [calc\\_phi\\_step](#) (variable Theta\_in, variable th, variable Theta\_st, variable Phi\_ran, variable Phi\_ref, string Holomode)  
*calculate delta-phi for a given theta*
- static variable [Calc\\_The\\_step](#) (variable th, variable Theta\_st, string Holomode)  
*calculate delta-theta for a given theta*
- static variable [CalcN\\_Theta](#) (string HoloMode, variable Theta\_in, variable Theta\_ran, variable Theta\_st)  
*calculate the number of thetas for a pattern*
- variable [make\\_hemi\\_grid](#) (variable npol, string nickname, variable xpdplot=defaultValue)  
*create a hemispherical, constant solid angle grid*
- string [get\\_hemi\\_nickname](#) (wave w)  
*finds the nick name given any hemi wave*
- string [get\\_hemi\\_prefix](#) (wave w)  
*finds the prefix given any hemi wave*
- dfr [find\\_hemi\\_data](#) (string nickname, string \*prefix, string \*intwave)  
*finds the folder, prefix and name of holo waves given their nick name*

- variable `clear_hemi_grid` (string nickname)  
*clear a hemispherical scan grid*
- variable `duplicate_hemi_scan` (string source\_nickname, dfref dest\_folder, string dest\_nickname, variable xpdplot=defaultValue)  
*duplicate a hemispherical scan dataset.*
- variable `rotate_hemi_scan` (string nickname, variable angle)  
*azimuthally rotate a hemispherical scan dataset.*
- string `display_hemi_scan` (string nickname, variable projection=defaultValue, variable graphtype=default↵  
Value, variable do\_ticks=defaultValue, variable do\_grids=defaultValue, string graphname=defaultValue)  
*display a plot of a hemispherical angle scan.*
- static string `display_polar_graph` (string graphname, variable angle\_offset=defaultValue, variable do\_↵  
ticks=defaultValue)  
*displays an empty polar graph*
- static string `draw_hemi_axes` (string graphname, variable do\_grids=defaultValue)  
*draw polar and azimuthal grids in an existing polar graph.*
- variable `draw_diffraction_cone` (string graphname, string groupname, variable theta\_axis, variable theta\_↵  
inner, variable phi)  
*draw the circle of a diffraction cone in a stereographic polar graph.*
- string `display_scanlines` (string nickname, variable alpha\_lo, variable alpha\_hi, wave m\_theta, wave m\_tilt,  
wave m\_phi, variable folding=defaultValue, variable projection=defaultValue)  
*display a polar graph with lines indicating the angles covered by an angle scan.*
- threadsafe variable `calc_graph_radius` (variable polar, variable projection=defaultValue)  
*calculate the projected polar angle*
- threadsafe variable `calc_graph_polar` (variable x, variable y, variable projection=defaultValue)  
*calculate polar angle from Cartesian coordinate*
- threadsafe variable `calc_graph_azi` (variable x, variable y, variable projection=defaultValue, variable zero\_↵  
Angle=defaultValue)  
*calculate azimuthal angle from Cartesian coordinate*
- static variable `update_polar_info` (string graphname)  
*update the angles info based on cursors A and B of a given polar graph window*
- static variable `polar_graph_hook` (WMWinHookStruct \*s)  
*polar graph window hook*
- variable `set_polar_graph_cursor` (string nickname, string cursorname, variable polar\_angle, variable azim\_↵  
angle, string graphname=defaultValue)
- variable `hemi_add_anglescan` (string nickname, wave values, wave polar, wave azi, wave weights=default\_↵  
Value)  
*add an arbitrary angle scan to a hemispherical scan grid.*
- variable `hemi_add_aziscan` (string nickname, wave values, variable polar, wave azi, wave weights=default\_↵  
Value)  
*add an azimuthal scan to a hemispherical scan grid.*
- variable `interpolate_hemi_scan` (string nickname)  
*interpolate a hemispherical scan onto a rectangular grid*
- variable `quick_pizza_image` (wave data, string nickname, variable theta\_offset, variable tilt\_offset, variable  
phi\_offset, variable npolar=defaultValue, variable nograph=defaultValue, variable folding=defaultValue)  
*map angle scan data onto a rectangular grid in stereographic projection*
- variable `save_hemi_scan` (string nickname, string pathname, string filename)  
*save a hemispherical scan to an Igor text file*
- variable `load_hemi_scan` (string nickname, string pathname, string filename)  
*load a hemispherical scan from an Igor text file*
- variable `import_tpi_scan` (string nickname, wave theta, wave phi, wave intensity, variable folding=defaultValue,  
variable npolar=defaultValue, variable nograph=defaultValue, variable xpdplot=defaultValue)  
*import a hemispherical scan from theta-phi-intensity waves and display it*

- variable `trim_hemi_scan` (string nickname, variable `theta_max`)  
*trim a hemispherical scan at grazing angle*
- wave `hemi_polar_cut` (string nickname, variable `azim`)  
*extract a polar cut from a hemispherical scan.*
- wave `hemi_azi_cut` (string nickname, variable `pol`)  
*extract an azimuthal cut from a hemispherical scan*
- static variable `check_contrast` (wave values, variable `pcmin`, variable `pcmax`, variable `*vmin`, variable `*vmax`)
- variable `set_contrast` (variable `pcmin`, variable `pcmax`, string `graphname=defaultValue`, string `colortable=defaultValue`)  
*set the pseudocolor contrast by percentile.*

## Variables

- const variable `kProjDist` = 0
- const variable `kProjStereo` = 1
- const variable `kProjArea` = 2
- const variable `kProjGnom` = 3
- const variable `kProjOrtho` = 4
- static const variable `kProjScaleDist` = 2
- static const variable `kProjScaleStereo` = 2
- static const variable `kProjScaleArea` = 2
- static const variable `kProjScaleGnom` = 0.06744519021
- static const variable `kProjScaleOrtho` = 2

### 11.3.1 Detailed Description

processing and holographic mapping of angle scanned XPD data.

the functions in this file map angle scanned data measured at PEARL onto a hemispherical angle grid which is compatible with XPDplot. the resulting data are in a canonical polar coordinate system (normal emission  $\leftarrow$   $\rightarrow$  polar angle = 0, azimuthal axis right-handed) which is anchored in the sample surface. the orientation of polar graphs (phi = 0 at 3 o'clock) created by this procedure corresponds to a top view of the sample surface at normal emission, and the handle of the sample plate pointing to the left (phi = 180). this is the canonical orientation of a spherical coordinate system where phi = 0 corresponds to the positive part of the x axis.

#### Note

the orientation of the sample coordinate system has changed in version 1.6. the change was necessary for compatibility with other data analysis software and calculation programs.

data imported with version 1.5 and earlier, must be offset by 180 deg in phi to be compatible with the new version. data imported and displayed by the same code version will give the same picture but with different azimuthal axis. the new graph functions shows a warning if they are applied to code imported with earlier versions.

the measurement geometry is hard-coded but may be parametrized in the future. the theta rotation axis is perpendicular to the scattering plane. the angle dispersive axis of the analyser is parallel to the theta rotation axis. the tilt rotation axis is in the scattering plane. it rotates with theta. at normal emission it is perpendicular to the axis of the lens stack of the analyser. the phi rotation axis corresponds to the surface normal of the sample. it rotates with theta and with tilt. at normal emission it is parallel to the axis of the lens stack of the analyser.

coordinate transformations: (to be revised - v1.6)

```
*      theta_sample = theta_manipulator - theta_offset
*      phi_sample = phi_manipulator - phi_offset
*
```

valid for `theta_manipulator` = normal emission only: (to be revised - v1.6)

```
*      theta_sample = | -(tilt_manipulator - tilt_offset) |
*      phi_sample = 270 if tilt_manipulator - tilt_offset > 0
*      phi_sample = 90 if tilt_manipulator - tilt_offset < 0
*
```

#### Author

matthias muntwiler, [matthias.muntwiler@psi.ch](mailto:matthias.muntwiler@psi.ch)

#### Copyright

2013-17 Paul Scherrer Institut

Licensed under the Apache License, Version 2.0 (the "License");

you may not use this file except in compliance with the License.

You may obtain a copy of the License at <http://www.apache.org/licenses/LICENSE-2.0>

#### Version

1.8 canonical orientation of spherical coordinate system.

### 11.3.2 Function Documentation

**11.3.2.1** `threadsafe variable calc_graph_azi ( variable x, variable y, variable projection = defaultValue, variable zeroAngle = defaultValue )`

calculate azimuthal angle from Cartesian coordinate

#### Parameters

<i>x,y</i>	projected Cartesian coordinate
<i>projection</i>	mapping function from polar to cartesian coordinates. all supported projections are azimuthal, they have no effect on the azimuthal coordinate. see <a href="#">Projections</a> for details. <ul style="list-style-type: none"> <li>• kProjDist = 0 azimuthal equidistant</li> <li>• kProjStereo = 1 stereographic (default)</li> <li>• kProjArea = 2 azimuthal equal-area</li> <li>• kProjGnom = 3 gnomonic (0 &lt;= polar &lt; 90)</li> <li>• kProjOrtho = 4 orthographic</li> </ul>
<i>zeroAngle</i>	zeroAngleWhere parameter of polar graphs <ul style="list-style-type: none"> <li>• 0 (default) zero is at the 3 o'clock position</li> <li>• 180 zero is at the 9 o'clock position</li> <li>• other values not tested</li> </ul>

#### Returns

polar angle in degrees

Definition at line 2129 of file pearl-anglescan-process.ipf.

### 11.3.2.2 threadsafe variable `calc_graph_polar` ( variable *x*, variable *y*, variable *projection* = defaultValue )

calculate polar angle from Cartesian coordinate

this is the reverse mapping to [calc\\_graph\\_radius\(\)](#)

#### Parameters

<i>x,y</i>	projected Cartesian coordinate
<i>projection</i>	mapping function from polar to cartesian coordinates. see <a href="#">Projections</a> for details. <ul style="list-style-type: none"> <li>• kProjDist = 0 azimuthal equidistant</li> <li>• kProjStereo = 1 stereographic (default)</li> <li>• kProjArea = 2 azimuthal equal-area</li> <li>• kProjGnom = 3 gnomonic (<math>0 \leq \text{polar} &lt; 90</math>)</li> <li>• kProjOrtho = 4 orthographic</li> </ul>

#### Returns

polar angle in degrees

Definition at line 2076 of file `pearl-anglescan-process.ipf`.

### 11.3.2.3 threadsafe variable `calc_graph_radius` ( variable *polar*, variable *projection* = defaultValue )

calculate the projected polar angle

#### Parameters

<i>polar</i>	polar angle in degrees
<i>projection</i>	mapping function from polar to cartesian coordinates. see <a href="#">Projections</a> for details. <ul style="list-style-type: none"> <li>• kProjDist = 0 azimuthal equidistant</li> <li>• kProjStereo = 1 stereographic (default)</li> <li>• kProjArea = 2 azimuthal equal-area</li> <li>• kProjGnom = 3 gnomonic (<math>0 \leq \text{polar} &lt; 90</math>)</li> <li>• kProjOrtho = 4 orthographic</li> </ul>

#### Returns

projected radius. the radius is scaled such that grazing emission maps to 2.

Definition at line 2031 of file `pearl-anglescan-process.ipf`.

### 11.3.2.4 static variable `calc_nth` ( variable *Theta\_st*, variable *Theta\_in*, variable *th*, variable *Phi\_ran*, variable *Phi\_ref*, string *Holomode* ) [static]

calculate the number of phis for a given theta

adapted from XPDplot 8.03

Definition at line 935 of file pearl-anglescan-process.ipf.

**11.3.2.5** `static variable calc_phi_step ( variable Theta_in, variable th, variable Theta_st, variable Phi_ran, variable Phi_ref, string Holomode ) [static]`

calculate delta-phi for a given theta

adapted from XPDplot 8.03

Definition at line 964 of file pearl-anglescan-process.ipf.

**11.3.2.6** `static variable Calc_The_step ( variable th, variable Theta_st, string Holomode ) [static]`

calculate delta-theta for a given theta

adapted from XPDplot 8.03

Definition at line 1008 of file pearl-anglescan-process.ipf.

**11.3.2.7** `static variable CalcN_Theta ( string HoloMode, variable Theta_in, variable Theta_ran, variable Theta_st ) [static]`

calculate the number of thetas for a pattern

adapted from XPDplot 8.03

Definition at line 1039 of file pearl-anglescan-process.ipf.

**11.3.2.8** `static variable check_contrast ( wave values, variable pcmin, variable pcmax, variable * vmin, variable * vmax ) [static]`

Definition at line 2925 of file pearl-anglescan-process.ipf.

**11.3.2.9** `variable clear_hemi_grid ( string nickname )`

clear a hemispherical scan grid

values and weights waves are set to zero. the intensity wave is set to NaN.

#### Parameters

<i>nickname</i>	folder name or name prefix of holo waves. may be empty.
-----------------	---

Definition at line 1318 of file pearl-anglescan-process.ipf.

**11.3.2.10** `variable convert_angles_ttpa2polar ( wave theta, wave tilt, wave phi, wave analyser, wave polar, wave azi )`

convert angles from TTPA (theta-tilt-phi-analyser) scheme to polar coordinates.

the angles are in the manipulator coordinate system.

#### Parameters

in	<i>theta</i>	offset-corrected theta angle, normal emission = 0, grazing emission = 90. one dimensional wave.
in	<i>tilt</i>	offset-corrected tilt angle, normal emission = 0 same dimension size and scale as theta

## Parameters

in	<i>phi</i>	phi angle, range $-360 < \phi < +360$ offset correction is optional as long as the angles lie in the accepted range. same dimension size and scale as theta
in	<i>analyser</i>	analyser angle scale corresponding to the slices scale of Scienta. one dimensional wave. this values are constant regardless of manipulator angle.
out	<i>polar</i>	wave to receive the polar coordinates.
out	<i>azi</i>	wave to receive the azimuthal coordinates.

for the output parameters polar and azi, you need to pass in existing numeric waves. dimension size does not matter, the waves are redimensioned by the function so that they have the same dimensions as the intensity data set. X dimension = analyser scale, Y dimension = manipulator scan.

Definition at line 874 of file pearl-anglescan-process.ipf.

#### 11.3.2.11 variable `convert_angles_ttpd2polar ( wave theta, wave tilt, wave phi, wave data, wave polar, wave azi )`

convert angles from TTPA (theta-tilt-phi-analyser) scheme to polar coordinates.

similar to `convert_angles_ttpa2polar()` but reads the analyser angles from the X scale of data

Definition at line 835 of file pearl-anglescan-process.ipf.

#### 11.3.2.12 variable `crop_strip ( wave strip, variable xlo, variable xhi )`

crop a strip at the sides.

the strip is cropped in place, data outside the region of interest is lost.

## Parameters

in, out	<i>strip</i>	2D data, X-axis = analyser angle, Y-axis = arbitrary manipulator scan
in	<i>xlo</i>	lowest analyser angle to keep (will be rounded to nearest existing point)
in	<i>xhi</i>	highest analyser angle to keep (will be rounded to nearest existing point)

## Remarks

cropping should be done after smoothing and normalization operations to reduce artefacts.

Definition at line 495 of file pearl-anglescan-process.ipf.

#### 11.3.2.13 string `display_hemi_scan ( string nickname, variable projection = defaultValue, variable graphtype = defaultValue, variable do_ticks = defaultValue, variable do_grids = defaultValue, string graphname = defaultValue )`

display a plot of a hemispherical angle scan.

the scan data must exist in the current data folder. azimuth = 0 should be at 9 o'clock. then the orientation is the same as the sample at normal emission and phi = 0, the handle of the sample plate pointing to the left.

## Parameters

<i>nickname</i>	name prefix of holo waves. may be empty.
-----------------	--

## Parameters

<i>projection</i>	mapping function from polar to cartesian coordinates <ul style="list-style-type: none"> <li>• 0 linear</li> <li>• 1 stereographic (default)</li> <li>• 2 azimuthal</li> <li>• 3 gnomonic (<math>0 \leq \text{polar} &lt; 90</math>).</li> </ul>
<i>graphtype</i>	type of graph <ul style="list-style-type: none"> <li>• 1 (pol, az) trace in Igor "New Polar" (default).</li> <li>• 2 XPDplot (reserved, not implemented).</li> <li>• 3 matrix in Igor "New Polar". the matrix wave is a 2D wave with X and Y scaling corresponding to the selected projection. matrix waves can be created by <a href="#">interpolate_hemi_scan()</a>. note: the pol and az waves are required as well.</li> </ul>
<i>do_ticks</i>	select which ticks to draw. value must be the arithmetic OR of all selected items. default: 3 <ul style="list-style-type: none"> <li>• 0 none</li> <li>• 1 major azimuthal</li> <li>• 2 minor azimuthal S</li> </ul>
<i>do_grids</i>	select which grids to draw. value must be the arithmetic OR of all selected items. default: 3 <ul style="list-style-type: none"> <li>• 0 none</li> <li>• 1 radius at 0 and 90 degree azimuth</li> <li>• 2 circle at 30 and 60 degree polar</li> </ul>
<i>graphname</i>	name of graph window. default: nickname if empty, a default name is assigned. if a window with this name is existing, the function brings it to the front, and does nothing else.

## Returns

the name of the graph window

Definition at line 1528 of file pearl-anglescan-process.ipf.

**11.3.2.14** `static string display_polar_graph ( string graphname, variable angle_offset = defaultvalue, variable do_ticks = defaultvalue ) [static]`

displays an empty polar graph

the graph is drawn using Wavemetrics "New Polar Graphs.ipf".

initially the graph is empty. hemispherical scans are displayed by adding a trace that covers the whole plot area, and setting the trace color to a function of the intensity. traces are added by calling `WMPolarAppendTrace`.

the following items of the graph > packages menu might be useful:

- modify polar graph
- color table control
- show polar cursors

- polar graph legend

parameters can be changed programmatically as shown in the code of this function. after programmatic parameter changes, call `WMPolarAxesRedrawGraphNow(graphname)`.

#### Parameters

<i>graphname</i>	requested name of new graph window. if empty, a default name is assigned. if a window with this name is existing, the function brings it to the front, and does nothing else.
<i>angle_offset</i>	azimuth (on screen) where angle 0 is plotted ( <code>zeroAngleWhere</code> parameter of polar graphs). starting with version 1.6, the default is 0. for hemi grids created with earlier versions, it should be set to 180 for correct orientation.
<i>do_ticks</i>	select which ticks to draw. value must be the arithmetic OR of all selected items. default: 3 <ul style="list-style-type: none"> <li>• 0 none</li> <li>• 1 major azimuthal</li> <li>• 2 minor azimuthal</li> </ul>

#### Returns

the name of the graph window.

#### Version

1.7 interface change: the trace drawing code is moved to `display_hemi_scan`, so that this function can be reused by other graph types, e.g. `display_scanlines`.

Definition at line 1665 of file `pearl-anglescan-process.ipf`.

**11.3.2.15** `string display_scanlines ( string nickname, variable alpha_lo, variable alpha_hi, wave m_theta, wave m_tilt, wave m_phi, variable folding = defaultValue, variable projection = defaultValue )`

display a polar graph with lines indicating the angles covered by an angle scan.

#### Parameters

<i>nickname</i>	nick name for output data. this will become the name of a child folder containing the output.
<i>alpha_lo</i>	low limit of the analyser angle.
<i>alpha_hi</i>	high limit of the analyser angle.
<i>m_theta</i>	manipulator theta angles, 0 = normal emission. size = <code>dimsize(data, 1)</code>
<i>m_tilt</i>	manipulator tilt angles, 0 = normal emission. size = <code>dimsize(data, 1)</code>
<i>m_phi</i>	manipulator phi angles, 0 = azimuthal origin. size = <code>dimsize(data, 1)</code>
<i>folding</i>	rotational averaging, default = 1
<i>projection</i>	mapping function from polar to cartesian coordinates. see <a href="#">calc_graph_radius()</a> .

#### Remarks

this function is extremely slow.

Definition at line 1916 of file `pearl-anglescan-process.ipf`.

11.3.2.16 `variable draw_diffraction_cone ( string graphname, string groupname, variable theta_axis, variable theta_inner, variable phi )`

draw the circle of a diffraction cone in a stereographic polar graph.

the diffraction cone consists of a circle marking the diffraction ring, and a dot marking the axis. the cone is drawn as a group of draw objects on the UserFront layer. the objects can be edited interactively.

#### Parameters

<i>graphname</i>	name of graph window (not implemented yet).
<i>groupname</i>	name of a drawing group. if the group exists (from a previous cone) it is replaced. if the group doesn't exist, a new one is created.
<i>theta_axis</i>	polar angle of the cone axis in degrees.
<i>theta_inner</i>	polar angle of the innermost point of the circle in degrees.
<i>phi</i>	azimuthal angle of the cone axis in degrees.

#### Warning

EXPERIMENTAL! this function is under development. the interface and behaviour of this function may change significantly in future versions.

Definition at line 1852 of file pearl-anglescan-process.ipf.

11.3.2.17 `static string draw_hemi_axes ( string graphname, variable do_grids = defaultValue ) [static]`

draw polar and azimuthal grids in an existing polar graph.

the function adds the following draw objects to a polar graph:

- concentric circles at polar angles 0, 30, and 60 degrees with labels.
- radial axes at 0 and 90 degree azimuth.

the objects are added to the ProgFront drawing layer and will appear in front of the data trace. in interactive drawing mode, you can select the active drawing layer by clicking the tree icon while holding the Alt key.

the graph must have been created by [display\\_polar\\_graph\(\)](#). the function reads the projection mode from the window user data "projection".

#### Parameters

<i>graphname</i>	name of graph window.
<i>do_grids</i>	select which optional grids to draw. value must be the arithmetic OR of all selected items. default: 3 <ul style="list-style-type: none"> <li>• 0 none</li> <li>• 1 radius at 0 and 90 degree azimuth</li> <li>• 2 circle at 30 and 60 degree polar</li> </ul>

#### Warning

EXPERIMENTAL! this function is under development. the interface and behaviour of this function may change significantly in future versions.

Definition at line 1780 of file pearl-anglescan-process.ipf.

**11.3.2.18** `variable duplicate_hemi_scan ( string source_nickname, dfref dest_folder, string dest_nickname, variable xpdplot = defaultvalue )`

duplicate a hemispherical scan dataset.

this function works only for hemi scans created by [make\\_hemi\\_grid\(\)](#) (or compatible functions). the angle grid is recreated rather than copied point-by-point. the new dataset is independent from the original one.

if the version of the source dataset is pre 1.6, it is converted to version 1.6.

#### Parameters

<i>source_nickname</i>	name prefix for waves. source data must be in current data folder.
<i>dest_folder</i>	destination folder. folder must exist.
<i>dest_nickname</i>	name prefix for destination waves. must be unique in the current data folder. otherwise existing waves get overwritten. may be empty.
<i>xpdplot</i>	XPDplot compatibility <ul style="list-style-type: none"> <li>• 0 (default) create the data structures required by this module</li> <li>• 1 create additional waves and notebook required by XPDplot</li> </ul>

Definition at line 1365 of file pearl-anglescan-process.ipf.

**11.3.2.19** `dfr find_hemi_data ( string nickname, string * prefix, string * intwave )`

finds the folder, prefix and name of holo waves given their nick name

the function looks for holo waves in the following order:

1. if nickname is empty, check for prefix-less waves in current folder.
2. if nickname is the name of a child folder in the current data folder, clear the (prefix-less) waves in the child folder.
3. nickname is prefix of waves in current folder.
4. nickname is prefix of waves in root folder.

#### Parameters

in	<i>nickname</i>	folder name or name prefix of holo waves. may be empty.
out	<i>prefix</i>	name prefix of waves. may be empty.
out	<i>intwave</i>	name of intensity/values wave

#### Returns

reference of the data folder which contains the waves

Definition at line 1284 of file pearl-anglescan-process.ipf.

**11.3.2.20** `string get_hemi_nickname ( wave w )`

finds the nick name given any hemi wave

the nick name is either the name of a child folder in the current data folder (PEARL specification), or a prefix of the hemi wave names (XPDplot specification).

**Returns**

the nick name

Definition at line 1228 of file pearl-anglescan-process.ipf.

**11.3.2.21 string get\_hemi\_prefix ( wave w )**

finds the prefix given any hemi wave

the prefix is the part of the wave name before the first underscore. the prefix is used by XPDplot where it is identical to the nick name. the prefix is empty in the PEARL specification.

**Returns**

the prefix

Definition at line 1253 of file pearl-anglescan-process.ipf.

**11.3.2.22 variable hemi\_add\_anglescan ( string nickname, wave values, wave polar, wave azi, wave weights = defaultValue )**

add an arbitrary angle scan to a hemispherical scan grid.

the hemi grid must have been created in the current data folder by the make\_hemi\_grid function. the function determines the bin size at the given polar angle, and adds all data points which fall into a bin. a point which lies exactly on the upper boundary falls into the next bin. this function does not clear previous values before adding new data. values are added to the \_tot wave, weights to the \_wt wave. the intensity (\_i) wave is calculated as \_tot / \_wt.

Definition at line 2275 of file pearl-anglescan-process.ipf.

**11.3.2.23 variable hemi\_add\_aziscan ( string nickname, wave values, variable polar, wave azi, wave weights = defaultValue )**

add an azimuthal scan to a hemispherical scan grid.

the hemi grid must have been created in the current data folder by the make\_hemi\_grid function. the function determines the bin size at the given polar angle, and calculates the mean values of the data points which fall into a bin. a point which lies exactly on the upper boundary falls into the next bin.

Definition at line 2348 of file pearl-anglescan-process.ipf.

**11.3.2.24 wave hemi\_azi\_cut ( string nickname, variable pol )**

extract an azimuthal cut from a hemispherical scan

the function extracts all azimuthal angles that are present for the given polar angle.

the hemi grid must have been created in the current data folder by the make\_hemi\_grid function. correct ordering is required.

**Parameters**

<i>nickname</i>	name of the scan dataset. can be empty if no prefix is used. the dataset must be in the current datafolder.
<i>pol</i>	polar angle in degrees

**Returns**

reference of the created wave. the wave has the same name as the intensity wave of the dataset with the suffix "\_azi" and the azimuthal angle rounded to integer. it is created in the same datafolder as the original data.

Definition at line 2874 of file pearl-anglescan-process.ipf.

**11.3.2.25 wave hemi\_polar\_cut ( string *nickname*, variable *azim* )**

extract a polar cut from a hemispherical scan.

for each polar angle, the function first extracts all azimuthal angles. the intensity is then interpolated between the nearest neighbours of the given azimuth.

the hemi grid must have been created in the current data folder by the make\_hemi\_grid function. correct ordering is required.

**Parameters**

<i>nickname</i>	name of the scan dataset. can be empty if no prefix is used. the dataset must be in the current datafolder.
<i>azim</i>	azimuthal angle in degrees

**Returns**

reference of the created wave. the wave has the same name as the intensity wave of the dataset with the suffix "\_azi" and the azimuthal angle rounded to integer. it is created in the same datafolder as the original data.

Definition at line 2790 of file pearl-anglescan-process.ipf.

**11.3.2.26 variable import\_tpi\_scan ( string *nickname*, wave *theta*, wave *phi*, wave *intensity*, variable *folding* = defaultvalue, variable *npolar* = defaultvalue, variable *nograph* = defaultvalue, variable *xpdplot* = defaultvalue )**

import a hemispherical scan from theta-phi-intensity waves and display it

in the tpi format, the hemi scan data is represented by a triple of flat one-dimensional waves corresponding to the polar angle (theta), azimuthal angle (phi) and intensity. no specific sort order is required.

**Parameters**

<i>nickname</i>	nick name for output data <ul style="list-style-type: none"> <li>• in default mode, this will become the name of a child folder containing the output.</li> <li>• in XPDplot mode, this will become a prefix of the generated data in the root folder.</li> </ul>
<i>theta</i>	theta angles, 0 = normal emission.
<i>phi</i>	phi angles, 0 = azimuthal origin. size = dimsize(data, 1)
<i>intensity</i>	intensity wave, see requirements above.
<i>npolar</i>	number of polar angles, determines polar and azimuthal step size. default = 91 (1 degree steps)
<i>folding</i>	rotational averaging. example: 3 = average to 3-fold symmetry. default = 1.
<i>nograph</i>	display a new graph window? <ul style="list-style-type: none"> <li>• 0 (default) display a new polar graph</li> <li>• 1 don't display a new graph</li> </ul>

## Parameters

<i>xpdplot</i>	XPDplot compatibility <ul style="list-style-type: none"> <li>• 0 (default) create waves in child folder \$nickname</li> <li>• 1 create waves in root folder (compatible with XPDplot)</li> </ul>
----------------	--

Definition at line 2699 of file pearl-anglescan-process.ipf.

#### 11.3.2.27 variable interpolate\_hemi\_scan ( string *nickname* )

interpolate a hemispherical scan onto a rectangular grid

## Warning

experimental this function has been tested for one specific set of scan parameters. the interface and code may change at any time. the function depends on the *ster\_x* and *ster\_y* waves that are created by `display_hemi_scan`.

Definition at line 2449 of file pearl-anglescan-process.ipf.

#### 11.3.2.28 static variable line\_average ( wave *source*, wave *dest* ) [static]

Definition at line 917 of file pearl-anglescan-process.ipf.

#### 11.3.2.29 variable load\_hemi\_scan ( string *nickname*, string *pathname*, string *filename* )

load a hemispherical scan from an Igor text file

**Todo** function not implemented

Definition at line 2650 of file pearl-anglescan-process.ipf.

#### 11.3.2.30 variable make\_hemi\_grid ( variable *npol*, string *nickname*, variable *xpdplot* = defaultValue )

create a hemispherical, constant solid angle grid

all necessary waves are created in the current data folder with step size  $90 / (npol - 1)$

adapted from XPDplot 8.03

## Parameters

<i>npol</i>	number of polar angles, determines polar and azimuthal step size. recommended 91 for 1-degree steps.
<i>nickname</i>	name prefix for waves. nick name must be unique in the current data folder. otherwise existing waves get overwritten. may be empty.
<i>xpdplot</i>	XPDplot compatibility <ul style="list-style-type: none"> <li>• 0 (default) create the data structures required by this module</li> <li>• 1 create additional waves and notebook required by XPDplot</li> </ul>

Definition at line 1077 of file pearl-anglescan-process.ipf.

**11.3.2.31** `variable normalize_strip_2d ( wave strip, wave theta, variable theta_offset = defaultValue, variable smooth_method = defaultValue, variable smooth_factor = defaultValue, variable check = defaultValue )`

divide the strip by a two-dimensional normalization function.

#### Warning

experimental. this function is under development.

#### Parameters

<i>check</i>	enable output of intermediate results <ul style="list-style-type: none"> <li>• 0 (default) don't create additional waves</li> <li>• 1 create check waves in the current folder</li> <li>• 2 calculate check waves only, do not modify strip</li> </ul>
--------------	--

#### Returns

if check waves are enabled, the following waves are created (overwritten if existing):

- *check\_dist* average theta distribution
- *check\_smoo* smoothed distribution used to normalize the strip

Definition at line 435 of file pearl-anglescan-process.ipf.

**11.3.2.32** `variable normalize_strip_phi ( wave strip, wave theta, wave phi, variable theta_offset = defaultValue, variable theta_range = defaultValue, variable check = defaultValue )`

divide the strip by a sine function in phi (wobble correction).

the sine function is a curve fit to the intensity integrated over detector angle with a period of 360.

this normalization may be useful if the intensity varies with a 360 periodicity in the azimuthal angle, e.g. due to misalignment of the surface normal and the azimuthal rotation axis of the manipulator (wobble). note, however, that this function does not correct other effects of wobble such as angle shifts.

the strip is normalized in place, previous data is overwritten.

#### Parameters

in, out	<i>strip</i>	2D data, X-axis = analyser angle, Y-axis = arbitrary manipulator scan
in	<i>theta</i>	polar manipulator angle.
in	<i>phi</i>	azimuthal manipulator angle, arbitrary offset.
in	<i>theta_offset</i>	theta value corresponding to normal emission (default 0).
in	<i>theta_range</i>	maximum (offset corrected) theta to consider in the sine fit (default 10).
	<i>check</i>	enable output of intermediate results <ul style="list-style-type: none"> <li>• 0 (default) don't create additional waves</li> <li>• 1 create check waves in the current folder</li> <li>• 2 calculate check waves only, do not modify strip</li> </ul>

**Returns**

if check waves are enabled, the following waves are created (overwritten if existing):

- check\_dist average theta distribution
- check\_smoo smoothed distribution used to normalize the strip

Definition at line 270 of file pearl-anglescan-process.ipf.

**11.3.2.33** `variable normalize_strip_theta ( wave strip, wave theta, variable theta_offset = defaultValue, variable smooth_method = defaultValue, variable smooth_factor = defaultValue, variable check = defaultValue )`

divide the strip by the average polar distribution.

this is a simple way to remove the polar angle dependence. the strip is normalized in place, previous data is overwritten.

**Parameters**

in, out	<i>strip</i>	2D data, X-axis = analyser angle, Y-axis = arbitrary manipulator scan
in	<i>theta</i>	polar manipulator angle, 0 = normal emission, 90 = grazing emission
in	<i>theta_offset</i>	
in	<i>smooth_method</i>	smoothing method <ul style="list-style-type: none"> <li>• 0 none</li> <li>• 1 binomial (requires monotonic theta), see Igor's Smooth operation</li> <li>• 2 boxcar (requires monotonic theta), see Igor's Smooth operation</li> <li>• 3 polynomial fit per slice</li> <li>• 4 (default) Loess, see Igor's Loess operation</li> </ul> caution: binomial and boxcar smoothing are not aware of theta. this may give unpredictable results if theta is non-monotonic.
in	<i>smooth_factor</i>	smoothing parameter, depends on smooth_method <ul style="list-style-type: none"> <li>• binomial/boxcar: see Igor's Smooth operation</li> <li>• loess: see Igor's Loess operation, 0 &lt;= smooth_factor &lt;= 1, default 0.5</li> <li>• polynomial fit: polynomial degree, 1 = linear (default), 2 = quadratic</li> </ul>
	<i>check</i>	enable output of intermediate results <ul style="list-style-type: none"> <li>• 0 (default) don't create additional waves</li> <li>• 1 create check waves in the current folder</li> <li>• 2 calculate check waves only, do not modify strip</li> </ul>

**Returns**

if check waves are enabled, the following waves are created (overwritten if existing):

- check\_dist average theta distribution
- check\_smoo smoothed distribution used to normalize the strip

Definition at line 353 of file pearl-anglescan-process.ipf.

**11.3.2.34** `variable normalize_strip_x ( wave strip, variable smooth_method = defaultValue, variable smooth_factor = defaultValue, variable check = defaultValue )`

divide the strip by the average X distribution.

this is a simple way to remove the effect of the angle-dependence of the analyser transmission function. the strip is normalized in place, previous data is overwritten.

#### Parameters

<i>in, out</i>	<i>strip</i>	2D data, X-axis = analyser angle, Y-axis = arbitrary manipulator scan
<i>in</i>	<i>smooth_method</i>	smoothing method <ul style="list-style-type: none"> <li>• 0 none</li> <li>• 1 (default) binomial, see Igor's Smooth operation</li> <li>• 2 boxcar, see Igor's Smooth operation</li> <li>• 3 <a href="#">scientia_ang_transm()</a> function fit</li> <li>• 4 LOESS smoothing, see Igor's Loess operation</li> </ul>
<i>in</i>	<i>smooth_factor</i>	num parameter of Igor's Smooth operation. the default value depends on <i>smooth_method</i> . it is 0.5 for LOESS smoothing, 2 otherwise.
<i>in</i>	<i>check</i>	enable output of intermediate results <ul style="list-style-type: none"> <li>• 0 (default) don't create additional waves</li> <li>• 1 create check waves in the current folder</li> <li>• 2 calculate check waves only, do not modify strip</li> </ul>

#### Returns

if check waves are enabled, the following waves are created (overwritten if existing):

- *check\_dist* average X distribution
- *check\_smoo* smoothed distribution used to normalize the strip

Definition at line 185 of file *pearl-anglescan-process.ipf*.

**11.3.2.35** `variable pizza_service ( wave data, string nickname, variable theta_offset, variable tilt_offset, variable phi_offset, variable npolar = defaultValue, variable nograph = defaultValue, variable folding = defaultValue, variable xpdplot = defaultValue )`

create a pizza plot from a measured (energy-integrated) data strip

accepts angle-scan data as returned by [adh5\\_load\\_reduced\(\)](#), maps them onto a hemispherical scan grid, and displays a polar graph.

#### Parameters

<i>data</i>	2D intensity wave, see requirements above <ul style="list-style-type: none"> <li>• X-axis analyser angle</li> <li>• Y-axis manipulator scan. no specific ordering required. manipulator angle waves (ManipulatorTheta, ManipulatorTilt, ManipulatorPhi) must be in the subfolder <i>attr</i> below the data wave.</li> </ul>
-------------	--

## Parameters

<i>nickname</i>	nick name for output data <ul style="list-style-type: none"> <li>• in default mode, this will become the name of a child folder containing the output.</li> <li>• in XPDplot mode, this will become a prefix of the generated data in the root folder.</li> </ul>
<i>theta_offset</i>	manipulator theta angle corresponding to normal emission. the offset is subtracted from the ManipulatorTheta wave before processing.
<i>tilt_offset</i>	manipulator tilt angle corresponding to normal emission the offset is subtracted from the ManipulatorTilt wave before processing.
<i>phi_offset</i>	manipulator phi angle corresponding to phi_result = 0 the offset is subtracted from the ManipulatorPhi wave before processing.
<i>npolar</i>	number of polar angles, determines polar and azimuthal step size. default = 91 (1 degree steps)
<i>folding</i>	rotational averaging, default = 1
<i>nograph</i>	display a new graph window? <ul style="list-style-type: none"> <li>• 0 (default) display a new polar graph</li> <li>• 1 don't display a new graph</li> </ul>
<i>xpdplot</i>	XPDplot compatibility <ul style="list-style-type: none"> <li>• 0 (default) create waves in child folder \$nickname</li> <li>• 1 create waves in root folder (compatible with XPDplot)</li> </ul>

## Attention

if you modify the structure of the data wave, e.g. delete some angles, this function cannot be used because the manipulator settings do not correspond to the original manipulator waves! instead, create your own manipulator waves and use [pizza\\_service\\_2\(\)](#).

Definition at line 555 of file pearl-anglescan-process.ipf.

**11.3.2.36** `variable pizza_service_2 ( wave data, string nickname, wave m_theta, wave m_tilt, wave m_phi, variable npolar = defaultValue, variable nograph = defaultValue, variable folding = defaultValue, variable xpdplot = defaultValue )`

create a pizza plot from a measured (energy-integrated) data strip

accepts angle-scan data as returned by [adh5\\_load\\_reduced\(\)](#), maps them onto a hemispherical scan grid, and displays a polar graph.

the behaviour of this function is the same as [pizza\\_service\(\)](#) except that the manipulator waves are specified explicitly.

## Parameters

<i>data</i>	2D intensity wave, see requirements above <ul style="list-style-type: none"> <li>• X-axis analyser angle</li> <li>• Y-axis manipulator scan. no specific ordering required. manipulator angle waves (ManipulatorTheta, ManipulatorTilt, ManipulatorPhi) must be in the subfolder <i>attr</i> below the data wave.</li> </ul>
-------------	--

## Parameters

<i>nickname</i>	nick name for output data <ul style="list-style-type: none"> <li>• in default mode, this will become the name of a child folder containing the output.</li> <li>• in XPDplot mode, this will become a prefix of the generated data in the root folder.</li> </ul>
<i>m_theta</i>	manipulator theta angles, 0 = normal emission. size = <code>dimsize(data, 1)</code>
<i>m_tilt</i>	manipulator tilt angles, 0 = normal emission. size = <code>dimsize(data, 1)</code>
<i>m_phi</i>	manipulator phi angles, 0 = azimuthal origin. size = <code>dimsize(data, 1)</code>
<i>npolar</i>	number of polar angles, determines polar and azimuthal step size. default = 91 (1 degree steps)
<i>folding</i>	rotational averaging, default = 1
<i>nograph</i>	display a new graph window? <ul style="list-style-type: none"> <li>• 0 (default) display a new polar graph</li> <li>• 1 don't display a new graph</li> </ul>
<i>xpdplot</i>	XPDplot compatibility <ul style="list-style-type: none"> <li>• 0 (default) create waves in child folder \$nickname</li> <li>• 1 create waves in root folder (compatible with XPDplot)</li> </ul>

Definition at line 648 of file `pearl-anglescan-process.ipf`.

11.3.2.37 `static variable polar_graph_hook ( WMWinHookStruct * s ) [static]`

polar graph window hook

this hook converts the cursor positions to polar coordinates and displays them in a text box on the graph. the text box is visible while the cursor info box is visible.

Definition at line 2209 of file `pearl-anglescan-process.ipf`.

11.3.2.38 `variable quick_pizza_image ( wave data, string nickname, variable theta_offset, variable tilt_offset, variable phi_offset, variable npolar = defaultValue, variable nograph = defaultValue, variable folding = defaultValue )`

map angle scan data onto a rectangular grid in stereographic projection

accepts angle-scan data as returned by `adh5_load_reduced`, maps them onto a rectangular grid in stereographic projection

## Parameters

<i>data</i>	2D data wave, X-axis = analyser angle, Y-axis = manipulator scan (no specific ordering required)
-------------	--

## Precondition

manipulator angles as attributes in `attr` folder next to the data wave

## Warning

EXPERIMENTAL

Definition at line 2503 of file `pearl-anglescan-process.ipf`.

11.3.2.39 variable rotate\_hemi\_scan ( string *nickname*, variable *angle* )

azimuthally rotate a hemispherical scan dataset.

this function works only for hemi scans created by [make\\_hemi\\_grid\(\)](#) (or compatible functions).

## Parameters

<i>nickname</i>	name prefix for waves. source data must be in current data folder.
<i>angle</i>	azimuthal rotation angle in degrees.

Definition at line 1449 of file pearl-anglescan-process.ipf.

11.3.2.40 variable save\_hemi\_scan ( string *nickname*, string *pathname*, string *filename* )

save a hemispherical scan to an Igor text file

Definition at line 2614 of file pearl-anglescan-process.ipf.

11.3.2.41 variable set\_contrast ( variable *pcmin*, variable *pcmax*, string *graphname* = defaultValue, string *colortable* = defaultValue )

set the pseudocolor contrast by percentile.

set the minimum and maximum values of the pseudocolor scale such that a specified percentile of the distribution lies outside the limits.

the new contrast is applied to traces and images of the selected graph that have pseudocolor tables.

the function is not specific to angle scans. it can be used for any pseudocolor trace or image plots except contour plots.

## Parameters

<i>pcmin</i>	percentile below the minimum color (0-100).
<i>pcmax</i>	percentile above the maximum color (0-100).
<i>graphname</i>	name of graph. default: top graph.
<i>colortable</i>	name of new colortable. default: keep current table.

Definition at line 2961 of file pearl-anglescan-process.ipf.

11.3.2.42 variable set\_polar\_graph\_cursor ( string *nickname*, string *cursorname*, variable *polar\_angle*, variable *azim\_angle*, string *graphname* = defaultValue )

Definition at line 2229 of file pearl-anglescan-process.ipf.

11.3.2.43 variable show\_analyser\_line ( variable *theta*, variable *tilt*, variable *phi*, variable *theta\_offset*, variable *tilt\_offset*, variable *phi\_offset*, variable *npolar* = defaultValue, variable *nograph* = defaultValue, variable *xpdplot* = defaultValue )

calculate and display the line seen by the analyser for a specific emission angle

this can be used to compare to an hemispherical plot and check the manipulator angle.

## Parameters

## Parameters

<i>theta</i>	manipulator theta angle
<i>tilt</i>	manipulator tilt angle
<i>phi</i>	manipulator phi angle
<i>theta_offset</i>	manipulator theta angle corresponding to normal emission
<i>tilt_offset</i>	manipulator tilt angle corresponding to normal emission
<i>phi_offset</i>	manipulator phi angle corresponding to phi_result = 0
<i>npolar</i>	number of polar angles, determines polar and azimuthal step size. default = 91 (1 degree steps)
<i>nograph</i>	display a new graph window? <ul style="list-style-type: none"> <li>• 0 (default) display a new polar graph</li> <li>• 1 don't display a new graph</li> </ul>
<i>xpdplot</i>	XPDplot compatibility <ul style="list-style-type: none"> <li>• 0 (default) create waves in child folder \$nickname</li> <li>• 1 create waves in root folder (compatible with XPDplot)</li> </ul>

## Remarks

the function creates angle scan data under the nickname *analyser*.

Definition at line 765 of file pearl-anglescan-process.ipf.

#### 11.3.2.44 variable strip\_delete\_frames ( wave *strip*, variable *qlo*, variable *qhi*, wave *theta*, wave *tilt*, wave *phi* )

delete a contiguous range of frames from a strip.

this can be used to remove a region of bad frames due to, e.g., measurement problems. the function operates on 2D intensity data and manipulator coordinates at the same time.

## Parameters

in, out	<i>strip</i>	2D data, X-axis = analyser angle, Y-axis = arbitrary manipulator scan. the result is written to the original wave.
in, out	<i>theta</i>	1D data, manipulator scan. the result is written to the original wave.
in, out	<i>tilt</i>	1D data, manipulator scan. the result is written to the original wave.
in, out	<i>phi</i>	1D data, manipulator scan. the result is written to the original wave.
in	<i>qlo</i>	point index of first frame to delete.
in	<i>qhi</i>	point index of last frame to delete. qhi must be greater or equal than qlo.

Definition at line 105 of file pearl-anglescan-process.ipf.

#### 11.3.2.45 variable trim\_hemi\_scan ( string *nickname*, variable *theta\_max* )

trim a hemispherical scan at grazing angle

the function recalculates the values wave from totals and weights but sets elements above a given polar angle to nan.

## Parameters

<i>nickname</i>	name of the scan dataset. can be empty if no prefix is used. the dataset must be in the current datafolder.
<i>theta_max</i>	highest polar angle to keep (0...90 degrees).

Definition at line 2748 of file pearl-anglescan-process.ipf.

#### 11.3.2.46 static variable update\_polar\_info ( string *graphname* ) [static]

update the angles info based on cursors A and B of a given polar graph window

the function reads the projection mode from the user data of the graph window and the zeroAngleWhere variable from the associated WMPolarGraph data folder.

the calculated angles are written to the csrA\_theta, csrA\_phi, csrB\_theta, and csrB\_phi global variables in the polar graph data folder. the angles text box of the graph updates from to these variables dynamically.

## Parameters

<i>graphname</i>	name of polar graph window
------------------	----------------------------

Definition at line 2174 of file pearl-anglescan-process.ipf.

### 11.3.3 Variable Documentation

#### 11.3.3.1 const variable kProjArea = 2

Definition at line 2006 of file pearl-anglescan-process.ipf.

#### 11.3.3.2 const variable kProjDist = 0

Definition at line 2004 of file pearl-anglescan-process.ipf.

#### 11.3.3.3 const variable kProjGnom = 3

Definition at line 2007 of file pearl-anglescan-process.ipf.

#### 11.3.3.4 const variable kProjOrtho = 4

Definition at line 2008 of file pearl-anglescan-process.ipf.

#### 11.3.3.5 const variable kProjScaleArea = 2 [static]

Definition at line 2012 of file pearl-anglescan-process.ipf.

#### 11.3.3.6 const variable kProjScaleDist = 2 [static]

Definition at line 2010 of file pearl-anglescan-process.ipf.

11.3.3.7 `const variable kProjScaleGnom = 0.06744519021` `[static]`

Definition at line 2014 of file `pearl-anglescan-process.ipf`.

11.3.3.8 `const variable kProjScaleOrtho = 2` `[static]`

Definition at line 2015 of file `pearl-anglescan-process.ipf`.

11.3.3.9 `const variable kProjScaleStereo = 2` `[static]`

Definition at line 2011 of file `pearl-anglescan-process.ipf`.

11.3.3.10 `const variable kProjStereo = 1`

Definition at line 2005 of file `pearl-anglescan-process.ipf`.

## 11.4 `pearl-anglescan-tracker.ipf` File Reference

```
#include "pearl-area-profiles"
#include "pearl-area-import"
#include "pearl-scienta-preprocess"
#include "pearl-anglescan-process"
#include <New Polar Graphs>
```

### Functions

- static variable `AfterCompiledHook` ()  
*initialize package data once when the procedure is first loaded*
- static variable `init_package` ()
- static variable `save_prefs` ()  
*save persistent package data to the preferences file.*
- static variable `load_prefs` ()  
*load persistent package data from the preferences file.*
- static variable `IgorQuitHook` (string app)  
*disconnect EPICS channels before Igor quits.*
- variable `ast_setup` ()  
*set up data structures, display graph, and try to connect to analyser.*
- variable `ast_prepare` (variable theta\_offset=defaultValue, variable tilt\_offset=defaultValue, variable phi\_offset=defaultValue)  
*prepare for new measurement and clear the data buffer.*
- variable `ast_set_processing` (string reduction\_func, string reduction\_params)  
*set the data processing parameters*
- variable `ast_add_image` (wave image, variable theta, variable tilt, variable phi)  
*process and add a detector image to the tracker scan.*
- variable `ast_export` (dfref folder, string nickname, variable xpdplot=defaultValue)  
*export tracker data to a separate, independent data set.*
- variable `ast_import` (string nickname)  
*import tracker data from an existing angle scan dataset.*
- variable `ast_update_detector` (variable theta, variable tilt, variable phi, variable range)

- update the current position indicator.*
- variable `ast_close` ()
  - stop tracker, close graph, release data structures.*
- static variable `setup_data` ()
- static variable `extend_data` (variable `num_slices`)
  - extend the data buffer for the next polar scan*
- static variable `setup_detector` ()
- static variable `add_image_data` (wave image, variable `theta`, variable `tilt`, variable `phi`)
  - reduce a detector image and add the result to the data buffer.*
- static variable `process_image_data` ()
  - process the data buffer to generate the tracker dataset.*
- static variable `update_detector` (variable `theta`, variable `tilt`, variable `phi`, variable `range`)
  - update the current position indicator.*
- static variable `setup_graph` ()
  - create the graph window.*
- static variable `update_data_graph` ()
- static variable `update_detector_graph` ()
- static variable `epics_connect` ()
  - connect the angle scan tracker to EPICS*
- static variable `epics_disconnect_chid` (string `chid_var_name`)
- static variable `epics_disconnect` ()
- static variable `ast_window_hook` (WMWinHookStruct \*s)
  - window hook*
- variable `ast_callback_data` (variable `chan`)
  - callback function for new analyser data from EPICS.*
- variable `ast_callback_detector` (variable `chan`)
  - callback function for new detector state from EPICS.*
- variable `ast_callback_manip` (variable `chan`)
  - callback function for new manipulator position from EPICS.*
- static variable `bp_capture` (WMButtonAction \*ba)
- static variable `toggle_capture` ()
- static variable `update_capture` ()
- static variable `pmp_data` (WMPopupAction \*pa)
- static variable `pmp_data_mouseup` (WMPopupAction \*pa)
- static variable `export_tracker_data` ()
  - export tracker data (with prompt)*
- static variable `import_tracker_data` ()
  - import tracker data (with prompt)*
- static variable `save_tracker_data` ()
  - save tracker data to file (with prompt)*
- static variable `load_tracker_data` ()
  - import tracker data from file (with prompt)*
- static variable `pmp_parameters` (WMPopupAction \*pa)
- static variable `pmp_parameters_mouseup` (WMPopupAction \*pa)
- static variable `edit_reduction_params` ()
- static variable `edit_offsets` ()

## Variables

- [version](#)
- static const string [package\\_path](#) = "root:packages:pearl\_anglescan\_tracker:"  
*data folder path*
- static const string [prefs\\_objects](#) = "projection;theta\_offset;tilt\_offset;phi\_offset;reduction\_func;reduction\_↔  
params"  
*semicolon-separated list of persistent variable, string, and wave names*

### 11.4.1 Function Documentation

#### 11.4.1.1 static variable `add_image_data ( wave image, variable theta, variable tilt, variable phi )` [static]

reduce a detector image and add the result to the data buffer.

##### Parameters

<i>image</i>	detector image with correct X (energy) and Y (angle) scaling
<i>theta</i>	polar angle of manipulator
<i>tilt</i>	tilt angle of manipulator
<i>phi</i>	azimuthal angle of manipulator

the manipulator angles are corrected by the preset offsets internally.

Definition at line 453 of file `pearl-anglescan-tracker.ipf`.

#### 11.4.1.2 static variable `AfterCompiledHook ( )` [static]

initialize package data once when the procedure is first loaded

Definition at line 73 of file `pearl-anglescan-tracker.ipf`.

#### 11.4.1.3 variable `ast_add_image ( wave image, variable theta, variable tilt, variable phi )`

process and add a detector image to the tracker scan.

##### Parameters

<i>image</i>	detector image with correct X (energy) and Y (angle) scaling
<i>theta</i>	polar angle of manipulator
<i>tilt</i>	tilt angle of manipulator
<i>phi</i>	azimuthal angle of manipulator

the manipulator angles are corrected by the preset offsets internally.

Definition at line 285 of file `pearl-anglescan-tracker.ipf`.

#### 11.4.1.4 variable `ast_callback_data ( variable chan )`

callback function for new analyser data from EPICS.

Definition at line 846 of file `pearl-anglescan-tracker.ipf`.

**11.4.1.5 variable `ast_callback_detector` ( variable *chan* )**

callback function for new detector state from EPICS.

save the manipulator position at the beginning of image acquisition. it is used by [ast\\_callback\\_data\(\)](#).

Definition at line 951 of file `pearl-anglescan-tracker.ipf`.

**11.4.1.6 variable `ast_callback_manip` ( variable *chan* )**

callback function for new manipulator position from EPICS.

Definition at line 980 of file `pearl-anglescan-tracker.ipf`.

**11.4.1.7 variable `ast_close` ( )**

stop tracker, close graph, release data structures.

Definition at line 360 of file `pearl-anglescan-tracker.ipf`.

**11.4.1.8 variable `ast_export` ( *dfref folder*, *string nickname*, *variable xpdplot* = `defaultValue` )**

export tracker data to a separate, independent data set.

the exported data can then be used for further processing. the data is exported to the current data folder, or root if XPDplot compatibility is requested.

**Parameters**

<i>folder</i>	destination folder path
<i>nickname</i>	name prefix for waves
<i>xpdplot</i>	xpdplot compatibility, see <a href="#">make_hemi_grid()</a> for details <ul style="list-style-type: none"> <li>• 0 (default)</li> <li>• 1 create additional waves and notebook required by XPDplot</li> </ul>

Definition at line 307 of file `pearl-anglescan-tracker.ipf`.

**11.4.1.9 variable `ast_import` ( *string nickname* )**

import tracker data from an existing angle scan dataset.

**Parameters**

<i>nickname</i>	name prefix for waves. data must be in current data folder.
-----------------	---

Definition at line 329 of file `pearl-anglescan-tracker.ipf`.

**11.4.1.10 variable `ast_prepare` ( *variable theta\_offset* = `defaultValue`, *variable tilt\_offset* = `defaultValue`, *variable phi\_offset* = `defaultValue` )**

prepare for new measurement and clear the data buffer.

optionally, set new manipulator offsets. the offsets are the manipulator readback coordinates where the sample surface is oriented in normal emission, and the handle of the sample plate points horizontally to the left (9 o'clock).

## Parameters

<i>theta_offset</i>	set new theta offset. default: no change.
<i>tilt_offset</i>	set new tilt offset. default: no change.
<i>phi_offset</i>	set new phi offset. default: no change.

Definition at line 211 of file pearl-anglescan-tracker.ipf.

#### 11.4.1.11 variable `ast_set_processing ( string reduction_func, string reduction_params )`

set the data processing parameters

the parameters will be effective for subsequent measurements only. previously acquired data is not affected. the processing parameters are saved with the preferences.

## Parameters

<i>reduction_func</i>	name of custom reduction function, e.g. "int_linbg_reduction". any user-defined function with the same signature as <a href="#">adh5_default_reduction()</a> is allowed.
<i>reduction_params</i>	parameter string for the reduction function. the format depends on the actual function. for <code>int_linbg_reduction</code> , e.g., "Lcrop=0.1;Hcrop=0.1;Lsize=0.2;Hsize=0.2;Cpos=0.5;Csize=0.4".

Definition at line 261 of file pearl-anglescan-tracker.ipf.

#### 11.4.1.12 variable `ast_setup ( )`

set up data structures, display graph, and try to connect to analyser.

Definition at line 194 of file pearl-anglescan-tracker.ipf.

#### 11.4.1.13 variable `ast_update_detector ( variable theta, variable tilt, variable phi, variable range )`

update the current position indicator.

## Parameters

<i>theta</i>	polar angle of manipulator
<i>tilt</i>	tilt angle of manipulator
<i>phi</i>	azimuthal angle of manipulator
<i>range</i>	angle range (60 or 45)

the manipulator angles are corrected by the preset offsets internally.

Definition at line 349 of file pearl-anglescan-tracker.ipf.

#### 11.4.1.14 static variable `ast_window_hook ( WMWinHookStruct * s ) [static]`

window hook

disconnects from EPICS when the window is closed.

Definition at line 831 of file pearl-anglescan-tracker.ipf.

11.4.1.15 `static variable bp_capture ( WMBUTTONACTION * ba ) [static]`

Definition at line 1013 of file pearl-anglescan-tracker.ipf.

11.4.1.16 `static variable edit_offsets ( ) [static]`

Definition at line 1234 of file pearl-anglescan-tracker.ipf.

11.4.1.17 `static variable edit_reduction_params ( ) [static]`

Definition at line 1218 of file pearl-anglescan-tracker.ipf.

11.4.1.18 `static variable epics_connect ( ) [static]`

connect the angle scan tracker to EPICS

the tracker uses channels of the analyser and the manipulator.

if the EPICS XOP is not loaded, the function does nothing. if channels are not available, the function exits with an error code after a timeout of 5 seconds. the Igor run-time error status is reset to suppress the error dialog.

#### Returns

zero if successful, non-zero if an error occurred

**Todo** the X03DA channel names are hard-coded.

Definition at line 675 of file pearl-anglescan-tracker.ipf.

11.4.1.19 `static variable epics_disconnect ( ) [static]`

Definition at line 801 of file pearl-anglescan-tracker.ipf.

11.4.1.20 `static variable epics_disconnect_chid ( string chid_var_name ) [static]`

Definition at line 787 of file pearl-anglescan-tracker.ipf.

11.4.1.21 `static variable export_tracker_data ( ) [static]`

export tracker data (with prompt)

Definition at line 1095 of file pearl-anglescan-tracker.ipf.

11.4.1.22 `static variable extend_data ( variable num_slices ) [static]`

extend the data buffer for the next polar scan

call this function if the buffer is full.

#### Parameters

<i>num_slices</i>	number of slices that the measurement contains
-------------------	--

Definition at line 402 of file pearl-anglescan-tracker.ipf.

**11.4.1.23 static variable IgorQuitHook ( string *app* ) [static]**

disconnect EPICS channels before Igor quits.

Definition at line 188 of file pearl-anglescan-tracker.ipf.

**11.4.1.24 static variable import\_tracker\_data ( ) [static]**

import tracker data (with prompt)

Definition at line 1125 of file pearl-anglescan-tracker.ipf.

**11.4.1.25 static variable init\_package ( ) [static]**

Definition at line 98 of file pearl-anglescan-tracker.ipf.

**11.4.1.26 static variable load\_prefs ( ) [static]**

load persistent package data from the preferences file.

the preferences file is an Igor packed experiment file in a special preferences folder.

this function is called automatically when the procedure is first compiled, or whenever the user clicks the corresponding button.

Definition at line 163 of file pearl-anglescan-tracker.ipf.

**11.4.1.27 static variable load\_tracker\_data ( ) [static]**

import tracker data from file (with prompt)

Definition at line 1165 of file pearl-anglescan-tracker.ipf.

**11.4.1.28 static variable pmp\_data ( WMPopupAction \* *pa* ) [static]**

Definition at line 1061 of file pearl-anglescan-tracker.ipf.

**11.4.1.29 static variable pmp\_data\_mouseup ( WMPopupAction \* *pa* ) [static]**

Definition at line 1075 of file pearl-anglescan-tracker.ipf.

**11.4.1.30 static variable pmp\_parameters ( WMPopupAction \* *pa* ) [static]**

Definition at line 1184 of file pearl-anglescan-tracker.ipf.

**11.4.1.31 static variable pmp\_parameters\_mouseup ( WMPopupAction \* *pa* ) [static]**

Definition at line 1198 of file pearl-anglescan-tracker.ipf.

**11.4.1.32 static variable process\_image\_data ( ) [static]**

process the data buffer to generate the tracker dataset.

Definition at line 502 of file pearl-anglescan-tracker.ipf.

**11.4.1.33 static variable save\_prefs ( ) [static]**

save persistent package data to the preferences file.

this function is called when the user clicks the corresponding button.

Definition at line 139 of file pearl-anglescan-tracker.ipf.

**11.4.1.34 static variable save\_tracker\_data ( ) [static]**

save tracker data to file (with prompt)

Definition at line 1154 of file pearl-anglescan-tracker.ipf.

**11.4.1.35 static variable setup\_data ( ) [static]**

Definition at line 372 of file pearl-anglescan-tracker.ipf.

**11.4.1.36 static variable setup\_detector ( ) [static]**

Definition at line 434 of file pearl-anglescan-tracker.ipf.

**11.4.1.37 static variable setup\_graph ( ) [static]**

create the graph window.

Definition at line 589 of file pearl-anglescan-tracker.ipf.

**11.4.1.38 static variable toggle\_capture ( ) [static]**

Definition at line 1027 of file pearl-anglescan-tracker.ipf.

**11.4.1.39 static variable update\_capture ( ) [static]**

Definition at line 1045 of file pearl-anglescan-tracker.ipf.

**11.4.1.40 static variable update\_data\_graph ( ) [static]**

Definition at line 640 of file pearl-anglescan-tracker.ipf.

**11.4.1.41 static variable update\_detector ( variable *theta*, variable *tilt*, variable *phi*, variable *range* ) [static]**

update the current position indicator.

**Parameters**

<i>theta</i>	polar angle of manipulator
<i>tilt</i>	tilt angle of manipulator
<i>phi</i>	azimuthal angle of manipulator
<i>range</i>	angle range (60 or 45)

the manipulator angles are corrected by the preset offsets internally.

Definition at line 554 of file pearl-anglescan-tracker.ipf.

#### 11.4.1.42 static variable `update_detector_graph` ( ) [static]

Definition at line 652 of file pearl-anglescan-tracker.ipf.

### 11.4.2 Variable Documentation

#### 11.4.2.1 const string `package_path` = "root:packages:pearl\_anglescan\_tracker:" [static]

data folder path

Definition at line 68 of file pearl-anglescan-tracker.ipf.

#### 11.4.2.2 const string `prefs_objects` = "projection;theta\_offset;tilt\_offset;phi\_offset;reduction\_func;reduction\_params" [static]

semicolon-separated list of persistent variable, string, and wave names

Definition at line 70 of file pearl-anglescan-tracker.ipf.

#### 11.4.2.3 version

**Initial value:**

= 1.6

```
static const string package_name = "pearl_anglescan_tracker"
```

Definition at line 5 of file pearl-anglescan-tracker.ipf.

## 11.5 pearl-area-display.ipf File Reference

visualization tools for 2D and 3D data.

### Namespaces

- [PearlAreaDisplay](#)

*instant visualization of angle scan and manipulator position.*

## Functions

- static string [graphname\\_from\\_dfref](#) (dfref df, string prefix)  
*compose a valid and unique graph name from a data folder reference*
- string [ad\\_display](#) (wave image)  
*open a new graph window with a 2D image.*
- string [ad\\_display\\_histogram](#) (wave image)  
*display the histogram of a 2D image.*
- string [ad\\_display\\_profiles](#) (wave image, string filter=defaultValue)  
*open a new profiles graph window.*
- variable [ad\\_update\\_profiles](#) (wave image)  
*update a profiles graph with new data.*
- variable [ad\\_profiles\\_cursor\\_mode](#) (wave image, variable mode)  
*switch cursors on a profiles graph*
- variable [ad\\_profiles\\_set\\_cursor](#) (wave image, string cursorname, variable xa, variable ya, variable pscale=defaultValue)  
*move a cursor to the specified position in a profiles graph.*
- variable [ad\\_profiles\\_crosshairs](#) (wave image, variable clear=defaultValue)  
*draw permanent crosshairs in a profiles graph.*
- static wave [get\\_source\\_image](#) (wave view)  
*find the source image wave corresponding to the given view.*
- static dfr [make\\_view\\_folder](#) (wave source)  
*create a view data folder.*
- static dfr [get\\_view\\_folder](#) (wave source)  
*find the view data folder corresponding to the given source.*
- static wave [get\\_view\\_image](#) (wave source)  
*find the view image wave corresponding to the given source.*
- static variable [bp\\_reset\\_cursors](#) (WMButtonAction \*ba)
- static variable [svp\\_smoothing](#) (WMSetVariableAction \*sva)
- static variable [pmp\\_export](#) (WMPopupAction \*pa)
- variable [ad\\_profiles\\_hook](#) (WMWinHookStruct \*s)  
*hook function for user events in the profiles window.*
- variable [ad\\_calc\\_cursor\\_profiles](#) (wave image)  
*calculate profiles, statistics, and histogram of a cross-hair delimited region of interest.*
- variable [ad\\_calc\\_profiles](#) (wave image, variable pa, variable qa, variable pb, variable qb)  
*calculate profiles, statistics, and histogram of a rectangular region of interest.*
- variable [ad\\_export\\_profile](#) (wave view\_image, variable dim, variable trace=defaultValue, variable show=defaultValue, variable overwrite=defaultValue)  
*export a profile from a profiles graph to the source data folder.*
- static variable [set\\_trace\\_colors](#) (string graphname)
- variable [ad\\_calc\\_histogram](#) (wave image)  
*calculate the histogram.*
- variable [ad\\_default\\_image\\_filter](#) (wave image, string options)  
*abstract filter function for image display.*
- variable [ad\\_box\\_filter](#) (wave image, string options)  
*boxcar smoothing filter.*
- variable [ad\\_transpose\\_filter](#) (wave image, string options)  
*transpose image filter.*
- string [ad\\_display\\_brick](#) (wave data)  
*open a new "gizmo" window with three-dimensional data.*
- variable [ad\\_brick\\_slicer](#) (wave data)

- open a slicer panel for 3D data.*
- string `ad_display_slice` (wave data)
  - display three-dimensional data by 2D slice.*
- static variable `update_slice_info` ()
  - update controls with data scale limits.*
- variable `ad_gizmo_set_plane` (wave brick, variable dim, variable value)
  - set the position of a slicing plane of a 3D brick in a Gizmo window.*
- variable `ad_profiles_set_slice` (wave brick, variable dim, variable value)
  - set the position of the slicing plane of a 3D brick in a profiles window.*
- static variable `slp_slice_position` (WMSliderAction \*sa)
  - set slice coordinate (slider procedure).*
- static variable `svp_slice_position` (WMSetVariableAction \*sva)
  - set slice coordinate (button procedure).*
- static variable `bp_move_slice` (WMButtonAction \*ba)
  - move slice (button procedure).*
- static variable `bp_extract_slice` (WMButtonAction \*ba)
  - export a slice (button procedure).*
- static variable `bp_move_slice_center` (wave brick, variable dim, string posvariable)
  - move the slice to the center of the dimension (button procedure).*
- static variable `ad_slicer_move_bg` (WMBackgroundStruct \*s)
  - move a slice by one step (background task).*
- variable `ad_slicer_init_bg` ()
  - initialize the slice animation background task.*
- variable `ad_slicer_start_bg` (wave brick, variable dimension, string posvariable, variable delta)
  - start the animation.*
- variable `ad_slicer_stop_bg` (string posvariable)
  - stop the animation.*

### 11.5.1 Detailed Description

visualization tools for 2D and 3D data.

these tools were initially developed for monitoring output from EPICS area detector software. they are, however, useful for any kind of intensity versus x,y(z) data.

### 11.5.2 2D data

TO DO...

### 11.5.3 3D data

### 11.5.4 Function Documentation

#### 11.5.4.1 variable `ad_box_filter` ( wave image, string options )

boxcar smoothing filter.

filters the image in X and Y directions using Igor's Smooth operation.

#### Parameters

<i>image</i>	image to be filtered: original data and filter result.
--------------	--

## Parameters

<i>options</i>	smoothing factors in <code>key1=value1;key2=value2; . . . format.</code> <ul style="list-style-type: none"> <li>• SmoothingX</li> <li>• SmoothingY</li> </ul>
----------------	---

Definition at line 1044 of file pearl-area-display.ipf.

11.5.4.2 variable `ad_brick_slicer ( wave data )`

open a slicer panel for 3D data.

if a panel exists, bring it to the front.

## Parameters

<i>data</i>	three-dimensional wave.
-------------	-------------------------

Definition at line 1197 of file pearl-area-display.ipf.

11.5.4.3 variable `ad_calc_cursor_profiles ( wave image )`

calculate profiles, statistics, and histogram of a cross-hair delimited region of interest.

## Parameters

<i>image</i>	wave which contains the image data from the detector.
--------------	---

the function expects further objects as created by [ad\\_display\\_profiles\(\)](#) in the same data folder as the image wave. the most recent profiles graph of the image must exist, and the cursors A and B must be set on the image.

Definition at line 726 of file pearl-area-display.ipf.

11.5.4.4 variable `ad_calc_histogram ( wave image )`

calculate the histogram.

## Parameters

<i>image</i>	wave which contains the image data from the detector. the function expects further objects as created by <a href="#">ad_display_histogram()</a> in the same data folder as the image wave.
--------------	--

Definition at line 1007 of file pearl-area-display.ipf.

11.5.4.5 variable `ad_calc_profiles ( wave image, variable pa, variable qa, variable pb, variable qb )`

calculate profiles, statistics, and histogram of a rectangular region of interest.

the region of interest a rectangle spanned by the two points A and B. pixels at these coordinates are included.

**Parameters**

<i>image</i>	wave which contains the image data.
<i>pa</i>	first point coordinate of A.
<i>qa</i>	second point coordinate of A.
<i>pb</i>	first point coordinate of B.
<i>qb</i>	second point coordinate of B.

the function expects further objects as created by [ad\\_display\\_profiles\(\)](#) in the same data folder as the image wave.

this function does not require that the graph exists as long as the data folder is complete.

Definition at line 773 of file pearl-area-display.ipf.

#### 11.5.4.6 variable `ad_default_image_filter` ( wave *image*, string *options* )

abstract filter function for image display.

this is a function prototype for filtering two-dimensional data for preview. to write your own filter, define a new function which has the same signature.

**Parameters**

<i>image</i>	image to be filtered: original data and filter result.
<i>options</i>	filter options in <code>key1=value1;key2=value2; . . .</code> format.

**Returns**

the result must be written to the incoming image wave.

Definition at line 1030 of file pearl-area-display.ipf.

#### 11.5.4.7 string `ad_display` ( wave *image* )

open a new graph window with a 2D image.

this is essentially `display; appendimage`. the graph is directly linked to the image wave. it is, thus, updated automatically.

**Parameters**

<i>image</i>	wave which contains the image data.
--------------	-------------------------------------

**Returns**

(string) name of the graph window

Definition at line 84 of file pearl-area-display.ipf.

#### 11.5.4.8 string `ad_display_brick` ( wave *data* )

open a new "gizmo" window with three-dimensional data.

**Parameters**

<i>data</i>	three-dimensional wave.
-------------	-------------------------

**Returns**

name of the gizmo window.

Definition at line 1082 of file pearl-area-display.ipf.

**11.5.4.9 string ad\_display\_histogram ( wave *image* )**

display the histogram of a 2D image.

the function will create additional objects in the same data folder as the image. this objects are displayed in the graph and are updated by calling [ad\\_calc\\_profiles\(\)](#). see the code.

**Parameters**

<i>image</i>	wave which contains the image data from the detector.
--------------	---

**Returns**

(string) name of the graph window

Definition at line 114 of file pearl-area-display.ipf.

**11.5.4.10 string ad\_display\_profiles ( wave *image*, string *filter* = defaultValue )**

open a new profiles graph window.

opens an extended graph window with profiles for the specified image. the function copies/creates all necessary data structures in a subfolder of the one which contains the image wave. the data folder name is derived from the image wave name by prefixing with "view\_". there can be at most one profiles window of each image wave. the original wave must not be renamed while the graph window is used. to update the graph after modifying the original wave, call [ad\\_update\\_profiles\(\)](#).

**Parameters**

<i>image</i>	wave which contains the image data.
<i>filter</i>	name of a filter function which maps the original data to the displayed data. the function must have the same parameters as <a href="#">ad_default_image_filter()</a> . default: boxcar average ( <a href="#">ad_box_filter()</a> ) using parameters view_filter_smoothing_x and _y.

**Returns**

name of the graph window

Definition at line 165 of file pearl-area-display.ipf.

**11.5.4.11 string ad\_display\_slice ( wave *data* )**

display three-dimensional data by 2D slice.

to select the slice data to display, call [ad\\_profiles\\_set\\_slice\(\)](#), or open a [ad\\_brick\\_slicer\(\)](#) panel. do not modify the content of the created view\_data folder.

#### Parameters

<i>data</i>	three-dimensional wave.
-------------	-------------------------

#### Returns

name of the graph window.

Definition at line 1345 of file pearl-area-display.ipf.

**11.5.4.12** variable `ad_export_profile ( wave view_image, variable dim, variable trace = defaultValue, variable show = defaultValue, variable overwrite = defaultValue )`

export a profile from a profiles graph to the source data folder.

this function does not require that the show exists as long as the view data folder is complete.

#### Parameters

<i>view_image</i>	wave which contains the view image (image wave on display in profiles window). the function expects further objects as created by <a href="#">ad_display_profiles()</a> in the same data folder as the view_image wave.
<i>dim</i>	dimension index (0 = x, 1 = y).
<i>trace</i>	select profile trace: <ul style="list-style-type: none"> <li>• 0 = cursor A</li> <li>• 1 = cursor B</li> <li>• 2 = average between cursors (default)</li> </ul>
<i>show</i>	display mode: <ul style="list-style-type: none"> <li>• 0 = do not show (default)</li> <li>• 1 = display in new graph, or append to existing graph</li> <li>• 2 = collate: common graph for all profiles of a dimension. rename graph manually to detach it from future additions.</li> </ul>
<i>overwrite</i>	overwrite mode: <ul style="list-style-type: none"> <li>• 0 = create new wave (default). wave name may get a suffix to be unique.</li> <li>• 1 = overwrite existing wave</li> </ul>

Definition at line 875 of file pearl-area-display.ipf.

**11.5.4.13** variable `ad_gizmo_set_plane ( wave brick, variable dim, variable value )`

set the position of a slicing plane of a 3D brick in a Gizmo window.

#### Parameters

<i>brick</i>	original data wave.
--------------	---------------------

## Parameters

<i>dim</i>	dimension index: 0, 1, or 2.
<i>value</i>	new coordinate of the slicing plane (axis scaling).

## Returns

0 if successful, non-zero otherwise

Definition at line 1422 of file pearl-area-display.ipf.

11.5.4.14 variable `ad_profiles_crosshairs` ( wave *image*, variable *clear* = defaultValue )

draw permanent crosshairs in a profiles graph.

adds dash-dotted horizontal and vertical crosshairs to a profiles graph. for each active cursor A and/or B, a pair of lines crossing at the cursor position is added. existing crosshairs are moved to the current cursor positions. optionally, existing crosshairs are removed from the graph.

in contrast to the cursors, these crosshairs will be exported and printed with the graph. they are drawn using Igor's DrawLine operation. all lines drawn by this function are part of the "crosshairs" draw group.

## Parameters

<i>image</i>	image displayed in the graph. this is the original image, not the one in the view data folder.
<i>clear</i>	0 (default) = add/update lines. 1 = remove lines.

Definition at line 481 of file pearl-area-display.ipf.

11.5.4.15 variable `ad_profiles_cursor_mode` ( wave *image*, variable *mode* )

switch cursors on a profiles graph

the standard cursors allow to select the profiles to display in the profiles panes. additional cursors are shown in the profiles panes.

in the background selection mode, additional cursors allow the user to select the limits of the background and peak integration regions. the meaning of the cursors depends on the particular processing function.

## Parameters

<i>mode</i>	cursor mode. <ul style="list-style-type: none"> <li>• 0 (default) standard profile selection. cursors C-F on profile panes.</li> <li>• 1 background selection. cursors A-F on image.</li> </ul>
<i>image</i>	image displayed in the graph. this is the original image, not the one in the view data folder.

Definition at line 368 of file pearl-area-display.ipf.

11.5.4.16 variable `ad_profiles_hook` ( WMWinHookStruct \* s )

hook function for user events in the profiles window.

Definition at line 686 of file pearl-area-display.ipf.

**11.5.4.17** `variable ad_profiles_set_cursor ( wave image, string cursorname, variable xa, variable ya, variable pscale = defaultvalue )`

move a cursor to the specified position in a profiles graph.

this function can only set cursors in the image part of the profiles graph.

#### Parameters

<i>image</i>	image displayed in the graph. this is the original image, not the one in the view data folder.
<i>cursorname</i>	name of the cursor, e.g. "A" or "B". other cursors are allowed but need to be activated separately.
<i>xa</i>	x-coordinate to move the cursor to. the position is coerced to the image scale. +/-inf is allowed.
<i>ya</i>	y-coordinate to move the cursor to. the position is coerced to the image scale. +/-inf is allowed.
<i>pscale</i>	scaling of the position argument <ul style="list-style-type: none"> <li>• 0 (default) wave scaling</li> <li>• 1 point scaling</li> </ul>

Definition at line 429 of file pearl-area-display.ipf.

**11.5.4.18** `variable ad_profiles_set_slice ( wave brick, variable dim, variable value )`

set the position of the slicing plane of a 3D brick in a profiles window.

#### Parameters

<i>brick</i>	original data wave.
<i>dim</i>	dimension index: 0, 1, or 2.
<i>value</i>	new coordinate of the slicing plane (axis scaling).

#### Returns

0 if successful, non-zero otherwise

Definition at line 1458 of file pearl-area-display.ipf.

**11.5.4.19** `variable ad_slicer_init_bg ( )`

initialize the slice animation background task.

Definition at line 1717 of file pearl-area-display.ipf.

**11.5.4.20** `static variable ad_slicer_move_bg ( WMBackgroundStruct * s )` [static]

move a slice by one step (background task).

Definition at line 1678 of file pearl-area-display.ipf.

**11.5.4.21** `variable ad_slicer_start_bg ( wave brick, variable dimension, string posvariable, variable delta )`

start the animation.

## Parameters

<i>brick</i>	3D data wave
<i>dimension</i>	dimension to animate, 0, 1, or 2.
<i>posvariable</i>	full path to the global position variable.
<i>delta</i>	step increment, should be +/- dimdelta.

Definition at line 1741 of file pearl-area-display.ipf.

11.5.4.22 variable `ad_slicer_stop_bg` ( *string posvariable* )

stop the animation.

## Parameters

<i>posvariable</i>	full path to the global position variable.
--------------------	--

Definition at line 1783 of file pearl-area-display.ipf.

11.5.4.23 variable `ad_transpose_filter` ( *wave image*, *string options* )

transpose image filter.

transposes the image.

## Parameters

<i>image</i>	image to be transposed: original data and result.
<i>options</i>	not used. should be empty.

Definition at line 1066 of file pearl-area-display.ipf.

11.5.4.24 variable `ad_update_profiles` ( *wave image* )

update a profiles graph with new data.

## Parameters

<i>image</i>	wave which contains the image data. must be the same (by data folder and name) wave used with <a href="#">ad_display_profiles()</a> .
--------------	---

Definition at line 321 of file pearl-area-display.ipf.

11.5.4.25 static variable `bp_extract_slice` ( *WMBUTTONACTION \* ba* ) [static]

export a slice (button procedure).

extract a slice and saves it in a separate wave.

Definition at line 1612 of file pearl-area-display.ipf.

**11.5.4.26** static variable `bp_move_slice ( WMButtonAction * ba )` `[static]`

move slice (button procedure).

Definition at line 1563 of file `pearl-area-display.ipf`.

**11.5.4.27** static variable `bp_move_slice_center ( wave brick, variable dim, string posvariable )` `[static]`

move the slice to the center of the dimension (button procedure).

Definition at line 1666 of file `pearl-area-display.ipf`.

**11.5.4.28** static variable `bp_reset_cursors ( WMButtonAction * ba )` `[static]`

Definition at line 618 of file `pearl-area-display.ipf`.

**11.5.4.29** static wave `get_source_image ( wave view )` `[static]`

find the source image wave corresponding to the given view.

#### Returns

wave reference of the original data wave. the reference may be invalid if the source wave cannot be found.

Definition at line 547 of file `pearl-area-display.ipf`.

**11.5.4.30** static dfr `get_view_folder ( wave source )` `[static]`

find the view data folder corresponding to the given source.

the result data folder reference may be invalid if no view is currently open. use the built-in `DataFolderRefStatus` function to check for validity.

#### Parameters

<i>source</i>	wave which contains the image data. must be the same (by data folder and name) wave used with <a href="#">ad_display_profiles()</a> .
---------------	---

Definition at line 585 of file `pearl-area-display.ipf`.

**11.5.4.31** static wave `get_view_image ( wave source )` `[static]`

find the view image wave corresponding to the given source.

#### Parameters

<i>source</i>	wave which contains the image data. must be the same (by data folder and name) wave used with <a href="#">ad_display_profiles()</a> .
---------------	---

Definition at line 608 of file `pearl-area-display.ipf`.

**11.5.4.32** static string `graphname_from_dhref ( dhref df, string prefix )` `[static]`

compose a valid and unique graph name from a data folder reference

Definition at line 56 of file pearl-area-display.ipf.

**11.5.4.33** `static dfr make_view_folder ( wave source ) [static]`

create a view data folder.

Definition at line 561 of file pearl-area-display.ipf.

**11.5.4.34** `static variable pmp_export ( WMPopupAction * pa ) [static]`

Definition at line 662 of file pearl-area-display.ipf.

**11.5.4.35** `static variable set_trace_colors ( string graphname ) [static]`

Definition at line 986 of file pearl-area-display.ipf.

**11.5.4.36** `static variable slp_slice_position ( WMSliderAction * sa ) [static]`

set slice coordinate (slider procedure).

Definition at line 1499 of file pearl-area-display.ipf.

**11.5.4.37** `static variable svp_slice_position ( WMSetVariableAction * sva ) [static]`

set slice coordinate (button procedure).

Definition at line 1531 of file pearl-area-display.ipf.

**11.5.4.38** `static variable svp_smoothing ( WMSetVariableAction * sva ) [static]`

Definition at line 637 of file pearl-area-display.ipf.

**11.5.4.39** `static variable update_slice_info ( ) [static]`

update controls with data scale limits.

current folder must be slicer info

Definition at line 1386 of file pearl-area-display.ipf.

## 11.6 pearl-area-import.ipf File Reference

HDF5 file import from EPICS area detectors.

```
#include <HDF5 Browser>
#include "pearl-gui-tools"
```

### Namespaces

- [PearlAreaImport](#)

*HDF5 file import from EPICS area detectors.*

## Functions

- static variable [BeforeFileOpenHook](#) (variable refNum, string fileName, string path, string type, string creator, variable kind)  
*callback function for drag&drop of HDF5 files into Igor.*
- string [ad\\_suggest\\_foldername](#) (string filename, variable ignoredate=defaultValue, string sourcename=defaultValue, variable unique=defaultValue)  
*generate the name of a data folder based on a file name.*
- variable [ad\\_load\\_dialog](#) (string APathName)  
*load area detector data files selected in a file dialog window*
- string [adh5\\_load\\_complete](#) (string ANickName, string APathName, string AFileName, variable load\_data=defaultValue, variable load\_attr=defaultValue)  
*import everything from a HDF5 file created by the Area Detector software.*
- string [adh5\\_load\\_reduced](#) (string ANickName, string APathName, string AFileName, funcref reduction\_func, string reduction\_param, variable load\_data=defaultValue, variable load\_attr=defaultValue, variable progress=defaultValue)  
*load and reduce a dataset from a HDF5 file created by the Area Detector software.*
- string [adh5\\_load\\_preview](#) (string ANickName, string APathName, string AFileName, variable load\_data=defaultValue, variable load\_attr=defaultValue)  
*load a single image from a HDF5 file created by the Area Detector software.*
- string [adh5\\_load\\_info](#) (string APathName, string AFileName)  
*load descriptive info from a HDF5 file created by the Area Detector software.*
- variable [adh5\\_load\\_detector](#) (variable fileID, string detectorpath)  
*load the detector dataset from the open HDF5 file.*
- variable [adh5\\_redim](#) (wave data)  
*redimension a multi-dimensional area detector array loaded from HDF5.*
- static dfr [GetAttrDataFolderDFR](#) (wave data)  
*find the attributes data folder of an area detector dataset.*
- variable [adh5\\_scale](#) (wave data, string source=defaultValue)  
*set the dimension scales of an area detector dataset.*
- variable [adh5\\_load\\_detector\\_slabs](#) (variable fileID, string detectorpath, variable progress=defaultValue)  
*load the detector dataset from the open HDF5 file.*
- variable [adh5\\_load\\_detector\\_image](#) (variable fileID, string detectorpath, variable dim2start, variable dim2count, variable dim3start, variable dim3count)  
*load a single image from the detector dataset of the open HDF5 file*
- string [adh5\\_list\\_reduction\\_funcs](#) ()  
*get a list of functions which can be used as reduction functions.*
- threadsafe variable [adh5\\_default\\_reduction](#) (wave source, wave dest1, wave dest2, string \*param)  
*function prototype for adh5\_load\_reduced\_detector*
- threadsafe variable [adh5\\_setup\\_profile](#) (wave image, wave profile, variable dim)  
*set up a one-dimensional wave for a line profile based on a 2D original wave.*
- string [adh5\\_test\\_reduction\\_func](#) (wave source, wave dest1, wave dest2, funcref reduction\_func, string reduction\_param)  
*wrapper function for testing reduction functions from the command line.*
- variable [adh5\\_load\\_reduced\\_detector](#) (variable fileID, string detectorpath, funcref reduction\_func, string reduction\_param, variable progress=defaultValue, variable nthreads=defaultValue)  
*load a reduced detector dataset from the open HDF5 file.*
- static threadsafe variable [reduce\\_slab\\_worker](#) (funcref reduction\_func)
- static threadsafe variable [reduce\\_slab\\_image](#) (wave slabdata, wave image, wave profile1, wave profile2, funcref reduction\_func, string reduction\_param)
- variable [adh5\\_loadattr\\_all](#) (variable fileID, string attributespath)  
*load an NDAttributes group from an open HDF5 file into the current data folder.*

- static variable `read_attribute_info` (string datawavename, string source, variable \*idest)  
*sub-function of `adh5_loadattr_all`.*
- variable `adh5_scale_scienta` (wave data)  
*set the energy and angle scales of an area detector dataset from the Scienta analyser.*
- variable `adh5_scale_scan` (wave data)  
*scales the extra dimensions of an area detector dataset according to the EPICS scan*

### 11.6.1 Detailed Description

HDF5 file import from EPICS area detectors.

HDF5 file import from EPICS area detectors such as CCD cameras, 2D electron analysers

as of Igor 6.3, Igor can open datasets of up to rank 4. i.e. the extra dimension Y of the file plugin cannot be used. the extra dimensions N and X are supported.

### 11.6.2 Function Documentation

#### 11.6.2.1 variable `ad_load_dialog` ( string *APathName* )

load area detector data files selected in a file dialog window

##### Parameters

<i>APathName</i>	Igor symbolic path name. if empty, Igor will choose a folder on its own
------------------	---

Definition at line 159 of file `pearl-area-import.ipf`.

#### 11.6.2.2 string `ad_suggest_foldername` ( string *filename*, variable *ignoreddate* = defaultValue, string *sourcename* = defaultValue, variable *unique* = defaultValue )

generate the name of a data folder based on a file name.

if the file name follows the naming convention source-date-index.extension, the function tries to generate the nick name as `source_date_index`. otherwise it's just a cleaned up version of the file name.

date must be in `yyyymmdd` or `yymmdd` format and is clipped to the short `yymmdd` format. index should be a running numeric index of up to 6 digits, or the time encoded as `hhmmss`. however, in the current version index can be any string that can be a valid Igor folder name.

##### Parameters

<i>filename</i>	file name, including extension. can also include a folder path (which is ignored). the extension is currently ignored, but may be used to select the parent folder in a later version.
<i>ignoreddate</i>	if non-zero, the nick name will not include the date part. defaults to zero.
<i>sourcename</i>	nick name of the data source. by default, the function tries to detect the source from the file name. this option can be used to override auto-detection. the automatic source names are: <code>sci</code> (scienta by area detector), <code>psh</code> (pshell), <code>sl</code> (optics slit camera by area detector), <code>es</code> (end station camera by area detector), <code>xy</code> (unidentified).
<i>unique</i>	if non-zero, the resulting name is made a unique data folder name in the current data folder defaults to zero.

Definition at line 98 of file `pearl-area-import.ipf`.

### 11.6.2.3 threadsafe variable `adh5_default_reduction ( wave source, wave dest1, wave dest2, string * param )`

function prototype for `adh5_load_reduced_detector`

derived functions reduce a two-dimensional dataset to a one-dimensional dataset, e.g. by ROI-integration, curve fitting, etc. each destination wave is a one-dimensional intensity distribution. the function must redimension each of these waves to one of the image dimensions by calling the `adh5_setup_profile()` function. this function will also copy the scale information and dimension labels, which is important for the proper scaling of the result.

the meaning of the data in `dest1` and `dest2` is up to the particular function, e.g. `dest1` could hold the mean value and `dest2` the one-sigma error, or `dest1` could hold the X-profile, and `dest2` the Y-profile.

#### Parameters

<i>source</i>	source wave two-dimensional intensity distribution (image)
<i>dest1, dest2</i>	destination waves
<i>param</i>	string with optional parameters, shared between calls. this is a pass-by-reference argument, the function may modify the string

#### Returns

zero if successful, non-zero if an error occurs.

Definition at line 1093 of file `pearl-area-import.ipf`.

### 11.6.2.4 string `adh5_list_reduction_funcs ( )`

get a list of functions which can be used as reduction functions.

the function evaluates only the function arguments, it may thus include functions which are not suitable as reduction functions.

Definition at line 1033 of file `pearl-area-import.ipf`.

### 11.6.2.5 string `adh5_load_complete ( string ANickName, string APathName, string AFileName, variable load_data = defaultValue, variable load_attr = defaultValue )`

import everything from a HDF5 file created by the Area Detector software.

if the data is from the electron analyser driver and some special attributes are included, the function will set the scales of the image dimensions.

#### Parameters

<i>ANickName</i>	destination folder name (top level under root)
<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in <i>FileName</i> or a dialog box should be displayed
<i>AFileName</i>	if empty a dialog box shows up
<i>load_data</i>	1 (default): load data; 0: do not load data
<i>load_attr</i>	1 (default): load attributes; 0: do not load attributes for proper wave scaling, the attributes must be loaded

Definition at line 199 of file `pearl-area-import.ipf`.

#### 11.6.2.6 variable `adh5_load_detector` ( variable `fileID`, string `detectorpath` )

load the detector dataset from the open HDF5 file.

the function loads the whole dataset at once and redimensions it so that the image dimensions are X and Y

##### Parameters

<code>fileID</code>	ID of open HDF5 file from <code>HDF5OpenFile</code>
<code>detectorpath</code>	path to detector group in the HDF5 file

Definition at line 573 of file `pearl-area-import.ipf`.

#### 11.6.2.7 variable `adh5_load_detector_image` ( variable `fileID`, string `detectorpath`, variable `dim2start`, variable `dim2count`, variable `dim3start`, variable `dim3count` )

load a single image from the detector dataset of the open HDF5 file

the function can average over a region in the extra dimensions.

##### Parameters

<code>fileID</code>	ID of open HDF5 file from <code>HDF5OpenFile</code>
<code>detectorpath</code>	path to detector group in the HDF5 file
<code>dim2start</code>	2nd dimension coordinate of the first image note that the order of dimensions is reversed in the file 2nd dimension = N dimension in area detector = dimension 0 of the three-dimensional HDF dataset set to 0 if dimension may not be present
<code>dim2count</code>	number of subsequent images to average set to 1 if dimension may not be present
<code>dim3start</code>	3rd dimension coordinate of the first image note that the order of dimensions is reversed in the file 3rd dimension = extra X dimension in area detector = dimension 0 of the four-dimensional HDF dataset set to 0 if dimension may not be present
<code>dim3count</code>	number of subsequent images to average set to 1 if dimension may not be present

Definition at line 937 of file `pearl-area-import.ipf`.

#### 11.6.2.8 variable `adh5_load_detector_slabs` ( variable `fileID`, string `detectorpath`, variable `progress` = `defaultValue` )

load the detector dataset from the open HDF5 file.

the function loads the dataset image by image using the hyperslab option. this function gives the same result as `adh5_load_detector`. it is about 5% slower, and it depends on HDF5 Browser code. but it does not choke on large datasets (as long as the final wave fits into memory).

##### Parameters

<code>fileID</code>	ID of open HDF5 file from <code>HDF5OpenFile</code> .
<code>detectorpath</code>	path to detector group in the HDF5 file.
<code>progress</code>	1 (default): show progress window; 0: do not show progress window.

##### Returns

0 if successful, non-zero if an error occurred.

Definition at line 772 of file `pearl-area-import.ipf`.

### 11.6.2.9 string adh5\_load\_info ( string *APathName*, string *AFileName* )

load descriptive info from a HDF5 file created by the Area Detector software.

the information returned is the array size and active scans

#### Attention

EXPERIMENTAL this function should be merged with adh5\_load\_preview

#### Parameters

<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in <i>FileName</i> or a dialog box should be displayed
<i>AFileName</i>	if empty a dialog box shows up

Definition at line 489 of file pearl-area-import.ipf.

### 11.6.2.10 string adh5\_load\_preview ( string *ANickName*, string *APathName*, string *AFileName*, variable *load\_data* = default**Value**, variable *load\_attr* = default**Value** )

load a single image from a HDF5 file created by the Area Detector software.

the data wave is loaded into the current data folder. attributes are loaded into the attr subfolder. existing waves in attr are deleted.

#### Warning

EXPERIMENTAL this function uses the root:pearl\_area:preview data folder. existing data there may be deleted!

#### Parameters

<i>ANickName</i>	destination wave name. the wave is created in the current data folder.
<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in <i>FileName</i> or a dialog box should be displayed
<i>AFileName</i>	if empty a dialog box shows up
<i>load_data</i>	1 (default): load data; 0: do not load data
<i>load_attr</i>	1 (default): load attributes; 0: do not load attributes note: for correct scaling of the image, the attributes need to be loaded

Definition at line 379 of file pearl-area-import.ipf.

### 11.6.2.11 string adh5\_load\_reduced ( string *ANickName*, string *APathName*, string *AFileName*, funcref *reduction\_func*, string *reduction\_param*, variable *load\_data* = default**Value**, variable *load\_attr* = default**Value**, variable *progress* = default**Value** )

load and reduce a dataset from a HDF5 file created by the Area Detector software.

the resulting dataset is reduced in one image dimension by a user-defined reduction function, e.g. by region-of-interest integration, curve fitting, etc.

the function loads the dataset image by image using the hyperslab option and applies a custom reduction function to each image. the results from the reduction function are composed into one result wave. the raw data are discarded.

if the data is from the electron analyser driver and some special attributes are included, the function will set the

scales of the image dimensions.

#### Parameters

<i>ANickName</i>	destination folder name (top level under root)
<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in <i>FileName</i> or a dialog box should be displayed
<i>AFileName</i>	if empty a dialog box shows up
<i>reduction_func</i>	custom reduction function (any user-defined function which has the same parameters as <a href="#">adh5_default_reduction()</a> )
<i>reduction_param</i>	parameter string for the reduction function
<i>load_data</i>	1 (default): load data; 0: do not load data
<i>load_attr</i>	1 (default): load attributes; 0: do not load attributes for proper wave scaling, the attributes must be loaded
<i>progress</i>	1 (default): show progress window; 0: do not show progress window

Definition at line 294 of file `pearl-area-import.ipf`.

**11.6.2.12** `variable adh5_load_reduced_detector ( variable fileID, string detectorpath, funcref reduction_func, string reduction_param, variable progress = defaultValue, variable nthreads = defaultValue )`

load a reduced detector dataset from the open HDF5 file.

the function loads the dataset image by image using the hyperslab option and applies a custom reduction function to each image. the results from the reduction function are composed into one result wave. the raw data are discarded.

by default, the reduction function is called in separate threads to reduce the total loading time. (see the global variable `adh5_perf_secs` which reports the total run time of the function.) the effect varies depending on the balance between file loading (image size) and data processing (complexity of the reduction function). for debugging the reduction function, multi-threading can be disabled.

#### Parameters

<i>fileID</i>	ID of open HDF5 file from <code>HDF5OpenFile</code>
<i>detectorpath</i>	path to detector group in the HDF5 file
<i>reduction_func</i>	custom reduction function (any user-defined function which has the same parameters as <a href="#">adh5_default_reduction()</a> )
<i>reduction_param</i>	parameter string for the reduction function
<i>progress</i>	1 (default): show progress window; 0: do not show progress window
<i>nthreads</i>	-1 (default): use as many threads as there are processor cores (in addition to main thread) 0: use main thread only (e.g. for debugging the reduction function) $\geq 1$ : use a fixed number of (additional) threads

Definition at line 1162 of file `pearl-area-import.ipf`.

**11.6.2.13** `variable adh5_loadattr_all ( variable fileID, string attributespath )`

load an `NDAAttributes` group from an open HDF5 file into the current data folder.

datasets contained in the group are loaded as waves. if a dataset contains only one data point, it is added to the IN, ID, IV, IU waves, where IN = EPICS channel name, ID = attribute name, IV = value, IU = unit (units are left empty as they are not saved in HDF5). attributes of the `NDAAttributes` group are added to the IN, ID, IV, IU waves, however, IN and IU are left empty as this information is not saved in the HDF5 file.

## Parameters

<i>fileID</i>	ID of open HDF5 file from HDF5OpenFile
<i>attributespath</i>	path to NDAttributes group in the HDF5 file

Definition at line 1479 of file pearl-area-import.ipf.

11.6.2.14 variable `adh5_redim ( wave data )`

redimension a multi-dimensional area detector array loaded from HDF5.

so that the image dimensions are X and Y singleton dimensions are removed (most common cases only)

in the redimensioned array, the original dimension type is noted in the dimension label: AD\_Dim0 = first image dimension AD\_Dim1 = second image dimension AD\_DimN = frame sequence AD\_DimX = extra dimension X AD\_DimY = extra dimension Y (cannot be loaded in Igor)

## Parameters

<i>data</i>	area detector data loaded from HDF5 to be redimensioned
-------------	---

Definition at line 618 of file pearl-area-import.ipf.

11.6.2.15 variable `adh5_scale ( wave data, string source = defaultValue )`

set the dimension scales of an area detector dataset.

the intrinsic dimensions 0 and 1 are scaled according to the data source (currently supported: Prosilica cameras, Scienta electron analyser). the extra dimensions are scaled according to the scan. the latter requires that the positioner names and position values are available.

Definition at line 718 of file pearl-area-import.ipf.

11.6.2.16 variable `adh5_scale_scan ( wave data )`

scales the extra dimensions of an area detector dataset according to the EPICS scan

the scan positioner name and its values must be available

**Todo** incomplete

Definition at line 1719 of file pearl-area-import.ipf.

11.6.2.17 variable `adh5_scale_scienta ( wave data )`

set the energy and angle scales of an area detector dataset from the Scienta analyser.

the dimension labels of the energy and angle scales must be set correctly: AD\_Dim0 = energy dimension; AD\_Dim1 = angle dimension. these dimensions must be the first two dimensions of a multi-dimensional dataset. normally, AD\_Dim0 is the X dimension, and AD\_Dim1 the Y dimension.

Definition at line 1614 of file pearl-area-import.ipf.

11.6.2.18 threadsafe variable `adh5_setup_profile ( wave image, wave profile, variable dim )`

set up a one-dimensional wave for a line profile based on a 2D original wave.

redimensions the profile wave to the given dimension. copies the scale and dimension label of the given dimension.  
Definition at line 1113 of file pearl-area-import.ipf.

**11.6.2.19** `string adh5_test_reduction_func ( wave source, wave dest1, wave dest2, funcref reduction_func, string reduction_param )`

wrapper function for testing reduction functions from the command line.

Igor does not allow global variables as pass-by-reference parameter for reduction\_param.

Definition at line 1128 of file pearl-area-import.ipf.

**11.6.2.20** `static variable BeforeFileOpenHook ( variable refNum, string fileName, string path, string type, string creator, variable kind ) [static]`

callback function for drag&drop of HDF5 files into Igor.

Definition at line 34 of file pearl-area-import.ipf.

**11.6.2.21** `static dfr GetAttrDataFolderDFR ( wave data ) [static]`

find the attributes data folder of an area detector dataset.

since version 1.04 attributes should be stored in a subfolder named attr. earlier versions had the attributes in the same data folder as the actual dataset.

#### Parameters

<i>data</i>	wave containing the main dataset.
-------------	-----------------------------------

#### Returns

data folder reference of the attributes folder. the reference may be invalid (and default to root) if the folder cannot be found, cf. built-in DataFolderRefStatus function.

Definition at line 699 of file pearl-area-import.ipf.

**11.6.2.22** `static variable read_attribute_info ( string datawavename, string source, variable * idest ) [static]`

sub-function of adh5\_loadattr\_all.

reads one attribute from a wave which was loaded from an HDF5 file into the info waves IN, ID, IV, IU. the attribute is read only if the input wave contains exactly one item, i.e. either the measurement is a single image, or the attribute has string type.

#### Parameters

<i>datawavename</i>	name of the attribute wave in the current folder. can be text or numeric.
<i>source</i>	source identifier (EPICS name) of the attribute.
<i>idest</i>	destination index in IN, ID, IV, IU where the results are written. the variable is incremented if data was written, otherwise it is left unchanged. make sure IN, ID, IV, IU have at least idest + 1 elements.

Definition at line 1561 of file pearl-area-import.ipf.

11.6.2.23 `static threadsafe variable reduce_slab_image ( wave slabdata, wave image, wave profile1, wave profile2, funcref reduction_func, string reduction_param ) [static]`

Definition at line 1444 of file `pearl-area-import.ipf`.

11.6.2.24 `static threadsafe variable reduce_slab_worker ( funcref reduction_func ) [static]`

Definition at line 1405 of file `pearl-area-import.ipf`.

## 11.7 `pearl-area-profiles.ipf` File Reference

profile extraction for multi-dimensional datasets acquired from area detectors.

### Namespaces

- [PearlAreaProfiles](#)

*profile extraction for multi-dimensional datasets acquired from area detectors.*

### Functions

- threadsafe wave [ad\\_extract\\_rod](#) (wave dataset, variable x1, variable x2, variable y1, variable y2, variable z1, variable z2, string destname, variable noavg=defaultValue, variable sdev=defaultValue, variable pscale=defaultValue)  
*1D cut through 3D dataset, integrate in normal dimensions*
- threadsafe wave [ad\\_extract\\_rod\\_x](#) (wave dataset, variable q1, variable q2, variable r1, variable r2, string destname, variable noavg=defaultValue, variable sdev=defaultValue)  
*1D cut through 3D dataset along X dimension.*
- threadsafe wave [ad\\_extract\\_rod\\_y](#) (wave dataset, variable p1, variable p2, variable r1, variable r2, string destname, variable noavg=defaultValue, variable sdev=defaultValue)  
*1D cut through 3D dataset along Y dimension.*
- threadsafe wave [ad\\_extract\\_rod\\_z](#) (wave dataset, variable p1, variable p2, variable q1, variable q2, string destname, variable noavg=defaultValue, variable sdev=defaultValue)  
*1D cut through 3D dataset along Z dimension.*
- threadsafe wave [ad\\_extract\\_slab](#) (wave dataset, variable x1, variable x2, variable y1, variable y2, variable z1, variable z2, string destname, variable noavg=defaultValue, variable pscale=defaultValue)  
*2D cut through 3D dataset, integrate in normal dimension*
- threadsafe wave [ad\\_extract\\_slab\\_x](#) (wave dataset, variable p1, variable p2, string destname, variable noavg=defaultValue)
- threadsafe wave [ad\\_extract\\_slab\\_y](#) (wave dataset, variable q1, variable q2, string destname, variable noavg=defaultValue)
- threadsafe wave [ad\\_extract\\_slab\\_z](#) (wave dataset, variable r1, variable r2, string destname, variable noavg=defaultValue)
- threadsafe wave [ad\\_profile\\_x](#) (wave dataset, variable q1, variable q2, string destname, variable noavg=defaultValue)  
*1D cut through 2D dataset along X dimension, new destination wave.*
- threadsafe wave [ad\\_profile\\_x\\_w](#) (wave dataset, variable q1, variable q2, wave destwave, variable noavg=defaultValue)  
*1D cut through 2D dataset along X dimension, existing destination wave.*
- threadsafe wave [ad\\_profile\\_y](#) (wave dataset, variable p1, variable p2, string destname, variable noavg=defaultValue)  
*1D cut through 2D dataset along Y dimension, new destination wave.*

- threadsafe wave `ad_profile_y_w` (wave dataset, variable `p1`, variable `p2`, wave `destwave`, variable `noavg=defaultValue`)  
*1D cut through 2D dataset along X dimension, existing destination wave.*
- threadsafe variable `calc_y_profile_mins` (wave image)
- variable `ad_collect_multiscan_y` (wave dataset, wave positions, wave `destwave`, variable `noavg=defaultValue`)  
*collect profiles from a multi-scan.*

### 11.7.1 Detailed Description

profile extraction for multi-dimensional datasets acquired from area detectors.

#### Author

matthias muntwiler, [matthias.muntwiler@psi.ch](mailto:matthias.muntwiler@psi.ch)

#### Copyright

2013-15 Paul Scherrer Institut  
Licensed under the Apache License, Version 2.0 (the "License");  
you may not use this file except in compliance with the License.  
You may obtain a copy of the License at <http://www.apache.org/licenses/LICENSE-2.0>

### 11.7.2 Function Documentation

**11.7.2.1** variable `ad_collect_multiscan_y` ( wave *dataset*, wave *positions*, wave *destwave*, variable *noavg* = `defaultValue` )

collect profiles from a multi-scan.

#### Warning

experimental: name and interface of this function may change.

Definition at line 621 of file `pearl-area-profiles.ipf`.

**11.7.2.2** threadsafe wave `ad_extract_rod` ( wave *dataset*, variable *x1*, variable *x2*, variable *y1*, variable *y2*, variable *z1*, variable *z2*, string *destname*, variable *noavg* = `defaultValue`, variable *sdev* = `defaultValue`, variable *pscale* = `defaultValue` )

1D cut through 3D dataset, integrate in normal dimensions

#### Parameters

<i>dataset</i>	
<i>x1,x2,y1,y2,z1,z2</i>	coordinates of integration region by default, the coordinates use wave scaling coordinates of rod dimensions (to be preserved) must be nan coordinate pairs don't have to be ordered, i.e. both $x1 \leq x2$ and $x1 \geq x2$ are allowed.
<i>destname</i>	name of destination wave. to be created in current data folder. if empty, the function returns a free wave
<i>noavg</i>	non-zero = calculate the sum, default = 0 as of version 1.05, this option should rather be called "calc_sum" or similar. it is noavg for compatibility with older code.
<i>sdev</i>	non-zero = calculate the standard deviation, default = 0 by default, the function calculates the average of the integration region set either the noavg or sdev option to select the sum or the standard deviation, respectively. if both options are set, noavg (sum) takes precedence.
<i>pscale</i>	scaling of the slab coordinates $x1, x2, \dots, z2$ : zero or default = wave scaling, non-zero = print scaling

## Remarks

- version 1.02: the specification of the destination coordinates has changed
- version 1.04: the function returns an empty wave reference if an error occurred

Definition at line 54 of file pearl-area-profiles.ipf.

**11.7.2.3** `threadsafe wave ad_extract_rod_x ( wave dataset, variable q1, variable q2, variable r1, variable r2, string destname, variable noavg = defaultValue, variable sdev = defaultValue )`

1D cut through 3D dataset along X dimension.

see [ad\\_extract\\_rod\(\)](#) for descriptions of common parameters.

Definition at line 106 of file pearl-area-profiles.ipf.

**11.7.2.4** `threadsafe wave ad_extract_rod_y ( wave dataset, variable p1, variable p2, variable r1, variable r2, string destname, variable noavg = defaultValue, variable sdev = defaultValue )`

1D cut through 3D dataset along Y dimension.

see [ad\\_extract\\_rod\(\)](#) for descriptions of common parameters.

Definition at line 168 of file pearl-area-profiles.ipf.

**11.7.2.5** `threadsafe wave ad_extract_rod_z ( wave dataset, variable p1, variable p2, variable q1, variable q2, string destname, variable noavg = defaultValue, variable sdev = defaultValue )`

1D cut through 3D dataset along Z dimension.

see [ad\\_extract\\_rod\(\)](#) for descriptions of common parameters.

Definition at line 231 of file pearl-area-profiles.ipf.

**11.7.2.6** `threadsafe wave ad_extract_slab ( wave dataset, variable x1, variable x2, variable y1, variable y2, variable z1, variable z2, string destname, variable noavg = defaultValue, variable pscale = defaultValue )`

2D cut through 3D dataset, integrate in normal dimension

## Parameters

<i>dataset</i>	
<i>x1,x2,y1,y2,z1,z2</i>	coordinates of integration region. by default, the coordinates use wave scaling. coordinates of slab dimensions (to be preserved) must be nan. coordinate pairs don't have to be ordered, i.e. both $x1 \leq x2$ and $x1 \geq x2$ are allowed. coordinates can be out of range (-inf and +inf allowed) to select the whole range.
<i>destname</i>	name of destination wave. to be created in current data folder. if empty, the function returns a free wave.
<i>noavg</i>	zero or default = average, non-zero = sum.
<i>pscale</i>	scaling of the slab coordinates $x1, x2, \dots, z2$ : zero or default = wave scaling, non-zero = point scaling.

## Remarks

- version 1.02: the specification of the destination coordinates has changed
- version 1.04: the function returns an empty wave reference if an error occurred

Definition at line 313 of file pearl-area-profiles.ipf.

**11.7.2.7** `threadsafe wave ad_extract_slab_x ( wave dataset, variable p1, variable p2, string destname, variable noavg = defaultValue )`

Definition at line 358 of file pearl-area-profiles.ipf.

**11.7.2.8** `threadsafe wave ad_extract_slab_y ( wave dataset, variable q1, variable q2, string destname, variable noavg = defaultValue )`

Definition at line 398 of file pearl-area-profiles.ipf.

**11.7.2.9** `threadsafe wave ad_extract_slab_z ( wave dataset, variable r1, variable r2, string destname, variable noavg = defaultValue )`

Definition at line 438 of file pearl-area-profiles.ipf.

**11.7.2.10** `threadsafe wave ad_profile_x ( wave dataset, variable q1, variable q2, string destname, variable noavg = defaultValue )`

1D cut through 2D dataset along X dimension, new destination wave.

Definition at line 480 of file pearl-area-profiles.ipf.

**11.7.2.11** `threadsafe wave ad_profile_x_w ( wave dataset, variable q1, variable q2, wave destwave, variable noavg = defaultValue )`

1D cut through 2D dataset along X dimension, existing destination wave.

Definition at line 504 of file pearl-area-profiles.ipf.

**11.7.2.12** `threadsafe wave ad_profile_y ( wave dataset, variable p1, variable p2, string destname, variable noavg = defaultValue )`

1D cut through 2D dataset along Y dimension, new destination wave.

Definition at line 542 of file pearl-area-profiles.ipf.

**11.7.2.13** `threadsafe wave ad_profile_y_w ( wave dataset, variable p1, variable p2, wave destwave, variable noavg = defaultValue )`

1D cut through 2D dataset along X dimension, existing destination wave.

Definition at line 567 of file pearl-area-profiles.ipf.

**11.7.2.14** `threadsafe variable calc_y_profile_mins ( wave image )`

Definition at line 603 of file pearl-area-profiles.ipf.

## 11.8 pearl-arpes.ipf File Reference

data acquisition and analysis package for ARPES at PEARL.

```
#include "pearl-area-display"
#include "pearl-area-profiles"
#include "pearl-area-import"
#include "pearl-pshell-import"
#include "pearl-data-explorer"
#include "pearl-anglescan-process"
#include "pearl-anglescan-tracker"
#include "pearl-scienta-preprocess"
#include "pearl-elog"
```

## Namespaces

- [PearlArpes](#)

*data acquisition and analysis package for ARPES at PEARL.*

## Functions

- static variable [AfterCompiledHook](#) ()  
*initializes package data once when the procedure is first loaded*
- variable [UnloadPearlArpesPackage](#) ()

### 11.8.1 Detailed Description

data acquisition and analysis package for ARPES at PEARL.

this procedure defines the PEARL ARPES package the main purpose of this file is to load the necessary dependent procedures (see the include statements at the top)

#### Precondition

- HDF5 XOP must be loaded.
- on-line data acquisition functionality requires the EPICS XOP to be loaded

#### Author

matthias muntwiler, [matthias.muntwiler@psi.ch](mailto:matthias.muntwiler@psi.ch)

#### Copyright

2012-15 Paul Scherrer Institut  
Licensed under the Apache License, Version 2.0 (the "License");  
you may not use this file except in compliance with the License.  
You may obtain a copy of the License at <http://www.apache.org/licenses/LICENSE-2.0>

### 11.8.2 Function Documentation

#### 11.8.2.1 static variable [AfterCompiledHook](#) ( ) [static]

initializes package data once when the procedure is first loaded

Definition at line 81 of file `pearl-arpes.ipf`.

11.8.2.2 variable `UnloadPearlArpesPackage ( )`

Definition at line 88 of file `pearl-arpes.ipf`.

## 11.9 pearl-data-explorer.ipf File Reference

preview and import panel for PEARL data

```
#include "pearl-area-import"
#include "pearl-area-profiles"
#include "pearl-area-display"
#include "pearl-pshell-import"
```

### Namespaces

- [PearlDataExplorer](#)  
*preview and import panel for PEARL data*

### Functions

- variable [pearl\\_data\\_explorer \( \)](#)
- static variable [init\\_package \( \)](#)  
*initialize the global variables of the data explorer.*
- static variable [save\\_prefs \( \)](#)
- static variable [load\\_prefs \( \)](#)
- static variable [pearl\\_file\\_type](#) (string filename)  
*check whether a file can be imported by this module.*
- static variable [update\\_filelist \( \)](#)  
*read a list of PEARL files from the file system*
- static variable [update\\_datasets \( \)](#)
- static variable [preview\\_file](#) (string filename)
- static wave [preview\\_pshell\\_file](#) (string filename)  
*load the preview of a PShell HDF5 file.*
- static wave [preview\\_hdf\\_file](#) (string filename)  
*load the preview of a PEARL HDF5 file.*
- static wave [preview\\_itx\\_file](#) (string filename)
- static wave [preview\\_mtrx\\_file](#) (string filename)  
*load the preview of a Matrix STM file.*
- static variable [extract\\_preview\\_image](#) (wave data, wave preview)
- static variable [preview\\_dataset](#) (string datasetname)
- static variable [preview\\_datafolder \( \)](#)
- static variable [preview\\_setscale\\_x](#) (wave data, wave preview)
- static variable [preview\\_attributes](#) (dfref attr\_folder, dfref dest\_folder=defaultValue, wave attr\_filter=defaultValue, variable include\_datawaves=defaultValue, variable include\_infowaves=defaultValue)
- static variable [display\\_dataset](#) (string datasetname)
- variable [test\\_attributes\\_notebook \( \)](#)
- static variable [attributes\\_notebook](#) (wave attr\_names, wave attr\_values, string title)
- static variable [notebook\\_add\\_attributes](#) (string notebook\_name, wave attr\_filter, wave attr\_names, wave attr\_values)
- static string [show\\_preview\\_graph](#) (wave data, wave xdata=defaultValue)
- static string [display\\_preview\\_trace](#) (wave xtrace, wave ytrace)

- static variable `load_selected_files` (string options=defaultValue)
- static variable `load_file` (string filename, string options=defaultValue)
- static variable `prompt_hdf_options` (string \*options)
- variable `prompt_default_process` (string \*param)  
*prototype for prompting for processing function parameters.*
- variable `prompt_func_params` (string func\_name, string \*func\_param)
- static dfr `load_pshell_file` (string filename, string options=defaultValue)
- static dfr `load_hdf_file` (string filename, string options=defaultValue)
- static dfr `load_itx_file` (string filename, string options=defaultValue)
- static dfr `load_mtrx_file` (string filename, string options=defaultValue)  
*load a matrix (STM) data file*
- string `itx_suggest_foldername` (string filename, variable ignoredate=defaultValue, string sourcename=defaultValue, variable unique=defaultValue)
- void `PearlDataExplorer` ()
- static variable `bp_load_prefs` (WMButtonAction \*ba)
- static variable `bp_save_prefs` (WMButtonAction \*ba)
- static variable `bp_browse_filepath` (WMButtonAction \*ba)
- static variable `bp_update_filelist` (WMButtonAction \*ba)
- static variable `bp_load_files` (WMButtonAction \*ba)
- static variable `bp_load_files_opt` (WMButtonAction \*ba)
- static variable `bp_file_next` (WMButtonAction \*ba)
- static variable `bp_file_prev` (WMButtonAction \*ba)
- static variable `lbp_filelist` (WMListboxAction \*lba)
- static variable `bp_update_datasets` (WMButtonAction \*ba)
- static variable `bp_dataset_folder` (WMButtonAction \*ba)
- static variable `bp_dataset_display` (WMButtonAction \*ba)
- static variable `bp_dataset_next` (WMButtonAction \*ba)
- static variable `bp_dataset_prev` (WMButtonAction \*ba)
- static variable `lbp_datasets` (WMListboxAction \*lba)
- static variable `bp_attr_notebook` (WMButtonAction \*ba)

## Variables

- static const string `package_name` = "pearl\_explorer"
- static const string `package_path` = "root:packages:pearl\_explorer:"
- static const string `ks_filematch_adh5` = "\*.h5"
- static const string `ks_filematch_pshell` = "psh\*.h5"
- static const string `ks_filematch_itx` = "\*.itx"
- static const string `ks_filematch_mtrx` = "\*\_mtrx"

### 11.9.1 Detailed Description

preview and import panel for PEARL data

preview and import panel for PEARL data:

- area detector (HDF5) files from scienta analyser and prosilica cameras (if HDF5.xop is installed).
- igor text files from s-scans and otf-scans.
- pshell (HDF5) data files (if HDF5.xop is installed).
- matrix STM files (if MatrixFileReader.xop is installed).

## 11.9.2 Function Documentation

11.9.2.1 `static variable attributes_notebook ( wave attr_names, wave attr_values, string title ) [static]`

Definition at line 805 of file pearl-data-explorer.ipf.

11.9.2.2 `static variable bp_attr_notebook ( WMButtonAction * ba ) [static]`

Definition at line 1733 of file pearl-data-explorer.ipf.

11.9.2.3 `static variable bp_browse_filepath ( WMButtonAction * ba ) [static]`

Definition at line 1419 of file pearl-data-explorer.ipf.

11.9.2.4 `static variable bp_dataset_display ( WMButtonAction * ba ) [static]`

Definition at line 1624 of file pearl-data-explorer.ipf.

11.9.2.5 `static variable bp_dataset_folder ( WMButtonAction * ba ) [static]`

Definition at line 1598 of file pearl-data-explorer.ipf.

11.9.2.6 `static variable bp_dataset_next ( WMButtonAction * ba ) [static]`

Definition at line 1644 of file pearl-data-explorer.ipf.

11.9.2.7 `static variable bp_dataset_prev ( WMButtonAction * ba ) [static]`

Definition at line 1671 of file pearl-data-explorer.ipf.

11.9.2.8 `static variable bp_file_next ( WMButtonAction * ba ) [static]`

Definition at line 1485 of file pearl-data-explorer.ipf.

11.9.2.9 `static variable bp_file_prev ( WMButtonAction * ba ) [static]`

Definition at line 1517 of file pearl-data-explorer.ipf.

11.9.2.10 `static variable bp_load_files ( WMButtonAction * ba ) [static]`

Definition at line 1457 of file pearl-data-explorer.ipf.

11.9.2.11 `static variable bp_load_files_opt ( WMButtonAction * ba ) [static]`

Definition at line 1471 of file pearl-data-explorer.ipf.

11.9.2.12 `static variable bp_load_prefs ( WMButtonAction * ba ) [static]`

Definition at line 1391 of file pearl-data-explorer.ipf.



11.9.2.24 `static dfr load_hdf_file ( string filename, string options = defaultValue ) [static]`

Definition at line 1125 of file `pearl-data-explorer.ipf`.

11.9.2.25 `static dfr load_itx_file ( string filename, string options = defaultValue ) [static]`

Definition at line 1175 of file `pearl-data-explorer.ipf`.

11.9.2.26 `static dfr load_mtrx_file ( string filename, string options = defaultValue ) [static]`

load a matrix (STM) data file

Definition at line 1218 of file `pearl-data-explorer.ipf`.

11.9.2.27 `static variable load_prefs ( ) [static]`

Definition at line 116 of file `pearl-data-explorer.ipf`.

11.9.2.28 `static dfr load_pshell_file ( string filename, string options = defaultValue ) [static]`

Definition at line 1073 of file `pearl-data-explorer.ipf`.

11.9.2.29 `static variable load_selected_files ( string options = defaultValue ) [static]`

Definition at line 955 of file `pearl-data-explorer.ipf`.

11.9.2.30 `static variable notebook_add_attributes ( string notebook_name, wave attr_filter, wave attr_names, wave attr_values ) [static]`

Definition at line 844 of file `pearl-data-explorer.ipf`.

11.9.2.31 `variable pearl_data_explorer ( )`

Definition at line 44 of file `pearl-data-explorer.ipf`.

11.9.2.32 `static variable pearl_file_type ( string filename ) [static]`

check whether a file can be imported by this module.

the file type is determined by the extension of the file name.

#### Returns

file type

- 0 not a recognized file type
- 1 PShell file (HDF5, name starts with `psh_`)
- 2 area detector HDF5 file
- 3 Igor text (itx) file
- 4 Matrix STM file (`*_mtrx`)

Definition at line 161 of file `pearl-data-explorer.ipf`.

### 11.9.2.33 void PearlDataExplorer ( )

Definition at line 1300 of file pearl-data-explorer.ipf.

```
11.9.2.34 static variable preview_attributes ( dfref attr_folder, dfref dest_folder = default_value, wave attr_filter
      = default_value, variable include_datawaves = default_value, variable include_infowaves =
      default_value ) [static]
```

Definition at line 631 of file pearl-data-explorer.ipf.

```
11.9.2.35 static variable preview_datafolder ( ) [static]
```

Definition at line 560 of file pearl-data-explorer.ipf.

```
11.9.2.36 static variable preview_dataset ( string datasetname ) [static]
```

Definition at line 517 of file pearl-data-explorer.ipf.

```
11.9.2.37 static variable preview_file ( string filename ) [static]
```

Definition at line 255 of file pearl-data-explorer.ipf.

```
11.9.2.38 static wave preview_hdf_file ( string filename ) [static]
```

load the preview of a PEARL HDF5 file.

the preview is an arbitrary detector image extracted from the file, see [adh5\\_load\\_preview\(\)](#). the preview is loaded to the preview\_image wave in the pearl\_explorer data folder.

the s\_file\_info string is updated with information about the scan dimensions.

#### Parameters

<i>filename</i>	name of a file in the directory specified by the pearl_explorer_filepath path object.
-----------------	---

#### Returns

wave reference of the preview image

Definition at line 363 of file pearl-data-explorer.ipf.

```
11.9.2.39 static wave preview_itx_file ( string filename ) [static]
```

Definition at line 395 of file pearl-data-explorer.ipf.

```
11.9.2.40 static wave preview_mtrx_file ( string filename ) [static]
```

load the preview of a Matrix STM file.

the preview is loaded to the preview\_image wave in the pearl\_explorer data folder.

the s\_file\_info string is updated with information about the scan dimensions.

this function requires the MatrixFileReader.xop and [pearl-matrix-import.ipf](#) to be loaded. otherwise it will return an empty wave reference.

**Parameters**

<i>filename</i>	name of a file in the directory specified by the <code>pearl_explorer_filepath</code> path object.
-----------------	--

**Returns**

wave reference of the preview image. empty wave reference if the function failed.

Definition at line 435 of file `pearl-data-explorer.ipf`.

**11.9.2.41 static wave preview\_pshell\_file ( string filename ) [static]**

load the preview of a PShell HDF5 file.

the preview is an arbitrary detector image extracted from the file, see [adh5\\_load\\_preview\(\)](#). the preview is loaded to the `preview_image` wave in the `pear_explorer` data folder.

the `s_file_info` string is updated with information about the scan dimensions.

**Parameters**

<i>filename</i>	name of a file in the directory specified by the <code>pearl_explorer_filepath</code> path object.
-----------------	--

**Returns**

wave reference of the preview image

Definition at line 307 of file `pearl-data-explorer.ipf`.

**11.9.2.42 static variable preview\_setscale\_x ( wave data, wave preview ) [static]**

Definition at line 601 of file `pearl-data-explorer.ipf`.

**11.9.2.43 variable prompt\_default\_process ( string \* param )**

prototype for prompting for processing function parameters.

the function should prompt the user for function parameters, and update the `param` argument if the user clicked OK. returns 0 if the user clicked OK, 1 if the user cancelled.

prompt functions must have the same name as the corresponding reduction function with the prefix "prompt\_". be aware of the limited length of function names in Igor.

this function is a prototype. it does nothing but returns OK.

Definition at line 1053 of file `pearl-data-explorer.ipf`.

**11.9.2.44 variable prompt\_func\_params ( string func\_name, string \* func\_param )**

Definition at line 1059 of file `pearl-data-explorer.ipf`.

**11.9.2.45 static variable prompt\_hdf\_options ( string \* options ) [static]**

Definition at line 1014 of file `pearl-data-explorer.ipf`.

**11.9.2.46** static variable `save_prefs ( )` `[static]`

Definition at line 96 of file `pearl-data-explorer.ipf`.

**11.9.2.47** static string `show_preview_graph ( wave data, wave xdata = defaultValue )` `[static]`

Definition at line 871 of file `pearl-data-explorer.ipf`.

**11.9.2.48** variable `test_attributes_notebook ( )`

Definition at line 798 of file `pearl-data-explorer.ipf`.

**11.9.2.49** static variable `update_datasets ( )` `[static]`

Definition at line 212 of file `pearl-data-explorer.ipf`.

**11.9.2.50** static variable `update_filelist ( )` `[static]`

read a list of PEARL files from the file system

`wtFiles` and `wSelectedFiles` in the package data folder are updated. only files for which `pearl_file_type()` returns non-zero are listed.

Definition at line 184 of file `pearl-data-explorer.ipf`.

### 11.9.3 Variable Documentation

**11.9.3.1** const string `ks_filematch_adh5 = "*.h5"` `[static]`

Definition at line 39 of file `pearl-data-explorer.ipf`.

**11.9.3.2** const string `ks_filematch_itx = "*.itx"` `[static]`

Definition at line 41 of file `pearl-data-explorer.ipf`.

**11.9.3.3** const string `ks_filematch_mtrx = "*.mtrx"` `[static]`

Definition at line 42 of file `pearl-data-explorer.ipf`.

**11.9.3.4** const string `ks_filematch_pshell = "psh*.h5"` `[static]`

Definition at line 40 of file `pearl-data-explorer.ipf`.

**11.9.3.5** const string `package_name = "pearl_explorer"` `[static]`

Definition at line 36 of file `pearl-data-explorer.ipf`.

**11.9.3.6** const string `package_path = "root:packages:pearl_explorer:"` `[static]`

Definition at line 37 of file `pearl-data-explorer.ipf`.

## 11.10 pearl-elog.ipf File Reference

interface for writing ELOG entries with Igor graphs as attachment.

### Namespaces

- [PearlElog](#)  
*interface for writing ELOG entries with Igor graphs as attachment.*

### Functions

- variable [pearl\\_elog](#) (string logbook)  
*main function to initialize ELOG and to open an ELOG panel.*
- static variable [IgorBeforeNewHook](#) (string igorApplicationNameStr)  
*save preferences and recent values before Igor opens a new experiment.*
- static variable [IgorQuitHook](#) (string igorApplicationNameStr)  
*save preferences and recent values before Igor quits.*
- static variable [AfterFileOpenHook](#) (variable refNum, string file, string pathName, string type, string creator, variable kind)  
*initialize the package and reload preferences after an experiment is loaded.*
- static dfr [get\\_elog\\_df](#) (string name, variable category)  
*get the package, logbook, or template datafolder.*
- static variable [init\\_package](#) (variable clean=defaultValue)  
*initialize the package data folder.*
- variable [elog\\_init\\_pearl\\_templates](#) ()  
*setup PEARL template logbooks.*
- static variable [init\\_volatile\\_vars](#) ()  
*initialize volatile variables.*
- variable [elog\\_create\\_logbook](#) (string name, string template=defaultValue)  
*create a new logbook.*
- variable [elog\\_config](#) (string elog\_path=defaultValue, string hostname=defaultValue, variable port=defaultValue, string subdir=defaultValue)  
*set global module configuration parameters*
- variable [elog\\_login](#) (string logbook, string username, string password)  
*set username and password for login to a logbook*
- variable [elog\\_logout](#) (string logbook)  
*clear username and password of a logbook or all logbooks.*
- static variable [save\\_prefs](#) ()  
*save persistent package data to the preferences file.*
- static variable [load\\_prefs](#) ()  
*load persistent package data from the preferences file.*
- static string [list\\_logbooks](#) (variable templates=defaultValue)  
*get a list of configured logbooks or templates.*
- variable [elog\\_validate\\_attributes](#) (string logbook, string attributes)  
*validate attributes*
- variable [elog\\_create\\_entry](#) (string logbook, string attributes, string message, variable encoding=defaultValue, string graphs=defaultValue, variable replyto=defaultValue)  
*create a new entry in ELOG*
- variable [elog\\_add\\_attachment](#) (string logbook, variable id, string graphs)  
*add one or more graphs to an existing ELOG entry*
- static string [prepare\\_command\\_line](#) (string logbook)

- format the ELOG command and essential address arguments.*
- static string `format_url` (string logbook)
  - format the URL for display to the user*
- static string `prepare_graph_attachments` (string graphs)
  - prepare screenshots of graph windows for attachments*
- static string `get_timestamp` (string sep)
- static string `create_message_file` (string message)
  - save the message to a temporary text file*
- static string `create_graph_file` (string graphname, variable fileindex)
  - save a graph to a temporary graphics file*
- static string `create_cmd_file` (string cmd)
  - write the command line to a file.*
- static string `get_log_path` ()
- static variable `cleanup_temp_files` ()
  - delete temporary files created by the ELOG module.*
- static variable `parse_result` ()
  - parse the result file from an elog invocation.*
- string `elog_prompt_logbook` ()
  - prompt to open or create a logbook*
- variable `elog_prompt_login` (string logbook)
  - prompt the user for login to a logbook*
- string `PearlElogPanel` (string logbook)
  - open a new panel for submitting data to ELOG.*
- static variable `elog_panel_hook` (WMWinHookStruct \*s)
- static variable `update_attach_items` (string logbook)
  - update the list of attachments*
- static variable `move_attach_item` (string logbook, variable item, variable distance)
  - move an attachment item in the list of attachments*
- static variable `bp_attach_updown` (WMButtonAction \*ba)
  - button procedure for the attachment up and down buttons*
- static variable `bp_submit` (WMButtonAction \*ba)
  - button procedure for the Submit and Reply buttons*
- static variable `bp_attach_top` (WMButtonAction \*ba)
  - select top graph window for attachment*
- static variable `bp_attach_allnone` (WMButtonAction \*ba)
  - select/deselect all graph windows for attachment*
- static variable `bp_attach` (WMButtonAction \*ba)
- static variable `bp_save_graphs` (WMButtonAction \*ba)
- static variable `bp_clear` (WMButtonAction \*ba)
- static variable `bp_login` (WMButtonAction \*ba)
- static variable `bp_logout` (WMButtonAction \*ba)
- static string `get_default_panel_name` ()
- static string `get_panel_attributes` (string windowname)
  - get a list of attributes from the fields of the ELOG panel.*
- static string `set_panel_attributes` (string windowname, string attributes, variable clear=defaultValue)
  - set the fields of the ELOG panel*
- static string `get_panel_message` (string windowname)
  - get the message field of the ELOG panel*
- static string `set_panel_message` (string windowname, string message)
  - set the message field of the ELOG panel*
- static string `get_panel_graphs` (string windowname)
  - get the names of the graphs selected for attachment*
- static string `set_panel_graphs` (string windowname, string graphs)
  - update selection of graphs for attachment*

## Variables

- static const string `package_name` = "pearl\_elog"
- static const string `package_path` = "root:packages:pearl\_elog:"
- static const variable `kdfRoot` = 0
- static const variable `kdfVolatile` = 1
- static const variable `kdfPersistent` = 2
- static const variable `kdfTemplates` = 3
- static const string `elog_success_msg` = "Message successfully transmitted"
- static const string `elog_parse_id` = "ID=%u"
- static const variable `kAttachColSel` = 0
- static const variable `kAttachColTitle` = 1
- static const variable `kAttachColName` = 2

### 11.10.1 Detailed Description

interface for writing ELOG entries with Igor graphs as attachment.

the functions in this module support the following ELOG features:

- submit new entries and replies to existing entries.
- text field, list box, and check box attributes.
- attach any Igor graph to ELOG.
- configurable logbook templates for logbooks that share the same configuration.
- common server configurations available on the ELOG command line (hostname, port, SSL, username, password, sub-directory).
- not specific to the configuration at PEARL. PEARL code is concentrated in the `elog_init_pearl_templates()` function.
- the configuration of the ELOG server and logbooks as well as the most recently used attributes are persisted in the preference file.

usage:

1. the administrator of the ELOG server creates logbook templates according to the configuration of the logbooks. the templates are written in Igor code.
2. the user opens logbooks via the *Open ELOG panel* menu item. before first use, select a template and enter a name for the logbook. the new logbook is written to the preference file, and can afterwards be opened directly.
3. if the server requires a user name and password, click the login button.
4. edit the message, attributes and attachments as necessary, and submit to ELOG.
5. log out before saving the experiment to clear the password.

## Attention

the user name and password are stored in the global data tree of an experiment. it is not possible to handle passwords safely in Igor. they can be read by anyone having access to an open Igor experiment or a saved experiment file (unless the password is reset before saving). therefore:

- use a password for the ELOG server which is different from your other passwords.
- clear the password (logout button in the panel) before saving an experiment.

elog command line

```

*  elog -h <hostname> [-p port] [-d subdir]
*                                     Location where elogd is running
*  -l logbook                         Name of logbook
*  -s                                 Use SSL for communication
*  [-v]                              For verbose output
*  [-w password]                     Write password defined on server
*  [-u username password]            User name and password
*  [-f <attachment>]                 Up to 50 attachments
*  -a <attribute>=<value>             Up to 50 attributes
*  [-r <id>]                         Reply to existing message
*  [-q]                              Quote original text on reply
*  [-e <id>]                         Edit existing message
*  [-x]                              Suppress email notification
*  [-n 0|1|2]                       Encoding: 0:ELcode,1:plain,2:HTML
*  -m <textfile>] | <text>
*

```

#### Author

matthias muntwiler, [matthias.muntwiler@psi.ch](mailto:matthias.muntwiler@psi.ch)

#### Copyright

2013-17 Paul Scherrer Institut  
 Licensed under the Apache License, Version 2.0 (the "License");  
 you may not use this file except in compliance with the License.  
 You may obtain a copy of the License at <http://www.apache.org/licenses/LICENSE-2.0>

### 11.10.2 Function Documentation

**11.10.2.1 static variable AfterFileOpenHook ( variable refNum, string file, string pathName, string type, string creator, variable kind ) [static]**

initialize the package and reload preferences after an experiment is loaded.

Definition at line 143 of file pearl-elog.ipf.

**11.10.2.2 static variable bp\_attach ( WMBUTTONACTION \* ba ) [static]**

Definition at line 1634 of file pearl-elog.ipf.

**11.10.2.3 static variable bp\_attach\_allnone ( WMBUTTONACTION \* ba ) [static]**

select/deselect all graph windows for attachment

Definition at line 1613 of file pearl-elog.ipf.

**11.10.2.4 static variable bp\_attach\_top ( WMBUTTONACTION \* ba ) [static]**

select top graph window for attachment

Definition at line 1597 of file pearl-elog.ipf.

**11.10.2.5 static variable bp\_attach\_updown ( WMBUTTONACTION \* ba ) [static]**

button procedure for the attachment up and down buttons

Definition at line 1518 of file pearl-elog.ipf.

11.10.2.6 static variable `bp_clear ( WMBUTTONACTION * ba )` [static]

Definition at line 1695 of file `pearl-elog.ipf`.

11.10.2.7 static variable `bp_login ( WMBUTTONACTION * ba )` [static]

Definition at line 1711 of file `pearl-elog.ipf`.

11.10.2.8 static variable `bp_logout ( WMBUTTONACTION * ba )` [static]

Definition at line 1729 of file `pearl-elog.ipf`.

11.10.2.9 static variable `bp_save_graphs ( WMBUTTONACTION * ba )` [static]

Definition at line 1667 of file `pearl-elog.ipf`.

11.10.2.10 static variable `bp_submit ( WMBUTTONACTION * ba )` [static]

button procedure for the Submit and Reply buttons

Definition at line 1550 of file `pearl-elog.ipf`.

11.10.2.11 static variable `cleanup_temp_files ( )` [static]

delete temporary files created by the ELOG module.

this deletes all temporary graph files that are referenced by the volatile `temp_graph_files` list. `temp_graph_files` is a semicolon-delimited string. items are added by [create\\_graph\\_file\(\)](#).

this function should be called before a new experiment is loaded or igor quits.

Definition at line 1115 of file `pearl-elog.ipf`.

11.10.2.12 static string `create_cmd_file ( string cmd )` [static]

write the command line to a file.

the command script changes the working directory to the Temporary directory. it also deletes a previous `elog.log` file.

#### Note

somewhere the command line (even inside command files) is limited to 1024 bytes. for this reason all files should now be in the Temporary directory assigned by igor.

Definition at line 1059 of file `pearl-elog.ipf`.

11.10.2.13 static string `create_graph_file ( string graphname, variable fileindex )` [static]

save a graph to a temporary graphics file

the file is saved to the Temporary directory returned by igor's `SpecialDirPath` function. the file name contains a time stamp and the specified file index to make it unique. the function returns the name of the file (excluding path!)

the full path is added to the `temp_graph_files` global list. a hook function will delete the files listed there when igor quits.

**Parameters**

<i>graphname</i>	object name of the graph to save.
<i>fileindex</i>	incremental index of the file within one submission. the file name is made unique by a time stamp and this file index. submissions within the same second must have a unique file index.

**Returns**

(string) name of the created file. empty string if unsuccessful.

**Version**

1.41 the return value has changed from full path to file name only due to the limited length of the command line (1024 bytes).

Definition at line 1023 of file pearl-elog.ipf.

**11.10.2.14** `static string create_message_file ( string message ) [static]`

save the message to a temporary text file

the file is saved to the Temporary directory returned by igor's SpecialDirPath function under the file name "elog\_↵temp\_message.txt". the function returns the name of the file (excluding path!)

**Note**

percent characters (%) cannot be passed to elog. they are removed silently from the message.

**Parameters**

<i>message</i>	text message to save to the file.
----------------	-----------------------------------

**Returns**

(string) name of the created file. empty string if unsuccessful.

**Version**

1.41 the return value has changed from full path to file name only due to the limited length of the command line (1024 bytes).

Definition at line 982 of file pearl-elog.ipf.

**11.10.2.15** `variable elog_add_attachment ( string logbook, variable id, string graphs )`

add one or more graphs to an existing ELOG entry

**Parameters**

<i>logbook</i>	name of the target logbook
<i>id</i>	identification number of the existing entry
<i>graphs</i>	names of graph windows to be added as attachments, semicolon separated

**Warning**

this will delete all existing attachments of the entry!

Definition at line 792 of file pearl-elog.ipf.

**11.10.2.16** `variable elog_config ( string elog_path = defaultValue, string hostname = defaultValue, variable port = defaultValue, string subdir = defaultValue )`

set global module configuration parameters

Definition at line 473 of file pearl-elog.ipf.

**11.10.2.17** `variable elog_create_entry ( string logbook, string attributes, string message, variable encoding = defaultValue, string graphs = defaultValue, variable replyto = defaultValue )`

create a new entry in ELOG

this is the main function to create a new entry in a logbook.

**Parameters**

<i>logbook</i>	name of the target logbook.
<i>attributes</i>	key=value list of attributes, semicolon separated.
<i>message</i>	free text part of the entry.
<i>encoding</i>	encoding of message, 0:ELcode, 1:plain (default), 2:HTML.
<i>graphs</i>	names of graph windows to be added as attachments, semicolon separated.
<i>replyto</i>	existing message ID (> 1) to follow up on. 0 or default: start new thread.

**Returns**

ID number of the new entry (> 0), or error code (< 0).

- -1: failed to save temporary message file.
- -2: invalid/missing command line.
- -3: invalid/missing attributes.
- -4: elog returned error

Definition at line 685 of file pearl-elog.ipf.

**11.10.2.18** `variable elog_create_logbook ( string name, string template = defaultValue )`

create a new logbook.

create a new empty logbook or duplicate from a template.

**Parameters**

<i>name</i>	name of the new logbook. if the logbook exists, the existing logbook folder is killed and replaced by a new one. this may fail if a window is still open.
<i>template</i>	name of the template. if empty string, a new empty logbook is created.

Definition at line 414 of file pearl-elog.ipf.

**11.10.2.19 variable `elog_init_pearl_templates ( )`**

setup PEARL template logbooks.

template logbooks for PEARL.

**Remarks**

this function is specific to the setup at PEARL.

Definition at line 262 of file `pearl-elog.ipf`.

**11.10.2.20 variable `elog_login ( string logbook, string username, string password )`**

set username and password for login to a logbook

the username and password are stored (in plain text) in global strings under the selected logbook folder. this is necessary for sending data to the ELOG server.

call `elog_logout()` to clear the password variables and to avoid unintended use of your credentials.

**Warning**

igor does not have a built-in mechanism to protect passwords. user names and passwords are stored in plain text in the data folder tree. as such they are saved to experiment files and preferences.

**Parameters**

<i>logbook</i>	name of the target logbook.
----------------	-----------------------------

Definition at line 513 of file `pearl-elog.ipf`.

**11.10.2.21 variable `elog_logout ( string logbook )`**

clear username and password of a logbook or all logbooks.

the username and password are stored (in plain text) in global strings under the selected logbook folder. this function resets the username and password strings.

**Parameters**

<i>logbook</i>	name of the target logbook. if empty, the passwords of all logbooks are cleared.
----------------	--

Definition at line 533 of file `pearl-elog.ipf`.

**11.10.2.22 static variable `elog_panel_hook ( WMWinHookStruct * s )` [static]**

Definition at line 1394 of file `pearl-elog.ipf`.

**11.10.2.23 string `elog_prompt_logbook ( )`**

prompt to open or create a logbook

Definition at line 1184 of file `pearl-elog.ipf`.

**11.10.2.24** variable `elog_prompt_login` ( *string logbook* )

prompt the user for login to a logbook

Definition at line 1214 of file `pearl-elog.ipf`.

**11.10.2.25** variable `elog_validate_attributes` ( *string logbook*, *string attributes* )

validate attributes

**Returns**

0 if all required attributes are present and enumerated items are correct. non-zero if a violation is detected.

**Todo** function currently not implemented, always returns 0

Definition at line 653 of file `pearl-elog.ipf`.

**11.10.2.26** static `string format_url` ( *string logbook* ) [static]

format the URL for display to the user

**Parameters**

<i>logbook</i>	name of the target logbook
----------------	----------------------------

Definition at line 902 of file `pearl-elog.ipf`.

**11.10.2.27** static `string get_default_panel_name` ( ) [static]

Definition at line 1746 of file `pearl-elog.ipf`.

**11.10.2.28** static `dfr get_elog_dfr` ( *string name*, *variable category* ) [static]

get the package, logbook, or template datafolder.

**Parameters**

<i>name</i>	name of logbook or template, or empty string for respective parent folder.
<i>category</i>	parameter category: <ul style="list-style-type: none"> <li>• <code>kdfRoot</code> package root</li> <li>• <code>kdfVolatile</code> volatile</li> <li>• <code>kdfPersistent</code> persistent</li> <li>• <code>kdfTemplates</code> template</li> </ul>

**Returns**

data folder reference

Definition at line 170 of file `pearl-elog.ipf`.

**11.10.2.29** `static string get_log_path ( ) [static]`

Definition at line 1095 of file pearl-elog.ipf.

**11.10.2.30** `static string get_panel_attributes ( string windowname ) [static]`

get a list of attributes from the fields of the ELOG panel.

#### Parameters

<i>windowname</i>	window name of the ELOG panel if empty, use default name "PearlElogPanel"
-------------------	---

#### Returns

list of attributes to in the format "key1=value1;key2=value2".

Definition at line 1759 of file pearl-elog.ipf.

**11.10.2.31** `static string get_panel_graphs ( string windowname ) [static]`

get the names of the graphs selected for attachment

#### Parameters

<i>windowname</i>	panel window name
-------------------	-------------------

#### Returns

a semicolon-separated list, or the empty string if the selection is not valid.

Definition at line 1914 of file pearl-elog.ipf.

**11.10.2.32** `static string get_panel_message ( string windowname ) [static]`

get the message field of the ELOG panel

#### Parameters

<i>windowname</i>	window name of the ELOG panel if empty, use default name "PearlElogPanel"
-------------------	---

#### Returns

message text

Definition at line 1868 of file pearl-elog.ipf.

**11.10.2.33** `static string get_timestamp ( string sep ) [static]`

Definition at line 958 of file pearl-elog.ipf.

**11.10.2.34** static variable `IgorBeforeNewHook ( string igorApplicationNameStr )` [static]

save preferences and recent values before Igor opens a new experiment.

Definition at line 127 of file `pearl-elog.ipf`.

**11.10.2.35** static variable `IgorQuitHook ( string igorApplicationNameStr )` [static]

save preferences and recent values before Igor quits.

Definition at line 135 of file `pearl-elog.ipf`.

**11.10.2.36** static variable `init_package ( variable clean = defaultValue )` [static]

initialize the package data folder.

the data folder is initialized with a default, local configuration without any logbooks. the server configuration should be set in the preferences.

#### Parameters

<i>clean</i>	decides what to do if the package configuration exists. <ul style="list-style-type: none"><li>• 0 (default) keep existing configuration.</li><li>• 1 overwrite existing configuration.</li></ul>
--------------	--

Definition at line 217 of file `pearl-elog.ipf`.

**11.10.2.37** static variable `init_volatile_vars ( )` [static]

initialize volatile variables.

create and initialize all volatile variables for the configured notebooks. values of existing variables are not changed.

this function must be called after new logbooks have been configured, specifically by [elog\\_create\\_logbook\(\)](#) and [load\\_prefs\(\)](#).

Definition at line 340 of file `pearl-elog.ipf`.

**11.10.2.38** static string `list_logbooks ( variable templates = defaultValue )` [static]

get a list of configured logbooks or templates.

this is list of data folder names under persistent:logbooks (or persistent:templates). the function does not check whether the folders contain valid data.

#### Parameters

<i>templates</i>	select whether logbooks (0, default) or templates (1) are returned.
------------------	---

#### Returns

semicolon-separated list of logbooks

Definition at line 618 of file `pearl-elog.ipf`.

**11.10.2.39 static variable load\_prefs ( ) [static]**

load persistent package data from the preferences file.

the preferences file is an Igor packed experiment file in a special preferences folder

Definition at line 584 of file pearl-elog.ipf.

**11.10.2.40 static variable move\_attach\_item ( string *logbook*, variable *item*, variable *distance* ) [static]**

move an attachment item in the list of attachments

Definition at line 1494 of file pearl-elog.ipf.

**11.10.2.41 static variable parse\_result ( ) [static]**

parse the result file from an elog invocation.

**Returns**

the ID of the generated message, or a value  $\leq 0$  if an error occurred.

Definition at line 1141 of file pearl-elog.ipf.

**11.10.2.42 variable pearl\_elog ( string *logbook* )**

main function to initialize ELOG and to open an ELOG panel.

this function takes care of all necessary initialization, configuration, and preferences. if a panel exists, it will be moved to the front.

**Parameters**

<i>logbook</i>	name of the logbook if empty, the user is prompted to select or create a logbook by <a href="#">elog_prompt_logbook()</a> .
----------------	---

Definition at line 97 of file pearl-elog.ipf.

**11.10.2.43 string PearlElogPanel ( string *logbook* )**

open a new panel for submitting data to ELOG.

this function creates only the panel but not the necessary data folders. call [init\\_package\(\)](#) and [load\\_prefs\(\)](#) once before creating panels.

**Parameters**

<i>logbook</i>	name of the target logbook
----------------	----------------------------

Definition at line 1241 of file pearl-elog.ipf.

**11.10.2.44 static string prepare\_command\_line ( string *logbook* ) [static]**

format the ELOG command and essential address arguments.

the following arguments are included (from global variables) if applicable: host name, port, SSL, sub-dir, username,

password the result string does not include leading or trailing space

#### Parameters

<i>logbook</i>	name of the target logbook
----------------	----------------------------

Definition at line 853 of file pearl-elog.ipf.

#### 11.10.2.45 static string prepare\_graph\_attachments ( string *graphs* ) [static]

prepare screenshots of graph windows for attachments

prepares the attachment files from Igor graph windows and returns the arguments to the elog command to attach the files. file names are returned without path. the result string does not include leading or trailing space.

#### Parameters

<i>graphs</i>	names of graph windows to be added as attachments, semicolon separated
---------------	--

Definition at line 939 of file pearl-elog.ipf.

#### 11.10.2.46 static variable save\_prefs ( ) [static]

save persistent package data to the preferences file.

saves everything under the persistent folder of the package.

Definition at line 564 of file pearl-elog.ipf.

#### 11.10.2.47 static string set\_panel\_attributes ( string *windowname*, string *attributes*, variable *clear* = defaultValue ) [static]

set the fields of the ELOG panel

#### Parameters

<i>windowname</i>	window name of the ELOG panel if empty, use default name "PearlElogPanel"
<i>attributes</i>	list of attributes to set (format "key1=value1;key2=value2")
<i>clear</i>	what to do if a key is missing in attributes? <ul style="list-style-type: none"> <li>• 0 (default) leave the field unchanged</li> <li>• 1 clear the field</li> </ul>

Definition at line 1806 of file pearl-elog.ipf.

#### 11.10.2.48 static string set\_panel\_graphs ( string *windowname*, string *graphs* ) [static]

update selection of graphs for attachment

#### Parameters

<i>windowname</i>	panel window name. looks for default panel if empty.
<i>graphs</i>	semicolon-separated list of names of graph windows to select for attachment.

Definition at line 1954 of file `pearl-elog.ipf`.

**11.10.2.49** `static string set_panel_message ( string windowname, string message )` `[static]`

set the message field of the ELOG panel

#### Parameters

<i>windowname</i>	window name of the ELOG panel if empty, use default name "PearlElogPanel"
<i>message</i>	message text that can be passed to the <code>Notebook</code> operation.

#### Returns

original message (unchanged)

Definition at line 1894 of file `pearl-elog.ipf`.

**11.10.2.50** `static variable update_attach_items ( string logbook )` `[static]`

update the list of attachments

Definition at line 1433 of file `pearl-elog.ipf`.

### 11.10.3 Variable Documentation

**11.10.3.1** `const string elog_parse_id = "ID=%u"` `[static]`

Definition at line 1134 of file `pearl-elog.ipf`.

**11.10.3.2** `const string elog_success_msg = "Message successfully transmitted"` `[static]`

Definition at line 1133 of file `pearl-elog.ipf`.

**11.10.3.3** `const variable kAttachColName = 2` `[static]`

Definition at line 1430 of file `pearl-elog.ipf`.

**11.10.3.4** `const variable kAttachColSel = 0` `[static]`

Definition at line 1428 of file `pearl-elog.ipf`.

**11.10.3.5** `const variable kAttachColTitle = 1` `[static]`

Definition at line 1429 of file `pearl-elog.ipf`.

**11.10.3.6** `const variable kdfPersistent = 2` `[static]`

Definition at line 155 of file `pearl-elog.ipf`.

11.10.3.7 `const variable kdfRoot = 0` `[static]`

Definition at line 153 of file pearl-elog.ipf.

11.10.3.8 `const variable kdfTemplates = 3` `[static]`

Definition at line 156 of file pearl-elog.ipf.

11.10.3.9 `const variable kdfVolatile = 1` `[static]`

Definition at line 154 of file pearl-elog.ipf.

11.10.3.10 `const string package_name = "pearl_elog"` `[static]`

Definition at line 85 of file pearl-elog.ipf.

11.10.3.11 `const string package_path = "root:packages:pearl_elog:"` `[static]`

Definition at line 86 of file pearl-elog.ipf.

## 11.11 pearl-fitfuncs.ipf File Reference

```
#include "mm-physconst"
```

### Data Structures

- struct [DoniachSunjicStruct](#)

### Functions

- threadsafe variable [ds1\\_bg](#) (wave w, variable x)
- threadsafe variable [ds2\\_bg](#) (wave w, variable x)
- variable [ds4\\_bg](#) (wave w, variable x)
- variable [ds6\\_bg](#) (wave w, variable x)
- threadsafe variable [DoniachSunjicBroadS](#) ([DoniachSunjicStruct](#) \*s)
- variable [DoniachSunjicBroad](#) (wave pw, wave yw, wave xw)
- variable [Calc\\_DoniachSunjicBroad](#) (wave pw, wave yw)
- variable [Fit\\_DoniachSunjicBroad](#) (wave pw, wave yw, wave xw, wave ww)
- variable [Au4f](#) (wave w, variable x)
- variable [Au4f\\_2p2](#) (wave w, variable x)
- variable [ShowComponents\\_Au4f\\_2p2](#) (wave coef\_wave, wave fit\_wave)
- variable [Au4f\\_2p3](#) (wave w, variable x)
- variable [ShowComponents\\_Au4f\\_2p3](#) (wave coef\_wave, wave fit\_wave)
- variable [FermiGaussConv](#) (wave pw, wave yw, wave xw)

*convolution of Fermi-Dirac distribution and a Gaussian.*

### Variables

- [version](#)

### 11.11.1 Function Documentation

#### 11.11.1.1 variable `Au4f` ( wave *w*, variable *x* )

Definition at line 419 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.2 variable `Au4f_2p2` ( wave *w*, variable *x* )

Definition at line 464 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.3 variable `Au4f_2p3` ( wave *w*, variable *x* )

Definition at line 549 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.4 variable `Calc_DoniachSunjicBroad` ( wave *pw*, wave *yw* )

Definition at line 362 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.5 variable `DoniachSunjicBroad` ( wave *pw*, wave *yw*, wave *xw* )

Definition at line 301 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.6 threadsafe variable `DoniachSunjicBroadS` ( `DoniachSunjicStruct` \* *s* )

Definition at line 212 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.7 threadsafe variable `ds1_bg` ( wave *w*, variable *x* )

Definition at line 38 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.8 threadsafe variable `ds2_bg` ( wave *w*, variable *x* )

Definition at line 61 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.9 variable `ds4_bg` ( wave *w*, variable *x* )

Definition at line 89 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.10 variable `ds6_bg` ( wave *w*, variable *x* )

Definition at line 130 of file `pearl-fitfuncs.ipf`.

#### 11.11.1.11 variable `FermiGaussConv` ( wave *pw*, wave *yw*, wave *xw* )

convolution of Fermi-Dirac distribution and a Gaussian.

- `pw[0]` = constant background
- `pw[1]` = linear background
- `pw[2]` = amplitude

- `pw[3]` = Fermi level in eV
- `pw[4]` = temperature in K
- `pw[5]` = gaussian width = FWHM / 1.66511

Definition at line 672 of file `pearl-fitfuncs.ipf`.

11.11.1.12 `variable Fit_DoniachSunjicBroad ( wave pw, wave yw, wave xw, wave ww )`

Definition at line 383 of file `pearl-fitfuncs.ipf`.

11.11.1.13 `variable ShowComponents_Au4f_2p2 ( wave coef_wave, wave fit_wave )`

Definition at line 518 of file `pearl-fitfuncs.ipf`.

11.11.1.14 `variable ShowComponents_Au4f_2p3 ( wave coef_wave, wave fit_wave )`

Definition at line 617 of file `pearl-fitfuncs.ipf`.

## 11.11.2 Variable Documentation

11.11.2.1 `version`

Definition at line 5 of file `pearl-fitfuncs.ipf`.

## 11.12 pearl-gui-tools.ipf File Reference

### Functions

- variable `display_progress_panel` (string *title*, string *message*, variable *progress\_max*)
- variable `update_progress_panel` (variable *progress*, string *message*=defaultValue, variable *progress\_max*↔max=defaultValue)
- variable `kill_progress_panel` ()

### 11.12.1 Function Documentation

11.12.1.1 `variable display_progress_panel ( string title, string message, variable progress_max )`

Definition at line 18 of file `pearl-gui-tools.ipf`.

11.12.1.2 `variable kill_progress_panel ( )`

Definition at line 50 of file `pearl-gui-tools.ipf`.

11.12.1.3 `variable update_progress_panel ( variable progress, string message = defaultValue, variable progress_max = defaultValue )`

Definition at line 32 of file `pearl-gui-tools.ipf`.

## 11.13 `pearl-matrix-import.ipf` File Reference

data file import for omicron matrix (STM) files

### Data Structures

- struct `errorCode`  
*from matrixfilereader help*

### Namespaces

- `PearlMatrixImport`  
*data file import for omicron matrix (STM) files*

### Functions

- static variable `init_package` ()  
*initialize the package data folder.*
- static variable `check_package_folder` ()  
*check that the package data folder exists*
- static variable `AfterFileOpenHook` (variable refNum, string file, string pathName, string type, string creator, variable kind)  
*initialize the package and reload preferences after an experiment is loaded.*
- static variable `BeforeFileOpenHook` (variable refNum, string fileName, string path, string type, string creator, variable kind)  
*open a matrix file that was dropped into Igor.*
- string `matrix_format_elog_message` (wave metadata)  
*generate elog message from bricklet metadata*
- variable `matrix_preview_2d` (wave data, wave metadata)
- static wave `preview_matrix_file` (string filename)  
*load the preview of a Matrix data file*
- static variable `initStruct` (`errorCode` \*`errorCode`)  
*from matrixfilereader help*
- variable `mtrx_load_all` ()  
*load all data from a Matrix data file.*
- variable `mtrx_parse_filename` (string fileName, string \*resultFile, variable \*runCycle, variable \*scanCycle, string \*channel)  
*parse matrix file names*
- string `mtrx_split_filename` (string fileName, string \*prefix, string \*datepart, string \*timepart)  
*split a matrix filename and return the first three parts*
- dfr `mtrx_create_folder` (string fileName, dfref df\_base=defaultValue)  
*create or look up a data folder based on a matrix file name.*
- dfr `mtrx_get_cycle_folder` (dfref df\_base=defaultValue, variable runCycle=defaultValue, variable scanCycle=defaultValue)  
*create a data folder for bricklet data.*
- variable `mtrx_file_brickletID` (string resultFile, variable runCycle, variable scanCycle, string channel)  
*find out bricklet ID of a file*
- variable `mtrx_open_file` (string pathName, string fileNameOrPath)  
*open a matrix result or data file*
- string `mtrx_load_preview` (string destName, string pathName, string fileName, string traces=defaultValue)

- load a preview image from a Matrix data file.*
- string `mtrx_load_file` (string pathName, string fileName, string traces=defaultValue)
- load all data from a Matrix data file.*
- variable `mtrx_scale_dataset` (wave data)
- string `mtrx_load_info` (string APathName, string AFileName)
- load descriptive info from a Matrix data file.*
- variable `subtract_line_bg` (wave img)
- remove linear background line-by-line*

## Variables

- static const string `package_name` = "pearl\_matrix\_import"
- static const string `package_path` = "root:packages:pearl\_matrix\_import:"
- static const string `ks_filematch_mtrx` = "\*\_mtrx"

### 11.13.1 Detailed Description

data file import for omicron matrix (STM) files

the matrix file import requires the matrix file reader XOP by thomas braun (<http://www.igorexchange.com/project/matrixFileReader>) which in turn requires an installation of vernissage by omicron nanotechnology.

#### Warning

EXPERIMENTAL the matrix import module and its interface may change radically in future revisions!

#### Author

matthias muntwiler, [matthias.muntwiler@psi.ch](mailto:matthias.muntwiler@psi.ch)

#### Copyright

2016 Paul Scherrer Institut  
 Licensed under the Apache License, Version 2.0 (the "License");  
 you may not use this file except in compliance with the License.  
 You may obtain a copy of the License at <http://www.apache.org/licenses/LICENSE-2.0>

### 11.13.2 Function Documentation

**11.13.2.1** static variable `AfterFileOpenHook` ( variable `refNum`, string `file`, string `pathName`, string `type`, string `creator`, variable `kind` ) [static]

initialize the package and reload preferences after an experiment is loaded.

Definition at line 85 of file pearl-matrix-import.ipf.

**11.13.2.2** static variable `BeforeFileOpenHook` ( variable `refNum`, string `fileName`, string `path`, string `type`, string `creator`, variable `kind` ) [static]

open a matrix file that was dropped into Igor.

preliminary implementation. this should rather load the entire file and display a preview. graph windows should be reused by subsequent loads. also decide on a data saving location.

Definition at line 102 of file pearl-matrix-import.ipf.

**11.13.2.3 static variable check\_package\_folder ( ) [static]**

check that the package data folder exists

initialize the package if the folder does not exist.

Definition at line 72 of file pearl-matrix-import.ipf.

**11.13.2.4 static variable init\_package ( ) [static]**

initialize the package data folder.

Definition at line 45 of file pearl-matrix-import.ipf.

**11.13.2.5 static variable initStruct ( errorCode \* errorCode ) [static]**

from matrixfilereader help

Definition at line 212 of file pearl-matrix-import.ipf.

**11.13.2.6 string matrix\_format\_elog\_message ( wave metadata )**

generate elog message from bricklet metadata

**Parameters**

<i>metadata</i>	two-column text wave
-----------------	----------------------

Definition at line 120 of file pearl-matrix-import.ipf.

**11.13.2.7 variable matrix\_preview\_2d ( wave data, wave metadata )**

Definition at line 145 of file pearl-matrix-import.ipf.

**11.13.2.8 dfr mtrx\_create\_folder ( string fileName, dref df\_base = defaultValue )**

create or look up a data folder based on a matrix file name.

the name of the folder is mtrx\_date\_time, where date and time are parsed from the file name. for this to work, the file name must consist of at least three parts that are separated by dash or underscore. the second (third) part contains the date (time). date and time are copied as strings.

if the data folder exists, a reference to the existing folder is returned.

**Parameters**

<i>fileName</i>	name of the result or data file.
<i>df_base</i>	(optional) base data folder. default: current folder.

**Returns**

reference of the newly created or existing data folder.

Definition at line 360 of file pearl-matrix-import.ipf.

11.13.2.9 `variable mtrx_file_brickletID ( string resultFile, variable runCycle, variable scanCycle, string channel )`

find out bricklet ID of a file

**Warning**

EXPERIMENTAL the code of this function is inefficient. the function may be removed in a later version.

**Parameters**

<i>resultFile</i>	base name of result file without chain link number and extension. as returned by <code>mtrx_parse_filename</code> .
<i>runCycle</i>	requested run cycle. 0 = first available.
<i>scanCycle</i>	requested scan cycle. 0 = first available.
<i>channel</i>	channel name. for example: "I", "Z", "Aux(V)", etc. empty string: first available.

**Returns**

bricklet ID, or -1 if an error occurred.

Definition at line 461 of file `pearl-matrix-import.ipf`.

11.13.2.10 `dfr mtrx_get_cycle_folder ( dref df_base = defaultValue, variable runCycle = defaultValue, variable scanCycle = defaultValue )`

create a data folder for bricklet data.

the name of the folder is, for example "r23s2" where the first (second) number is the run (scan) cycle. run cycle and scan cycle numbers are taken from the open matrix file unless overridden by optional arguments.

if the data folder exists, a reference to the existing folder is returned. if one of the run or scan cycle numbers is lower than 1, the base folder is returned.

**Parameters**

<i>df_base</i>	(optional) base data folder. default: current folder.
<i>runCycle</i>	(optional) run cycle number. must be $\geq 1$ . default: from last <code>mtrx_open_file</code> call.
<i>scanCycle</i>	(optional) scan cycle number. must be $\geq 1$ . default: from last <code>mtrx_open_file</code> call.

**Returns**

reference of the newly created or existing data folder.

Definition at line 405 of file `pearl-matrix-import.ipf`.

11.13.2.11 `variable mtrx_load_all ( )`

load all data from a Matrix data file.

Definition at line 231 of file `pearl-matrix-import.ipf`.

11.13.2.12 `string mtrx_load_file ( string pathName, string fileName, string traces = defaultValue )`

load all data from a Matrix data file.

the data wave is loaded into a sub-subfolder the current data folder. the relative path has the format ":mtrx\_{date}\_{time}:r{run\_cycle}s{scan\_cycle}", where the parameters {date}, {time}, {run\_cycle} and {scan\_cycle} are copied from the file name. the file name must be formatted according to the specifications set out below.

#### Parameters

<i>pathName</i>	igor symbolic path name. can be empty if the path is specified in FileName or a dialog box should be displayed
<i>fileName</i>	if empty a dialog box shows up the file name must adhere to the format "{prefix}-{date}-{time}-{anything}-{run_cycle}_{scan_cycle}.{extension}". the first three separators can alternatively be underscores. it may be necessary to change the configuration of the Matrix application.
<i>traces</i>	(currently not used) semicolon-separated list of preferred traces. the items of the list are match strings for the Igor StringMatch function. only matching traces are loaded from the file. default: "*Up;*Down;*ReUp;*ReDown;"

#### Returns

semicolon-separated list of loaded waves including partial path from current data folder.

Definition at line 767 of file pearl-matrix-import.ipf.

#### 11.13.2.13 string mtrx\_load\_info ( string APathName, string AFileName )

load descriptive info from a Matrix data file.

the info string lists the following information for each scan contained in the file:

- path of the scan group inside the file.
- number of scan positions.
- dataset names of scan positioners.
- dataset names of detectors.

#### Parameters

<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in AFileName or a dialog box should be displayed
<i>AFileName</i>	if empty a dialog box shows up

#### Returns

newline terminated string.

Definition at line 863 of file pearl-matrix-import.ipf.

#### 11.13.2.14 string mtrx\_load\_preview ( string destName, string pathName, string fileName, string traces = defaultValue )

load a preview image from a Matrix data file.

the data wave is loaded into the current data folder.

## Parameters

<i>destName</i>	destination wave name. the wave is created in the current data folder.
<i>pathName</i>	igor symbolic path name. can be empty if the path is specified in <i>FileName</i> or a dialog box should be displayed
<i>fileName</i>	if empty a dialog box shows up the file name must adhere to the format "{prefix}-{date}-{time}-{anything}--{run_cycle}_{scan_cycle}.{extension}". the first three separators can alternatively be underscores. it may be necessary to change the configuration of the Matrix application.
<i>traces</i>	(currently not used) semicolon-separated list of preferred traces. the items of the list are match strings for the Igor StringMatch function. only the first matching trace is loaded from the file. default: "*Up;*Down;*ReUp;*ReDown;"

## Returns

semicolon-separated list of loaded waves including partial path from current data folder.

Definition at line 682 of file pearl-matrix-import.ipf.

11.13.2.15 variable mtrx\_open\_file ( string *pathName*, string *fileNameOrPath* )

open a matrix result or data file

this function opens a matrix result file (.mtrx) or data file (\*.mtrx).

if a data file is selected, the function locates the corresponding result file, opens it, and looks up the bricklet ID of the data file. if a result file is selected, the function opens it but does not look up bricklet IDs.

the result file remains open and can be accessed using the mtrx\_ functions or MFR\_ operations. once a result file is open, you can easily access any bricklets linked to it, i.e., any run cycle, scan cycle, and channel.

the function stores information about the opened file in a global package data folder. if the same result file is opened again later, the information is reused and the file not read again. this may cause problems if the file has been modified in the meantime, or if the cached data become corrupt for some reason. the function detects if a data file is not linked in the open result file, and updates the cache. in other situations it may be necessary to force a reload.

**Todo** fix possible cache issues, add an option to override the cache.

## Parameters

<i>pathName</i>	igor path name or empty string.
<i>fileName</i>	file name, with or without path, or empty string.

## Returns

file type

- 0 result file (logbook)
- 1 result data file (bricklet)
- -1 error, no data loaded
- -2 matrixfilereader.xop not installed

Definition at line 545 of file pearl-matrix-import.ipf.

**11.13.2.16** `variable mtrx_parse_filename ( string fileName, string * resultFile, variable * runCycle, variable * scanCycle, string * channel )`

parse matrix file names

parse matrix file names for result name, run cycle, scan cycle, and channel.

#### Parameters

<i>fileName</i>	matrix result or data file name (without path).
<i>resultFile</i>	(out) base name of the result file. append "_%04u.mtrx" to get the actual result file. we do not know the chain link number at this stage.
<i>runCycle</i>	(out) run cycle number. necessary to look up the bricklet ID.
<i>scanCycle</i>	(out) scan cycle number. necessary to look up the bricklet ID.
<i>channel</i>	(out) channel name.

#### Returns

file type

- 0 result file (logbook)
- 1 result data file (bricklet)

result file names look like: default\_2015Apr20-124353\_STM-STM\_AtomManipulation\_0001.mtrx, default\_↵  
2015Apr20-124353\_STM-STM\_AtomManipulation\_0002.mtrx, etc. the function returns the first part up to the  
experiment name ("AtomManipulation" in the examples). all other return values set to defaults and must not  
be regarded.

result data files look like: default\_2015Apr20-124353\_STM-STM\_AtomManipulation-136\_1.Aux1(V)\_mtrx,  
default\_2015Apr20-124353\_STM-STM\_AtomManipulation-136\_1.I(V)\_mtrx, default\_2015Apr20-124353\_STM-  
STM\_AtomManipulation-14\_1.I\_mtrx, default\_2015Apr20-124353\_STM-STM\_AtomManipulation-14\_1.Z\_mtrx,  
etc. the function returns all results as described in the parameter list.

Definition at line 294 of file pearl-matrix-import.ipf.

**11.13.2.17** `variable mtrx_scale_dataset ( wave data )`

Definition at line 829 of file pearl-matrix-import.ipf.

**11.13.2.18** `string mtrx_split_filename ( string fileName, string * prefix, string * datepart, string * timepart )`

split a matrix filename and return the first three parts

we assume that the second (third) part contains the date (time). the parts are separated by dash or underscore.

Definition at line 332 of file pearl-matrix-import.ipf.

**11.13.2.19** `static wave preview_matrix_file ( string filename ) [static]`

load the preview of a Matrix data file

the preview is loaded to the preview\_image wave in the pearl\_explorer data folder.

the s\_file\_info string is updated with information about the scan dimensions.

#### Parameters

<i>filename</i>	name of a file in the directory specified by the pearl_explorer_filepath path object.
-----------------	---

**Returns**

wave reference of the preview image

Definition at line 170 of file pearl-matrix-import.ipf.

**11.13.2.20 variable subtract\_line\_bg ( wave *img* )**

remove linear background line-by-line

Definition at line 886 of file pearl-matrix-import.ipf.

**11.13.3 Variable Documentation****11.13.3.1 const string ks\_filematch\_mtrx = " \*\_mtrx" [static]**

Definition at line 40 of file pearl-matrix-import.ipf.

**11.13.3.2 const string package\_name = "pearl\_matrix\_import" [static]**

Definition at line 37 of file pearl-matrix-import.ipf.

**11.13.3.3 const string package\_path = "root:packages:pearl\_matrix\_import:" [static]**

Definition at line 38 of file pearl-matrix-import.ipf.

**11.14 pearl-menu.ipf File Reference****Functions**

- string [PearlMenuEnableFunc](#) (string funcname)
- variable [LoadPearlOptics](#) ()
- variable [LoadPearlArpes](#) ()
- variable [LoadPearlPreparation](#) ()
- variable [Display2dProfiles](#) ()
- variable [Display3dSlicer](#) ()
- variable [DisplayGizmoSlicer](#) ()
- variable [PearlLiveDisplay](#) (string epicsname, string nickname, string wbRGB)
- variable [PearlAnglescanTracker](#) (string epicsname, string wbRGB)

**11.14.1 Function Documentation****11.14.1.1 variable Display2dProfiles ( )**

Definition at line 108 of file pearl-menu.ipf.

**11.14.1.2 variable Display3dSlicer ( )**

Definition at line 125 of file pearl-menu.ipf.

#### 11.14.1.3 variable DisplayGizmoSlicer ( )

Definition at line 144 of file pearl-menu.ipf.

#### 11.14.1.4 variable LoadPearlArpes ( )

Definition at line 96 of file pearl-menu.ipf.

#### 11.14.1.5 variable LoadPearlOptics ( )

Definition at line 89 of file pearl-menu.ipf.

#### 11.14.1.6 variable LoadPearlPreparation ( )

Definition at line 102 of file pearl-menu.ipf.

#### 11.14.1.7 variable PearlAnglescanTracker ( string *epicsname*, string *wbRGB* )

Definition at line 182 of file pearl-menu.ipf.

#### 11.14.1.8 variable PearlLiveDisplay ( string *epicsname*, string *nickname*, string *wbRGB* )

Definition at line 163 of file pearl-menu.ipf.

#### 11.14.1.9 string PearlMenuEnableFunc ( string *funcname* )

Definition at line 77 of file pearl-menu.ipf.

## 11.15 pearl-otf-import.ipf File Reference

### Functions

- variable [otf\\_load\\_itx\\_all](#) (string *pathname*)
- variable [otf\\_load\\_itx\\_match](#) (string *pathname*, string *matchstr*)
- variable [otf\\_load\\_itx](#) (string *pathname*, string *filename*)
- variable [otf\\_gather\\_iterator](#) (dfref *df*, string *\*sdata*)
- variable [otf\\_gather\\_batch](#) (string *ywavematch*, string *xwavematch*, string *destfolder*)
- variable [gather\\_batch](#) (string *foldermatch*, string *ywavematch*, string *xwavematch*, string *destfolder*)
- variable [otf\\_rename\\_folders\\_iterator](#) (dfref *df*, string *\*sdata*)
- variable [otf\\_rename\\_folders](#) (string *pattern*, variable *unique\_index*=defaultValue, string *new\_suffix*=defaultValue, string *match\_str*=defaultValue)
- variable [otf\\_interp](#) (variable *e1*, variable *e2*, variable *npts*, variable *smo*)
- variable [otf\\_smo\\_int](#) (wave *win*, wave *wout*, wave *wpe*, variable *smo*)

### 11.15.1 Function Documentation

#### 11.15.1.1 variable gather\_batch ( string *foldermatch*, string *ywavematch*, string *xwavematch*, string *destfolder* )

Definition at line 162 of file pearl-otf-import.ipf.

11.15.1.2 variable `otf_gather_batch` ( string *ywavematch*, string *xwavematch*, string *destfolder* )

Definition at line 152 of file `pearl-otf-import.ipf`.

11.15.1.3 variable `otf_gather_iterator` ( *dfref df*, string \* *sdata* )

Definition at line 106 of file `pearl-otf-import.ipf`.

11.15.1.4 variable `otf_interp` ( variable *e1*, variable *e2*, variable *npts*, variable *smo* )

Definition at line 271 of file `pearl-otf-import.ipf`.

11.15.1.5 variable `otf_load_itx` ( string *pathname*, string *filename* )

Definition at line 76 of file `pearl-otf-import.ipf`.

11.15.1.6 variable `otf_load_itx_all` ( string *pathname* )

Definition at line 31 of file `pearl-otf-import.ipf`.

11.15.1.7 variable `otf_load_itx_match` ( string *pathname*, string *matchstr* )

Definition at line 53 of file `pearl-otf-import.ipf`.

11.15.1.8 variable `otf_rename_folders` ( string *pattern*, variable *unique\_index* = `defaultValue`, string *new\_suffix* = `defaultValue`, string *match\_str* = `defaultValue` )

Definition at line 231 of file `pearl-otf-import.ipf`.

11.15.1.9 variable `otf_rename_folders_iterator` ( *dfref df*, string \* *sdata* )

Definition at line 185 of file `pearl-otf-import.ipf`.

11.15.1.10 variable `otf_smo_int` ( wave *win*, wave *wout*, wave *wpe*, variable *smo* )

Definition at line 304 of file `pearl-otf-import.ipf`.

## 11.16 pearl-polar-coordinates.ipf File Reference

### Functions

- variable `cart2polar` (variable *xx*, variable *yy*, variable *zz*, variable \**radius*, variable \**theta*, variable \**phi*)
- variable `cart2polar_wave` (wave *in*, wave *out*)
- variable `polar2cart` (variable *radius*, variable *theta*, variable *phi*, variable \**xx*, variable \**yy*, variable \**zz*)
- variable `polar2cart_wave` (wave *in*, wave *out*)
- variable `polar_distance` (variable *polar1*, variable *azim1*, variable *polar2*, variable *azim2*)

### 11.16.1 Function Documentation

11.16.1.1 `variable cart2polar ( variable xx, variable yy, variable zz, variable * radius, variable * theta, variable * phi )`

Definition at line 10 of file `pearl-polar-coordinates.ipf`.

11.16.1.2 `variable cart2polar_wave ( wave in, wave out )`

Definition at line 36 of file `pearl-polar-coordinates.ipf`.

11.16.1.3 `variable polar2cart ( variable radius, variable theta, variable phi, variable * xx, variable * yy, variable * zz )`

Definition at line 48 of file `pearl-polar-coordinates.ipf`.

11.16.1.4 `variable polar2cart_wave ( wave in, wave out )`

Definition at line 58 of file `pearl-polar-coordinates.ipf`.

11.16.1.5 `variable polar_distance ( variable polar1, variable azim1, variable polar2, variable azim2 )`

Definition at line 69 of file `pearl-polar-coordinates.ipf`.

## 11.17 `pearl-pshell-import.ipf` File Reference

import data from PShell

```
#include <HDF5 Browser>
#include "pearl-gui-tools"
#include "pearl-area-import"
```

### Namespaces

- [PearlPShellImport](#)  
*import data from PShell*

### Functions

- variable [psh5\\_open\\_file](#) (string ANickName, string APathName, string AFileName)  
*open a HDF5 file created by the PShell data acquisition program and prepare the data folder.*
- variable [psh5\\_close\\_file](#) (variable fileID)  
*close a HDF5 file opened by psh5\_open\_file.*
- string [psh5\\_load\\_complete](#) (string ANickName, string APathName, string AFileName, variable load\_↔  
data=defaultValue, variable load\_attr=defaultValue)  
*load everything from a PShell data file.*
- string [psh5\\_load\\_preview](#) (string APathName, string AFileName, variable load\_data=defaultValue, variable load\_attr=defaultValue, string pref\_scans=defaultValue, string pref\_datasets=defaultValue)  
*load a preview image from a PShell data file.*
- string [psh5\\_load\\_scan\\_complete](#) (variable fileID, string scanpath, variable load\_data=defaultValue, variable load\_attr=defaultValue)

- load all data of a selected scan from a PShell data file.*
- string `psh5_list_scans` (variable fileID)
  - list scan groups of a PShell data file.*
- string `psh5_list_scan_datasets` (variable fileID, string scanpath, variable include\_regions=defaultValue)
  - list datasets of a PShell scan group.*
- string `psh5_list_scan_regions` (variable fileID, string scanpath)
  - list regions of a PShell scan group.*
- string `psh5_load_scan_data` (variable fileID, string scanpath)
  - load all datasets of a PShell scan group.*
- string `psh5_load_scan_attrs` (variable fileID, string scanpath, variable attr\_sets=defaultValue)
  - load attributes of a PShell scan group.*
- string `psh5_load_scan_meta` (variable fileID, string scanpath)
  - load metadata of a PShell scan group.*
- string `psh5_load_dataset` (variable fileID, string scanpath, string datasetname, variable set\_scale=default↵Value)
  - load a dataset from an open PShell HDF5 file.*
- static string `select_dataset` (string file\_datasets, string pref\_datasets)
  - select the preferred dataset from a list of available datasets.*
- string `psh5_load_scan_preview` (variable fileID, string scanpath, variable set\_scale=defaultValue, string pref\_datasets=defaultValue)
  - load a preview dataset from an open PShell HDF5 file.*
- string `psh5_load_scan_section` (variable fileID, string scanpath, variable dim, variable set\_scale=default↵Value, string pref\_datasets=defaultValue)
  - load a longitudinal section of a scan from an open PShell HDF5 file.*
- variable `psh5_load_dataset_meta` (variable fileID, string datapath, string datasetname, wave datawave)
  - load metadata of a PShell dataset.*
- string `psh5_load_dataset_slabs` (variable fileID, string datapath, string datasetname, variable progress=default↵Value)
  - load a dataset slab-wise from the open PShell HDF5 file.*
- string `psh5_load_dataset_slab` (variable fileID, string datapath, string datasetname, variable dim2start, variable dim2count, variable dim3start, variable dim3count)
  - load a single image from the open PShell data file.*
- variable `ps_set_dimlabels` (wave data)
  - set dimension labels according to the axis type*
- variable `ps_set_dimlabels2` (wave data, string name)
  - set dimension labels according to the axis type*
- static dfr `find_scan_folder` (dfref dataDF)
  - find the scan folder*
- static dfr `find_attr_folder` (dfref dataDF)
  - find the attributes data folder*
- variable `ps_scale_datasets` ()
  - set the dimension scales of loaded PShell Scienta datasets according to attributes.*
- variable `ps_scale_dataset` (wave data)
  - set the dimension scales of a loaded PShell Scienta dataset according to attributes.*
- static wave `find_scale_wave` (string name, dfref dataDF, dfref scanDF, dfref attrDF)
- variable `ps_detect_scale` (wave ax, wave lo, wave hi, wave un)
  - detect the dimension scales from attributes.*
- variable `ps_scale_dataset_2` (wave data, wave ax, wave lo, wave hi, wave un)
  - set the dimension scales of a dataset.*
- string `psh5_load_reduced` (string ANickName, string APathName, string AFileName, funcref reduction\_func, string reduction\_param, variable progress=defaultValue)

*load and reduce the ScientaImage dataset of the first scan of a PShell data file.*

- string `psh5_load_dataset_reduced` (variable fileID, string scanpath, string datasetname, funcref reduction\_func, string reduction\_param, variable progress=defaultValue, variable nthreads=defaultValue)

*load a reduced dataset from the open PShell HDF5 file.*

- static threadsafe variable `reduce_slab_worker` (funcref reduction\_func)
- static threadsafe variable `reduce_slab_image` (wave slabdata, wave image, wave profile1, wave profile2, funcref reduction\_func, string reduction\_param)
- string `psh5_load_info` (string APathName, string AFileName)

*load descriptive info from a PShell data file.*

- string `psh5_load_scan_info` (variable fileID, string scanpath)

*load descriptive info from a PShell scan.*

- static string `twave2list` (wave wt, string sep)

*convert text wave to list.*

- static string `wave2list` (wave w, string format, string sep)

*convert numeric wave to list.*

## Variables

- const string `kEnergyDimLabel` = "energy"

*Dimension label for the energy dispersive dimension of multi-dimensional datasets.*

- const string `kAngleDimLabel` = "angle"

*Dimension label for the angle dispersive dimension of multi-dimensional datasets.*

- const string `kScanDimLabel` = "scan"

*Dimension label for the scan dimension of multi-dimensional datasets.*

- const string `kDataDimLabel` = "data"

*Dimension label for the data dimension.*

- const string `kPreviewDatasets` = "ScientaImage;ScientaSpectrum;ImageAngleDistribution;ImageEnergyDistribution;Counts;SampleCurrent;"

*List of preferred datasets to load for preview.*

- const string `kScientaScalingDatasets` = "LensMode;ScientaChannelBegin;ScientaChannelEnd;ScientaSliceBegin;ScientaSliceEnd;"

*List of datasets that must be loaded to determine the axis scaling of a Scienta image.*

- const string `kTransposedDatasets` = "ScientaImage;"

*List of datasets that should be transposed upon loading.*

- const variable `kDetectorSensitivity` = 4

*multiply scienta detector intensity by this value to get actual counts.*

### 11.17.1 Detailed Description

import data from PShell

HDF5 file import from the PShell data acquisition program. the main import functions are:

- `psh5_load_complete()` load all scans and datasets from a file.
- `psh5_load_reduced()` load the ScientaImage dataset of the first scan and reduce its dimensionality.
- `psh5_load_scan_complete()` load all datasets of a selected scan.
- `psh5_load_scan_preview()` load a preview of a selected scan.
- `psh5_load_dataset()` load a selected dataset.
- `psh5_load_dataset_reduced()` load a selected dataset and reduce its dimensionality.

the following helper functions are also needed:

- [psh5\\_open\\_file\(\)](#)
- [psh5\\_close\\_file\(\)](#)
- [psh5\\_list\\_scans\(\)](#)
- [psh5\\_list\\_scan\\_datasets\(\)](#)
- [psh5\\_load\\_scan\\_meta\(\)](#)
- [psh5\\_load\\_scan\\_attrs\(\)](#)

## 11.17.2 Function Documentation

### 11.17.2.1 static dfr find\_attr\_folder ( dfref dataDF ) [static]

find the attributes data folder

this is the :attr folder.

Definition at line 1451 of file pearl-pshell-import.ipf.

### 11.17.2.2 static wave find\_scale\_wave ( string name, dfref dataDF, dfref scanDF, dfref attrDF ) [static]

Definition at line 1539 of file pearl-pshell-import.ipf.

### 11.17.2.3 static dfr find\_scan\_folder ( dfref dataDF ) [static]

find the scan folder

the scan folder is the one that contains the :attr folder the data and scan folders may refer to the same folder.

Definition at line 1434 of file pearl-pshell-import.ipf.

### 11.17.2.4 variable ps\_detect\_scale ( wave ax, wave lo, wave hi, wave un )

detect the dimension scales from attributes.

the function checks the data , scan and attributes folders for scan parameters. the results are written to the provided waves. the function is normally called by [ps\\_scale\\_datasets\(\)](#) but can also be used independently.

the current datafolder must be the data or the scan folder. the data folder contains the waves that are to be scaled. the scan folder contains the scan positions and the :attr folder.

the provided waves are redimensioned by the function, and dimension labels are set. the scale parameters can then be extracted by keyword, e.g.,

- `lo[%energy]` analyser energy dimension.
- `lo[%angle]` analyser angle dimension.
- `lo[%scan]` scan dimension.
- `lo[%data]` data dimension.

the function tries to read the following waves, in the data, scan, and attributes folders, where the first folder in the list takes precedence. it may fall back to more or less reasonable default values if no data is not found.

- `LensMode`

- ScientaChannelBegin
- ScientaChannelEnd
- ScientaSliceBegin
- ScientaSliceEnd
- ScanWritables
- wave referenced by ScanWritables[0]

#### Parameters

<i>ax</i>	text wave to receive the axis labels.
<i>lo</i>	wave to receive the lower limits.
<i>hi</i>	wave to receive the upper limits.
<i>un</i>	text wave to receive the unit labels.

#### Returns

the function results are written to the lo, hi, un, and ax waves.

#### Version

this function supports regions from version 1.03. check that you're in the correct data folder!

Definition at line 1596 of file pearl-pshell-import.ipf.

#### 11.17.2.5 variable `ps_scale_dataset ( wave data )`

set the dimension scales of a loaded PShell Scienta dataset according to attributes.

the current datafolder must contain the :attr folder. the data wave can be in the current folder or a sub-folder.

the dimension labels of the dataset waves must have been set correctly, e.g. by [ps\\_set\\_dimlabels\(\)](#). this is implicitly done by the high-level load functions.

the function is useful if a single dataset is loaded and scaled. if multiple datasets are loaded, [ps\\_scale\\_datasets\(\)](#) is slightly more efficient.

#### Parameters

<i>data</i>	data wave to be scaled. dimension labels (index -1) must be set correctly, cf. <a href="#">ps_set_dimlabels()</a> .
-------------	---

#### Version

this function supports regions from version 1.03.

Definition at line 1525 of file pearl-pshell-import.ipf.

#### 11.17.2.6 variable `ps_scale_dataset_2 ( wave data, wave ax, wave lo, wave hi, wave un )`

set the dimension scales of a dataset.

the function is normally called by [ps\\_scale\\_datasets\(\)](#) but can also be used independently. the limits and units must be given as function arguments with proper dimension labels.

the provided limit and unit waves must have dimension labels matching the -1 index dimension labels of the data wave, such as set by the [ps\\_detect\\_scale\(\)](#) function. the scale parameters are extracted by keyword, e.g.,

- `lo[%energy]` analyser energy dimension.
- `lo[%angle]` analyser angle dimension.
- `lo[%scan]` scan dimension.
- `lo[%data]` data dimension.

#### Parameters

<i>data</i>	data wave to be scaled. dimension labels (index -1) must be set to match the limit waves.
<i>ax</i>	axis labels. the axis labels are written to the wave note in the format <code>AxisLabel%s=%s</code> where X, Y, Z, D is substituted for the first place holder and the label for the second one.
<i>lo</i>	lower limits. the lower limits are applied using the SetScale operation.
<i>hi</i>	upper limits. the upper limits are applied using the SetScale operation.
<i>un</i>	unit labels. the unit labels are applied using the SetScale operation.

#### Version

this function supports regions from version 1.03.

Definition at line 1750 of file pearl-pshell-import.ipf.

#### 11.17.2.7 variable `ps_scale_datasets ( )`

set the dimension scales of loaded PShell Scienta datasets according to attributes.

datasets listed in the ScanReadables waves are scaled according to the attribute waves in the data, scan, and attributes folders, whichever is found first.

the current datafolder must contain the ScanReadables wave and the :attr folder. the ScanReadables text wave contains names of the waves to scale. wave names can include a relative path to a sub-folder. the path separator is "/".

the dimension labels of the dataset waves must have been set correctly, e.g. by [ps\\_set\\_dimlabels\(\)](#). this is implicitly done by the high-level load functions.

#### Version

this function supports regions from version 1.03. check that you're in the correct data folder!

Definition at line 1479 of file pearl-pshell-import.ipf.

#### 11.17.2.8 variable `ps_set_dimlabels ( wave data )`

set dimension labels according to the axis type

this function asserts a particular ordering of dimensions types based on the name of the wave for ScientaImage, ScientaSpectrum, ImageAngleDistribution, ImageEnergyDistribution. all other waves must be one-dimensional, and the dimension must be the scan dimension.

dimension labels are required by scaling functions.

#### Parameters

<i>data</i>	data wave as loaded from PShell file
-------------	--------------------------------------

**Returns**

- 0 all labels set successfully.
- 1 unidentified data source.
- 2 wave does not contain data.

Definition at line 1358 of file pearl-pshell-import.ipf.

#### 11.17.2.9 variable `ps_set_dimlabels2` ( wave *data*, string *name* )

set dimension labels according to the axis type

same as [ps\\_set\\_dimlabels\(\)](#) except that the dimension labels are set according to a separate name argument instead of the wave name.

**Parameters**

<i>data</i>	data wave as loaded from PShell file.
<i>name</i>	original name of the dataset in the PShell file.

**Returns**

- 0 all labels set successfully.
- 1 unidentified data source.
- 2 wave does not contain data.

Definition at line 1377 of file pearl-pshell-import.ipf.

#### 11.17.2.10 variable `psh5_close_file` ( variable *fileID* )

close a HDF5 file opened by `psh5_open_file`.

this function just closes the HDF5 file. no change is made to the loaded data.

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
---------------	--

Definition at line 132 of file pearl-pshell-import.ipf.

#### 11.17.2.11 string `psh5_list_scan_datasets` ( variable *fileID*, string *scanpath*, variable *include\_regions* = `defaultValue` )

list datasets of a PShell scan group.

the function returns a list of all datasets of the selected scan. this does not include datasets from the attributes sub-group.

**Note**

in a future version, an option may be introduced to filter datasets by function (*Readable* and/or *Writable*).

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>scanpath</i>	path to the scan group in the HDF5 file, e.g. "/scan 1".

**Returns**

semicolon-separated list of dataset paths.

**Version**

since version 1.03 this function returns paths relative to scanpath.

Definition at line 432 of file pearl-pshell-import.ipf.

**11.17.2.12 string psh5\_list\_scan\_regions ( variable fileID, string scanpath )**

list regions of a PShell scan group.

the function returns a list of all region groups of the selected scan.

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>scanpath</i>	path to the scan group in the HDF5 file, e.g. "/scan 1".

**Returns**

semicolon-separated list of datagroup paths.

Definition at line 473 of file pearl-pshell-import.ipf.

**11.17.2.13 string psh5\_list\_scans ( variable fileID )**

list scan groups of a PShell data file.

the function returns a list of all top-level groups whose name starts with "scan".

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
---------------	--

**Returns**

semicolon-separated list of group paths.

Definition at line 397 of file pearl-pshell-import.ipf.

**11.17.2.14 string psh5\_load\_complete ( string ANickName, string APathName, string AFileName, variable load\_data = defaultvalue, variable load\_attr = defaultvalue )**

load everything from a PShell data file.

**Parameters**

<i>ANickName</i>	destination folder name (top level under root)
<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in FileName or a dialog box should be displayed
<i>AFileName</i>	if empty a dialog box shows up

## Parameters

<i>load_data</i>	select whether datasets (positioners and detectors) are loaded. <ul style="list-style-type: none"> <li>• 1 (default) load data.</li> <li>• 0 do not load data.</li> </ul>
<i>load_attr</i>	select whether attributes (auxiliary device readbacks) are loaded. for proper wave scaling, the attributes must be loaded. <ul style="list-style-type: none"> <li>• 1 (default) load attributes.</li> <li>• 0 do not load attributes.</li> </ul>

## Returns

complete path of the loaded file if successful. empty string otherwise.  
global string `s_filepath` in new data folder contains the full file path on disk.  
global string `s_scanpaths` in new data folder contains a list of scan groups inside the file.

Definition at line 162 of file `pearl-pshell-import.ipf`.

```
11.17.2.15 string psh5_load_dataset ( variable fileID, string scanpath, string datasetname, variable set_scale =
      defaultValue )
```

load a dataset from an open PShell HDF5 file.

if the dataset has a maximum of two dimensions, the function loads it at once. if it has more than two dimension, the function calls [psh5\\_load\\_dataset\\_slabs\(\)](#) to load the data slab by slab.

- the metadata (HDF5 attributes) are loaded into the wave note, cf. [psh5\\_load\\_dataset\\_meta\(\)](#).
- dimension labels are set according the dataset name, cf. [ps\\_set\\_dimlabels\(\)](#).
- wave scaling is set if the necessary scan attributes have been loaded and the `set_scale` option is selected (default). the attributes must be loaded by [psh5\\_load\\_scan\\_meta\(\)](#) and [psh5\\_load\\_scan\\_attrs\(\)](#) (`attr_sets=2`).

the dataset is loaded into the current data folder unless `datasetname` contains a region specifier. in the latter case, the dataset is loaded into sub-folder with the name of the region. the function returns from the original data folder.

## Parameters

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>scanpath</i>	path to the scan group in the HDF5 file, e.g. <code>"/scan 1"</code> .
<i>datasetname</i>	name of the dataset. the name of the loaded wave is a cleaned up version of the dataset name. the name can include the region name as a relative path, e.g. <code>"region1/ScientaSpectrum"</code> . in this case, the dataset is loaded into a sub-folder named <code>"region1"</code> .
<i>set_scale</i>	by default, the function tries to set the wave scaling if the attributes have been loaded. if multiple datasets are loaded from a file, it is more efficient to set the scaling of all loaded datasets at the end by calling <a href="#">ps_scale_datasets()</a> . <ul style="list-style-type: none"> <li>• 1 (default) set the wave scaling.</li> <li>• 0 do not set the wave scaling.</li> </ul>

**Returns**

name of loaded wave if successful. empty string otherwise.

**Version**

this function supports regions as of version 1.03.

Definition at line 681 of file pearl-pshell-import.ipf.

### 11.17.2.16 variable psh5\_load\_dataset\_meta ( variable *fileID*, string *datapath*, string *datasetname*, wave *datawave* )

load metadata of a PShell dataset.

"metadata" are the HDF5 attributes attached to the scan dataset.

data is added to the wave note.

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>datapath</i>	path to the containing group in the HDF5 file. path separator is the slash "/".
<i>datasetname</i>	name of the dataset. may include relative path.
<i>datawave</i>	metadata is added to the wave note of this wave.

**Returns**

0 if successful, non-zero if an error occurred.

Definition at line 1066 of file pearl-pshell-import.ipf.

### 11.17.2.17 string psh5\_load\_dataset\_reduced ( variable *fileID*, string *scanpath*, string *datasetname*, funcref *reduction\_func*, string *reduction\_param*, variable *progress* = defaultValue, variable *nthreads* = defaultValue )

load a reduced dataset from the open PShell HDF5 file.

the function loads the dataset image by image using the hyperslab option and applies a custom reduction function to each image. the results from the reduction function are written to the `ReducedData1` and `ReducedData2` waves. the raw data are discarded.

by default, the reduction function is called in separate threads to reduce the total loading time. (see the global variable `psh5_perf_secs` which reports the total run time of the function.) the effect varies depending on the balance between file loading (image size) and data processing (complexity of the reduction function). for debugging the reduction function, multi-threading can be disabled.

if the reduction function requires the image waves to be scaled properly, the attributes must have been loaded by [psh5\\_load\\_scan\\_attrs\(\)](#) before. in this case, the scales of the result waves are also set by the function. otherwise, the results can also be scaled by [ps\\_scale\\_dataset\(\)](#) later.

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>scanpath</i>	path to scan group in the HDF5 file.
<i>datasetname</i>	name of the dataset. this must currently be "ScientiaImage", other data is not supported. the name of the loaded wave is a cleaned up version of the dataset name. the name can include the region name as a relative path, e.g. "region1/ScientiaImage". in this case, the dataset is loaded into a sub-folder named "region1".

## Parameters

<i>reduction_func</i>	custom reduction function (any user-defined function which has the same parameters as <a href="#">adh5_default_reduction()</a> ).
<i>reduction_param</i>	parameter string for the reduction function.
<i>progress</i>	progress window. <ul style="list-style-type: none"> <li>• 1 (default) show progress window</li> <li>• 0 do not show progress window</li> </ul>
<i>nthreads</i>	<ul style="list-style-type: none"> <li>• -1 (default) use as many threads as there are processor cores (in addition to main thread).</li> <li>• 0 use main thread only (e.g. for debugging the reduction function).</li> <li>• <math>\geq 1</math> use a fixed number of (additional) threads.</li> </ul>

## Returns

semicolon-separated list of the loaded waves, `ReducedData1` and `ReducedData2` if successful. empty string if an error occurred. error messages are printed to the history.

## Version

this function supports regions as of version 1.03.

Definition at line 1970 of file `pearl-pshell-import.ipf`.

**11.17.2.18** `string psh5_load_dataset_slab ( variable fileID, string datapath, string datasetname, variable dim2start, variable dim2count, variable dim3start, variable dim3count )`

load a single image from the open PShell data file.

the function can average over a region in the extra dimensions.

## Parameters

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>datapath</i>	path to the containing group in the HDF5 file. path separator is the slash "/".
<i>dataset</i>	name of the dataset. also defines the name of the loaded wave.
<i>dim2start</i>	2nd dimension coordinate of the first image set to 0 if dimension may not be present
<i>dim2count</i>	number of subsequent images to average set to 1 if dimension may not be present
<i>dim3start</i>	3rd dimension coordinate of the first image set to 0 if dimension may not be present
<i>dim3count</i>	number of subsequent images to average set to 1 if dimension may not be present

## Returns

name of loaded wave if successful. empty string otherwise.

Definition at line 1259 of file `pearl-pshell-import.ipf`.

11.17.2.19 `string psh5_load_dataset_slabs ( variable fileID, string datapath, string datasetname, variable progress = defaultvalue )`

load a dataset slab-wise from the open PShell HDF5 file.

the function loads the dataset image by image using the hyperslab option.

#### Parameters

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>datapath</i>	path to the containing group in the HDF5 file. path separator is the slash "/".
<i>dataset</i>	name of the dataset. also defines the name of the loaded wave.
<i>progress</i>	select whether a progress window is displayed during the process. <ul style="list-style-type: none"> <li>• 1 (default) show progress window.</li> <li>• 0 do not show progress window.</li> </ul>

#### Returns

name of loaded wave if successful. empty string otherwise.

Definition at line 1123 of file pearl-pshell-import.ipf.

11.17.2.20 `string psh5_load_info ( string APathName, string AFileName )`

load descriptive info from a PShell data file.

the info string lists the following information for each scan contained in the file:

- path of the scan group inside the file.
- number of scan positions.
- dataset names of scan positioners.
- dataset names of detectors.

#### Parameters

<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in <i>AFileName</i> or a dialog box should be displayed
<i>AFileName</i>	if empty a dialog box shows up

#### Returns

newline terminated string.

Definition at line 2286 of file pearl-pshell-import.ipf.

11.17.2.21 `string psh5_load_preview ( string APathName, string AFileName, variable load_data = defaultvalue, variable load_attr = defaultvalue, string pref_scans = defaultvalue, string pref_datasets = defaultvalue )`

load a preview image from a PShell data file.

the data wave is loaded into the current data folder. attributes are loaded into the attr subfolder. existing waves in attr are deleted.

#### Parameters

<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in <i>FileName</i> or a dialog box should be displayed
<i>AFileName</i>	if empty a dialog box shows up
<i>load_data</i>	1 (default): load data; 0: do not load data
<i>load_attr</i>	1 (default): load attributes; 0: do not load attributes note: for correct scaling of the image, the attributes need to be loaded
<i>pref_scans</i>	semicolon-separated list of preferred scans. the items of the list are match strings for the Igor StringMatch function. the first matching scan (i.e. top-level HDF5 group with a matching name) is loaded from the file. if no match is found, the first scan is loaded.
<i>pref_datasets</i>	semicolon-separated list of preferred datasets. the items of the list are match strings for the Igor StringMatch function. the first matching dataset is loaded from the file. if no match is found, the first dataset listed in the file is loaded.

#### Returns

name of loaded preview wave.

Definition at line 243 of file pearl-pshell-import.ipf.

11.17.2.22 `string psh5_load_reduced ( string ANickName, string APathName, string AFileName, funcref reduction_func, string reduction_param, variable progress = defaultValue )`

load and reduce the ScientImage dataset of the first scan of a PShell data file.

the resulting dataset is reduced in one image dimension by a user-defined reduction function, e.g. by region-of-interest integration, curve fitting, etc.

the function loads the dataset image by image using the hyperslab option and applies a custom reduction function to each image. the results from the reduction function are composed into one result wave. the raw data are discarded.

if the data is from the electron analyser driver and some special attributes are included, the function will set the scales of the image dimensions.

#### Parameters

<i>ANickName</i>	destination folder name (top level under root).
<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in <i>FileName</i> or a dialog box should be displayed.
<i>AFileName</i>	if empty a dialog box shows up.
<i>reduction_func</i>	custom reduction function (any user-defined function which has the same parameters as <a href="#">adh5_default_reduction()</a> )
<i>reduction_param</i>	parameter string for the reduction function.
<i>progress</i>	progress window. <ul style="list-style-type: none"> <li>• 1 (default) show progress window</li> <li>• 0 do not show progress window</li> </ul>

**Returns**

semicolon-separated list of the loaded waves, `ReducedData1` and `ReducedData2` if successful. empty string if an error occurred. error messages are printed to the history.  
 global string `s_filepath` in new data folder contains the full file path on disk.  
 global string `s_scanpaths` in new data folder contains a list of scan groups inside the file.

**Todo** load scan positions.

Definition at line 1846 of file `pearl-pshell-import.ipf`.

**11.17.2.23** `string psh5_load_scan_attrs ( variable fileID, string scanpath, variable attr_sets = defaultValue )`

load attributes of a PShell scan group.

"attributes" are the auxiliary data inside the `attrs` group. do not confuse with HDF5 attributes! HDF5 attributes are loaded by the `psh5_load_scan_meta()` function.

data is loaded into the current data folder. this should normally be the `:attr` folder inside the respective scan folder.

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <code>psh5_open_file()</code> .
<i>scanpath</i>	path to the scan group in the HDF5 file, e.g. <code>"/scan 1"</code> .
<i>attr_sets</i>	specify the attribute sets to be loaded. this value can be an arithmetic OR of the following constants. by default, all attributes are loaded. <ul style="list-style-type: none"> <li>• 1 all datasets that are present in the file.</li> <li>• 2 datasets relevant for wave scaling of Scienta data.</li> </ul>

**Returns**

semicolon-separated list of the loaded waves.

Definition at line 546 of file `pearl-pshell-import.ipf`.

**11.17.2.24** `string psh5_load_scan_complete ( variable fileID, string scanpath, variable load_data = defaultValue, variable load_attr = defaultValue )`

load all data of a selected scan from a PShell data file.

data is loaded into the current data folder. attribute datasets are loaded into sub-folder `attr`. region datasets are loaded into region sub-folders. existing data, if present, is overwritten.

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <code>psh5_open_file()</code> .
<i>scanpath</i>	path to the scan group in the HDF5 file, e.g. <code>"/scan 1"</code> .
<i>load_data</i>	select whether datasets (positioners and detectors) are loaded. <ul style="list-style-type: none"> <li>• 1 (default) load data.</li> <li>• 0 do not load data.</li> </ul>

**Parameters**

<i>load_attr</i>	select whether attributes (auxiliary device readbacks) are loaded. for proper wave scaling, the attributes must be loaded. <ul style="list-style-type: none"> <li>• 1 (default) load attributes.</li> <li>• 0 do not load attributes.</li> </ul>
------------------	--

**Returns**

semicolon-separated list of the loaded data waves (excluding attributes).

Definition at line 353 of file `pearl-pshell-import.ipf`.

**11.17.2.25** `string psh5_load_scan_data ( variable fileID, string scanpath )`

load all datasets of a PShell scan group.

data is loaded into the current data folder. region datasets are loaded into the respective region sub-folders.

this function does not scale the datasets. call `ps_scale_datasets()` separately.

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <code>psh5_open_file()</code> .
<i>scanpath</i>	path to the scan group in the HDF5 file, e.g. <code>"/scan 1"</code> .

**Returns**

semicolon-separated list of the loaded waves.

Definition at line 506 of file `pearl-pshell-import.ipf`.

**11.17.2.26** `string psh5_load_scan_info ( variable fileID, string scanpath )`

load descriptive info from a PShell scan.

the info string contains up to three lines which are made up of the following information:

- number of scan positions.
- dataset names of scan positioners.
- dataset names of detectors (without region names).
- region names

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <code>psh5_open_file()</code> .
<i>scanpath</i>	path to scan group in the HDF5 file.

**Returns**

newline terminated string.

Definition at line 2333 of file pearl-pshell-import.ipf.

**11.17.2.27 string psh5\_load\_scan\_meta ( variable *fileID*, string *scanpath* )**

load metadata of a PShell scan group.

*metadata* are the HDF5 attributes attached to the scan group. the following attributes are loaded. the respective wave names under Igor are given in parentheses.

- Dimensions (ScanDimensions)
- Writables (ScanWritables)
- Readables (ScanReadables)
- Steps (ScanSteps)

if they are missing in the file, *ScanDimensions* and *ScanReadables* are set to default values assuming the file contains a single spectrum.

data is loaded into the current data folder.

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>scanpath</i>	path to the scan group in the HDF5 file, e.g. <code>"/scan 1"</code> .

**Returns**

semicolon-separated list of the loaded waves.

Definition at line 614 of file pearl-pshell-import.ipf.

**11.17.2.28 string psh5\_load\_scan\_preview ( variable *fileID*, string *scanpath*, variable *set\_scale* = defaultValue, string *pref\_datasets* = defaultValue )**

load a preview dataset from an open PShell HDF5 file.

if the dataset has a maximum of two dimensions, the function loads it at once. if it has more than two dimension, the function selects and loads one two-dimensional slab.

**Parameters**

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>scanpath</i>	path to the scan group in the HDF5 file, e.g. <code>"/scan 1"</code> .
<i>set_scale</i>	by default, the function tries to set the wave scaling if the attributes have been loaded. if multiple datasets are loaded from a file, it is more efficient to set the scaling of all loaded datasets at the end by calling <a href="#">ps_scale_datasets()</a> . <ul style="list-style-type: none"> <li>• 1 (default) set the wave scaling.</li> <li>• 0 do not set the wave scaling.</li> </ul>

## Parameters

<i>pref_datasets</i>	semicolon-separated list of preferred datasets. the items of the list are match strings for the Igor StringMatch function. the first matching dataset is loaded from the file. if no match is found, the first dataset listed in the file is loaded. if empty, a hard-coded default preference list is used.
----------------------	--

## Returns

name of loaded wave if successful. empty string otherwise.

Definition at line 815 of file pearl-pshell-import.ipf.

```
11.17.2.29 string psh5_load_scan_section ( variable fileID, string scanpath, variable dim, variable set_scale =
        defaultValue, string pref_datasets = defaultValue )
```

load a longitudinal section of a scan from an open PShell HDF5 file.

the dataset must have three dimensions.

## Parameters

<i>fileID</i>	ID of open HDF5 file from <a href="#">psh5_open_file()</a> .
<i>scanpath</i>	path to the scan group in the HDF5 file, e.g. "/scan 1".
<i>dim</i>	reserved, must be 0.
<i>set_scale</i>	by default, the function tries to set the wave scaling if the attributes have been loaded. if multiple datasets are loaded from a file, it is more efficient to set the scaling of all loaded datasets at the end by calling <a href="#">ps_scale_datasets()</a> . <ul style="list-style-type: none"> <li>• 1 (default) set the wave scaling.</li> <li>• 0 do not set the wave scaling.</li> </ul>
<i>pref_datasets</i>	semicolon-separated list of preferred datasets. the items of the list are match strings for the Igor StringMatch function. the first matching dataset is loaded from the file. if no match is found, the first dataset listed in the file is loaded. if empty, a hard-coded default preference list is used.

## Returns

name of loaded wave if successful. empty string otherwise.

## Warning

EXPERIMENTAL: this function is under development.

Definition at line 928 of file pearl-pshell-import.ipf.

```
11.17.2.30 variable psh5_open_file ( string ANickName, string APathName, string AFileName )
```

open a HDF5 file created by the PShell data acquisition program and prepare the data folder.

the function opens a specified or interactively selected HDF5 file, creates a data folder \$ANickName under root, and changes to the new data folder.

the file must be closed by [psh5\\_close\\_file\(\)](#) after use.

## Parameters

<i>ANickName</i>	destination folder name (top level under root).
<i>APathName</i>	igor symbolic path name. can be empty if the path is specified in <i>FileName</i> or a dialog box should be displayed
<i>AFileName</i>	if empty a dialog box shows up

## Returns

ID of open HDF5 file from `HDF5OpenFile`. zero if an error occurred.  
 global string `s_filepath` in new data folder contains the full file path on disk.  
 global string `s_scanpaths` in new data folder contains a list of scan groups inside the file.

Definition at line 102 of file `pearl-pshell-import.ipf`.

**11.17.2.31** `static threadsafe variable reduce_slab_image ( wave slabdata, wave image, wave profile1, wave profile2, funcref reduction_func, string reduction_param ) [static]`

Definition at line 2258 of file `pearl-pshell-import.ipf`.

**11.17.2.32** `static threadsafe variable reduce_slab_worker ( funcref reduction_func ) [static]`

Definition at line 2219 of file `pearl-pshell-import.ipf`.

**11.17.2.33** `static string select_dataset ( string file_datasets, string pref_datasets ) [static]`

select the preferred dataset from a list of available datasets.

## Parameters

<i>file_datasets</i>	semicolon-separated list of datasets that are available in the file. the items may include a path separated by slashes "/". only the last component of the path is checked.
<i>pref_datasets</i>	semicolon-separated list of preferred datasets. the items of the list are match strings for the <code>Igor StringMatch</code> function. the first matching dataset is loaded from the file. if no match is found, the first file dataset is selected.

## Returns

selected dataset.

Definition at line 753 of file `pearl-pshell-import.ipf`.

**11.17.2.34** `static string twave2list ( wave wt, string sep ) [static]`

convert text wave to list.

Definition at line 2392 of file `pearl-pshell-import.ipf`.

**11.17.2.35** `static string wave2list ( wave w, string format, string sep ) [static]`

convert numeric wave to list.

Definition at line 2409 of file `pearl-pshell-import.ipf`.

### 11.17.3 Variable Documentation

#### 11.17.3.1 `const string kAngleDimLabel = "angle"`

Dimension label for the angle dispersive dimension of multi-dimensional datasets.

Definition at line 60 of file `pearl-pshell-import.ipf`.

#### 11.17.3.2 `const string kDataDimLabel = "data"`

Dimension label for the data dimension.

This label may be used to store the parameters for the `set scale d` operation.

Definition at line 67 of file `pearl-pshell-import.ipf`.

#### 11.17.3.3 `const variable kDetectorSensitivity = 4`

multiply scienta detector intensity by this value to get actual counts.

Definition at line 79 of file `pearl-pshell-import.ipf`.

#### 11.17.3.4 `const string kEnergyDimLabel = "energy"`

Dimension label for the energy dispersive dimension of multi-dimensional datasets.

Definition at line 57 of file `pearl-pshell-import.ipf`.

#### 11.17.3.5 `const string kPreviewDatasets = "ScientaImage;ScientaSpectrum;ImageAngleDistribution;ImageEnergy↵ Distribution;Counts;SampleCurrent;"`

List of preferred datasets to load for preview.

Definition at line 70 of file `pearl-pshell-import.ipf`.

#### 11.17.3.6 `const string kScanDimLabel = "scan"`

Dimension label for the scan dimension of multi-dimensional datasets.

Definition at line 63 of file `pearl-pshell-import.ipf`.

#### 11.17.3.7 `const string kScientaScalingDatasets = "LensMode;ScientaChannelBegin;ScientaChannelEnd;ScientaSlice↵ Begin;ScientaSliceEnd;"`

List of datasets that must be loaded to determine the axis scaling of a Scienta image.

Definition at line 73 of file `pearl-pshell-import.ipf`.

#### 11.17.3.8 `const string kTransposedDatasets = "ScientaImage;"`

List of datasets that should be transposed upon loading.

Definition at line 76 of file `pearl-pshell-import.ipf`.

## 11.18 pearl-scienta-preprocess.ipf File Reference

preprocessing functions for Scienta detector images.

```
#include "mm-fitfuncs"
```

### Namespaces

- [PearlScientaPreprocess](#)  
*preprocessing functions for Scienta detector images.*

### Functions

- variable [prompt\\_int\\_linbg\\_reduction](#) (string \*param)
- string [capture\\_int\\_linbg\\_cursors](#) ()
- string [csr\\_int\\_linbg\\_reduction](#) (string win)
- variable [test\\_int\\_linbg](#) (wave image)
- threadsafe variable [int\\_linbg\\_reduction](#) (wave source, wave dest1, wave dest2, string \*param)
- variable [test\\_shockley\\_anglefit](#) (wave image, variable branch)
- variable [prompt\\_Shockley\\_anglefit](#) (string \*param)
- threadsafe variable [Shockley\\_anglefit](#) (wave source, wave dest1, wave dest2, string \*param)
- variable [prompt\\_int\\_quadbg\\_reduction](#) (string \*param)
- variable [test\\_int\\_quadbg](#) (wave image)
- threadsafe variable [int\\_quadbg\\_reduction](#) (wave source, wave dest1, wave dest2, string \*param)
- variable [scienta\\_norm](#) (wave w, variable x)
- wave [fit\\_scienta\\_ang\\_transm](#) (wave data, wave params)
- threadsafe variable [scienta\\_ang\\_transm](#) (wave w, variable x)
- wave [fit\\_scienta\\_poly\\_bg](#) (wave data, wave params, variable bgterms)
- variable [scienta\\_poly\\_bg](#) (wave w, variable e, variable a)
- variable [prompt\\_redim\\_linbg\\_reduction](#) (string \*param)  
*parameter dialog for the [redim\\_linbg\\_reduction\(\)](#) function*
- threadsafe variable [redim\\_linbg\\_reduction](#) (wave source, wave dest1, wave dest2, string \*param)  
*linear background reduction function for incorrectly dimensioned scienta image*
- variable [test\\_gauss2\\_reduction](#) (wave image)  
*apply the [gauss2\\_reduction](#) function to a single image*
- variable [prompt\\_gauss2\\_reduction](#) (string \*param)  
*prompt for the [gauss2\\_reduction](#) parameters*
- threadsafe variable [gauss2\\_reduction](#) (wave source, wave dest1, wave dest2, string \*param)  
*fit horizontal cuts of an image with two gaussian peaks on a linear background*

#### 11.18.1 Detailed Description

preprocessing functions for Scienta detector images.

this procedure contains functions for data reduction and instrument-specific normalization.

#### Author

matthias muntwiler, [matthias.muntwiler@psi.ch](mailto:matthias.muntwiler@psi.ch)

#### Copyright

2013-17 Paul Scherrer Institut

Licensed under the Apache License, Version 2.0 (the "License");

you may not use this file except in compliance with the License.

You may obtain a copy of the License at <http://www.apache.org/licenses/LICENSE-2.0>

## 11.18.2 Function Documentation

### 11.18.2.1 string capture\_int\_linbg\_cursors ( )

Definition at line 66 of file pearl-scienta-preprocess.ipf.

### 11.18.2.2 string csr\_int\_linbg\_reduction ( string win )

Definition at line 76 of file pearl-scienta-preprocess.ipf.

### 11.18.2.3 wave fit\_scienta\_ang\_transm ( wave data, wave params )

Definition at line 551 of file pearl-scienta-preprocess.ipf.

### 11.18.2.4 wave fit\_scienta\_poly\_bg ( wave data, wave params, variable bgterms )

Definition at line 610 of file pearl-scienta-preprocess.ipf.

### 11.18.2.5 threadsafe variable gauss2\_reduction ( wave source, wave dest1, wave dest2, string \* param )

fit horizontal cuts of an image with two gaussian peaks on a linear background

the function fits each horizontal profile (EDC) with two gaussian peaks on a linear background. the position and width of the peaks is kept fixed according to input parameters. the peak amplitude is constrained to positive value.

the width parameter is defined as in Igor's gauss curve fit function (standard deviation divided by the square root of two). the return value in dest1 is the integrated peak either of peak 1 or peak 2. dest2 returns the corresponding error estimate.

#### Parameters

<i>source</i>	source wave. two-dimensional distribution of counts. for correct weighting and error estimation it is important that the source wave contains actual counts (Poisson statistics).
<i>dest1</i>	(out) peak area
<i>dest2</i>	(out) error estimate of peak area
<i>param</i>	<p>(in, out) semicolon-separated key=value list of processing parameters. this is a pass-by-reference argument. the following parameters are required. position, width and limit parameters are on the x (energy) scale.</p> <ul style="list-style-type: none"> <li>• rngl low limit of fit interval</li> <li>• rngl high limit of fit interval</li> <li>• pos1 position of peak 1</li> <li>• wid1 width of peak 1</li> <li>• pos2 position of peak 2</li> <li>• wid2 width of peak 2</li> <li>• return select result to return, either "int1" or "int2".</li> <li>• ybox box size of averaging in y direction, must be 1 or 3. other values lead to corrupt data.</li> </ul>

**Returns**

zero if successful, non-zero if an error occurs.

Definition at line 903 of file pearl-scienta-preprocess.ipf.

**11.18.2.6 threadsafe variable int\_linbg\_reduction ( wave *source*, wave *dest1*, wave *dest2*, string \* *param* )**

Definition at line 197 of file pearl-scienta-preprocess.ipf.

**11.18.2.7 threadsafe variable int\_quadbg\_reduction ( wave *source*, wave *dest1*, wave *dest2*, string \* *param* )**

Definition at line 457 of file pearl-scienta-preprocess.ipf.

**11.18.2.8 variable prompt\_gauss2\_reduction ( string \* *param* )**

prompt for the gauss2\_reduction parameters

useful for testing or manual processing. to debug, (temporarily) remove the threadsafe attribute from the gauss2\_reduction function.

Definition at line 834 of file pearl-scienta-preprocess.ipf.

**11.18.2.9 variable prompt\_int\_linbg\_reduction ( string \* *param* )**

Definition at line 36 of file pearl-scienta-preprocess.ipf.

**11.18.2.10 variable prompt\_int\_quadbg\_reduction ( string \* *param* )**

Definition at line 410 of file pearl-scienta-preprocess.ipf.

**11.18.2.11 variable prompt\_redim\_linbg\_reduction ( string \* *param* )**

parameter dialog for the [redim\\_linbg\\_reduction\(\)](#) function

**Parameters**

<i>param</i>	parameter string in a key1=value1;key2=value2;... list. the parameter string is passed by reference. see <a href="#">redim_linbg_reduction()</a> for a description of parameters.
--------------	---

**Returns**

zero if the user clicked OK, non-zero if the user clicked Cancel.

Definition at line 716 of file pearl-scienta-preprocess.ipf.

**11.18.2.12 variable prompt\_Shockley\_anglefit ( string \* *param* )**

Definition at line 322 of file pearl-scienta-preprocess.ipf.

**11.18.2.13 threadsafe variable redim\_linbg\_reduction ( wave *source*, wave *dest1*, wave *dest2*, string \* *param* )**

linear background reduction function for incorrectly dimensioned scienta image

if the energy step size does not divide the energy range to an integer number, the scienta image is exported with the wrong array size. this can be fixed by redimensioning the array.

the current implementation works in the case where dimension 0 needs to be incremented. the function may be generalized to dimension 1 and/or decrementing by additional parameters. it is not known yet whether a generalization is needed or whether it can cover all cases.

background subtraction and peak integration is the same as by the [int\\_linbg\\_reduction\(\)](#) function.

#### Parameters

<i>source</i>	source wave Scienta detector image, energy axis along X, angle axis along Y
<i>dest1</i>	destination wave 1
<i>dest2</i>	destination wave 2 each wave is a one-dimensional intensity distribution the function may redimension these waves to one of the image dimensions (it must be clear to the user which dimension this is). the meaning of dest1 and dest2 is up to the particular function, e.g. dest1 could hold the mean value and dest2 the one-sigma error, or dest1 could hold the X-profile, and dest2 the Y-profile.
<i>param</i>	parameter string in a key1=value1;key2=value2;... list. the parameter string is passed by reference.

all region parameters are relative to the image size (0...1).

- Lcrop size of the lower cropping region
- Hcrop size of the upper cropping region
- Lsize size of the lower background integration region
- Hsize size of the upper background integration region
- Cpos center position of the of the peak integration region
- Csize size of the peak integration region

typical values (peak centered on detector, FWHM  $\sim$  20 % of image) Lcrop=0.11;Hcrop=0.11;Lsize=0.2;Hsize=0.2;Cpos=0.5;Csize=0.2

#### Returns

zero if successful, non-zero if an error occurs.

Definition at line 787 of file pearl-scienta-preprocess.ipf.

#### 11.18.2.14 threadsafe variable scienta\_ang\_transm ( wave *w*, variable *x* )

Definition at line 578 of file pearl-scienta-preprocess.ipf.

#### 11.18.2.15 variable scienta\_norm ( wave *w*, variable *x* )

Definition at line 544 of file pearl-scienta-preprocess.ipf.

#### 11.18.2.16 variable scienta\_poly\_bg ( wave *w*, variable *e*, variable *a* )

Definition at line 654 of file pearl-scienta-preprocess.ipf.

#### 11.18.2.17 threadsafe variable Shockley\_anglefit ( wave *source*, wave *dest1*, wave *dest2*, string \* *param* )

Definition at line 337 of file pearl-scienta-preprocess.ipf.

11.18.2.18 variable `test_gauss2_reduction ( wave image )`

apply the `gauss2_reduction` function to a single image

useful for testing or manual processing. to debug, (temporarily) remove the `threadsafe` attribute from the `gauss2_reduction` function.

Definition at line 810 of file `pearl-scienta-preprocess.ipf`.

11.18.2.19 variable `test_int_linbg ( wave image )`

Definition at line 180 of file `pearl-scienta-preprocess.ipf`.

11.18.2.20 variable `test_int_quadbg ( wave image )`

Definition at line 440 of file `pearl-scienta-preprocess.ipf`.

11.18.2.21 variable `test_shockley_anglefit ( wave image, variable branch )`

Definition at line 297 of file `pearl-scienta-preprocess.ipf`.

## 11.19 pearl-tools.ipf File Reference

```
#include "pearl-gui-tools"
```

### Functions

- variable `DefaultWaveIterator` (wave *w*, string *\*sdata*)
- variable `AppendToGraphIterator` (wave *w*, string *\*sdata*)
- variable `SumWavesIterator` (wave *w*, string *\*sdata*)
- string `IterateWaves` (string *matchStr*, funcref *iterator*, string *sdata*)
- variable `DefaultFolderIterator` (dfref *df*, string *\*sdata*)
- string `IterateDataFolders` (string *matchStr*, funcref *iterator*, string *sdata*, string *progress\_title*=defaultValue)

### 11.19.1 Function Documentation

11.19.1.1 variable `AppendToGraphIterator ( wave w, string * sdata )`

Definition at line 26 of file `pearl-tools.ipf`.

11.19.1.2 variable `DefaultFolderIterator ( dfref df, string * sdata )`

Definition at line 65 of file `pearl-tools.ipf`.

11.19.1.3 variable `DefaultWaveIterator ( wave w, string * sdata )`

Definition at line 18 of file `pearl-tools.ipf`.

11.19.1.4 `string IterateDataFolders ( string matchStr, funcref iterator, string sdata, string progress_title = defaultValue )`

Definition at line 79 of file `pearl-tools.ipf`.

11.19.1.5 `string IterateWaves ( string matchStr, funcref iterator, string sdata )`

Definition at line 45 of file `pearl-tools.ipf`.

11.19.1.6 `variable SumWavesIterator ( wave w, string * sdata )`

Definition at line 34 of file `pearl-tools.ipf`.

## 11.20 `pearl-vector-operations.ipf` File Reference

### Functions

- variable `rotate2d_x` (variable *xx*, variable *yy*, variable *angle*)
- variable `rotate2d_y` (variable *xx*, variable *yy*, variable *angle*)
- wave `create_rotation_matrix_free` ()
- wave `set_rotation_x` (wave matrix, variable *angle*)
- wave `set_rotation_y` (wave matrix, variable *angle*)
- wave `set_rotation_z` (wave matrix, variable *angle*)
- variable `rotate_x_wave` (wave *inout*, variable *angle*)
- variable `rotate_y_wave` (wave *inout*, variable *angle*)
- variable `rotate_z_wave` (wave *inout*, variable *angle*)

### 11.20.1 Function Documentation

11.20.1.1 `wave create_rotation_matrix_free ( )`

Definition at line 26 of file `pearl-vector-operations.ipf`.

11.20.1.2 `variable rotate2d_x ( variable xx, variable yy, variable angle )`

Definition at line 10 of file `pearl-vector-operations.ipf`.

11.20.1.3 `variable rotate2d_y ( variable xx, variable yy, variable angle )`

Definition at line 18 of file `pearl-vector-operations.ipf`.

11.20.1.4 `variable rotate_x_wave ( wave inout, variable angle )`

Definition at line 83 of file `pearl-vector-operations.ipf`.

11.20.1.5 `variable rotate_y_wave ( wave inout, variable angle )`

Definition at line 101 of file `pearl-vector-operations.ipf`.

11.20.1.6 variable rotate\_z\_wave ( wave *inout*, variable *angle* )

Definition at line 119 of file pearl-vector-operations.ipf.

11.20.1.7 wave set\_rotation\_x ( wave *matrix*, variable *angle* )

Definition at line 35 of file pearl-vector-operations.ipf.

11.20.1.8 wave set\_rotation\_y ( wave *matrix*, variable *angle* )

Definition at line 51 of file pearl-vector-operations.ipf.

11.20.1.9 wave set\_rotation\_z ( wave *matrix*, variable *angle* )

Definition at line 67 of file pearl-vector-operations.ipf.



# Index

ALREADY\_FILE\_OPEN

    errorCode, [23](#)

ARPES package, [15](#)

ad\_box\_filter

    pearl-area-display.ipf, [60](#)

ad\_brick\_slicer

    pearl-area-display.ipf, [61](#)

ad\_calc\_cursor\_profiles

    pearl-area-display.ipf, [61](#)

ad\_calc\_histogram

    pearl-area-display.ipf, [61](#)

ad\_calc\_profiles

    pearl-area-display.ipf, [61](#)

ad\_collect\_multiscan\_y

    pearl-area-profiles.ipf, [79](#)

ad\_default\_image\_filter

    pearl-area-display.ipf, [62](#)

ad\_display

    pearl-area-display.ipf, [62](#)

ad\_display\_brick

    pearl-area-display.ipf, [62](#)

ad\_display\_histogram

    pearl-area-display.ipf, [63](#)

ad\_display\_profiles

    pearl-area-display.ipf, [63](#)

ad\_display\_slice

    pearl-area-display.ipf, [63](#)

ad\_export\_profile

    pearl-area-display.ipf, [64](#)

ad\_extract\_rod

    pearl-area-profiles.ipf, [79](#)

ad\_extract\_rod\_x

    pearl-area-profiles.ipf, [80](#)

ad\_extract\_rod\_y

    pearl-area-profiles.ipf, [80](#)

ad\_extract\_rod\_z

    pearl-area-profiles.ipf, [80](#)

ad\_extract\_slab

    pearl-area-profiles.ipf, [80](#)

ad\_extract\_slab\_x

    pearl-area-profiles.ipf, [80](#)

ad\_extract\_slab\_y

    pearl-area-profiles.ipf, [81](#)

ad\_extract\_slab\_z

    pearl-area-profiles.ipf, [81](#)

ad\_gizmo\_set\_plane

    pearl-area-display.ipf, [64](#)

ad\_load\_dialog

    pearl-area-import.ipf, [71](#)

ad\_profile\_x

    pearl-area-profiles.ipf, [81](#)

ad\_profile\_x\_w

    pearl-area-profiles.ipf, [81](#)

ad\_profile\_y

    pearl-area-profiles.ipf, [81](#)

ad\_profile\_y\_w

    pearl-area-profiles.ipf, [81](#)

ad\_profiles\_crosshairs

    pearl-area-display.ipf, [65](#)

ad\_profiles\_cursor\_mode

    pearl-area-display.ipf, [65](#)

ad\_profiles\_hook

    pearl-area-display.ipf, [65](#)

ad\_profiles\_set\_cursor

    pearl-area-display.ipf, [65](#)

ad\_profiles\_set\_slice

    pearl-area-display.ipf, [66](#)

ad\_slicer\_init\_bg

    pearl-area-display.ipf, [66](#)

ad\_slicer\_move\_bg

    pearl-area-display.ipf, [66](#)

ad\_slicer\_start\_bg

    pearl-area-display.ipf, [66](#)

ad\_slicer\_stop\_bg

    pearl-area-display.ipf, [67](#)

ad\_suggest\_foldername

    pearl-area-import.ipf, [71](#)

ad\_transpose\_filter

    pearl-area-display.ipf, [67](#)

ad\_update\_profiles

    pearl-area-display.ipf, [67](#)

add\_image\_data

    pearl-anglescan-tracker.ipf, [52](#)

adh5\_default\_reduction

    pearl-area-import.ipf, [71](#)

adh5\_list\_reduction\_funcs

    pearl-area-import.ipf, [72](#)

adh5\_load\_complete

    pearl-area-import.ipf, [72](#)

adh5\_load\_detector

    pearl-area-import.ipf, [72](#)

adh5\_load\_detector\_image

    pearl-area-import.ipf, [73](#)

adh5\_load\_detector\_slabs

    pearl-area-import.ipf, [73](#)

adh5\_load\_info

    pearl-area-import.ipf, [73](#)

adh5\_load\_preview

- pearl-area-import.ipf, 74
- adh5\_load\_reduced
  - pearl-area-import.ipf, 74
- adh5\_load\_reduced\_detector
  - pearl-area-import.ipf, 75
- adh5\_loadattr\_all
  - pearl-area-import.ipf, 75
- adh5\_redim
  - pearl-area-import.ipf, 76
- adh5\_scale
  - pearl-area-import.ipf, 76
- adh5\_scale\_scan
  - pearl-area-import.ipf, 76
- adh5\_scale\_scienta
  - pearl-area-import.ipf, 76
- adh5\_setup\_profile
  - pearl-area-import.ipf, 76
- adh5\_test\_reduction\_func
  - pearl-area-import.ipf, 77
- AfterCompiledHook
  - pearl-anglescan-tracker.ipf, 52
  - pearl-arpes.ipf, 82
- AfterFileOpenHook
  - pearl-elog.ipf, 94
  - pearl-matrix-import.ipf, 109
- analyse\_curved\_edge
  - fermi-edge-analysis.ipf, 26
- analyser\_energy\_resolution
  - fermi-edge-analysis.ipf, 26
- AppendToGraphIterator
  - pearl-tools.ipf, 141
- ast\_add\_image
  - pearl-anglescan-tracker.ipf, 52
- ast\_callback\_data
  - pearl-anglescan-tracker.ipf, 52
- ast\_callback\_detector
  - pearl-anglescan-tracker.ipf, 52
- ast\_callback\_manip
  - pearl-anglescan-tracker.ipf, 53
- ast\_close
  - pearl-anglescan-tracker.ipf, 53
- ast\_export
  - pearl-anglescan-tracker.ipf, 53
- ast\_import
  - pearl-anglescan-tracker.ipf, 53
- ast\_prepare
  - pearl-anglescan-tracker.ipf, 53
- ast\_set\_processing
  - pearl-anglescan-tracker.ipf, 54
- ast\_setup
  - pearl-anglescan-tracker.ipf, 54
- ast\_update\_detector
  - pearl-anglescan-tracker.ipf, 54
- ast\_window\_hook
  - pearl-anglescan-tracker.ipf, 54
- attributes\_notebook
  - pearl-data-explorer.ipf, 85
- Au4f
  - pearl-fitfuncs.ipf, 106
- Au4f\_2p2
  - pearl-fitfuncs.ipf, 106
- Au4f\_2p3
  - pearl-fitfuncs.ipf, 106
- BeforeFileOpenHook
  - pearl-area-import.ipf, 77
  - pearl-matrix-import.ipf, 109
- bp\_attach
  - pearl-elog.ipf, 94
- bp\_attach\_allnone
  - pearl-elog.ipf, 94
- bp\_attach\_top
  - pearl-elog.ipf, 94
- bp\_attach\_updown
  - pearl-elog.ipf, 94
- bp\_attr\_notebook
  - pearl-data-explorer.ipf, 85
- bp\_browse\_filepath
  - pearl-data-explorer.ipf, 85
- bp\_capture
  - pearl-anglescan-tracker.ipf, 54
- bp\_clear
  - pearl-elog.ipf, 94
- bp\_dataset\_display
  - pearl-data-explorer.ipf, 85
- bp\_dataset\_folder
  - pearl-data-explorer.ipf, 85
- bp\_dataset\_next
  - pearl-data-explorer.ipf, 85
- bp\_dataset\_prev
  - pearl-data-explorer.ipf, 85
- bp\_extract\_slice
  - pearl-area-display.ipf, 67
- bp\_file\_next
  - pearl-data-explorer.ipf, 85
- bp\_file\_prev
  - pearl-data-explorer.ipf, 85
- bp\_load\_files
  - pearl-data-explorer.ipf, 85
- bp\_load\_files\_opt
  - pearl-data-explorer.ipf, 85
- bp\_load\_prefs
  - pearl-data-explorer.ipf, 85
- bp\_login
  - pearl-elog.ipf, 95
- bp\_logout
  - pearl-elog.ipf, 95
- bp\_move\_slice
  - pearl-area-display.ipf, 67
- bp\_move\_slice\_center
  - pearl-area-display.ipf, 68
- bp\_reset\_cursors
  - pearl-area-display.ipf, 68
- bp\_save\_graphs
  - pearl-elog.ipf, 95
- bp\_save\_prefs
  - pearl-data-explorer.ipf, 85

- bp\_submit
  - pearl-elog.ipf, 95
- bp\_update\_datasets
  - pearl-data-explorer.ipf, 86
- bp\_update\_filelist
  - pearl-data-explorer.ipf, 86
- broadening
  - DoniachSunjicStruct, 21
- Calc\_DoniachSunjicBroad
  - pearl-fitfuncs.ipf, 106
- Calc\_The\_step
  - pearl-anglescan-process.ipf, 33
- calc\_graph\_azi
  - pearl-anglescan-process.ipf, 31
- calc\_graph\_polar
  - pearl-anglescan-process.ipf, 31
- calc\_graph\_radius
  - pearl-anglescan-process.ipf, 32
- calc\_nth
  - pearl-anglescan-process.ipf, 32
- calc\_phi\_step
  - pearl-anglescan-process.ipf, 33
- calc\_y\_profile\_mins
  - pearl-area-profiles.ipf, 81
- CalcN\_Theta
  - pearl-anglescan-process.ipf, 33
- capture\_int\_linbg\_cursors
  - pearl-scienta-preprocess.ipf, 138
- cart2polar
  - pearl-polar-coordinates.ipf, 118
- cart2polar\_wave
  - pearl-polar-coordinates.ipf, 118
- check\_contrast
  - pearl-anglescan-process.ipf, 33
- check\_package\_folder
  - pearl-matrix-import.ipf, 109
- cleanup\_temp\_files
  - pearl-elog.ipf, 95
- clear\_hemi\_grid
  - pearl-anglescan-process.ipf, 33
- convert\_angles\_ttpa2polar
  - pearl-anglescan-process.ipf, 33
- convert\_angles\_ttpd2polar
  - pearl-anglescan-process.ipf, 34
- convolution
  - DoniachSunjicStruct, 21
- create\_cmd\_file
  - pearl-elog.ipf, 95
- create\_graph\_file
  - pearl-elog.ipf, 95
- create\_message\_file
  - pearl-elog.ipf, 96
- create\_rotation\_matrix\_free
  - pearl-vector-operations.ipf, 142
- crop\_strip
  - pearl-anglescan-process.ipf, 34
- csr\_int\_linbg\_reduction
  - pearl-scienta-preprocess.ipf, 138
- DefaultFolderIterator
  - pearl-tools.ipf, 141
- DefaultWavelterator
  - pearl-tools.ipf, 141
- Display2dProfiles
  - pearl-menu.ipf, 115
- Display3dSlicer
  - pearl-menu.ipf, 115
- display\_dataset
  - pearl-data-explorer.ipf, 86
- display\_hemi\_scan
  - pearl-anglescan-process.ipf, 34
- display\_polar\_graph
  - pearl-anglescan-process.ipf, 35
- display\_preview\_trace
  - pearl-data-explorer.ipf, 86
- display\_progress\_panel
  - pearl-gui-tools.ipf, 107
- display\_scanlines
  - pearl-anglescan-process.ipf, 36
- DisplayGizmoSlicer
  - pearl-menu.ipf, 115
- DoniachSunjicBroad
  - pearl-fitfuncs.ipf, 106
- DoniachSunjicBroadS
  - pearl-fitfuncs.ipf, 106
- DoniachSunjicStruct, 21
  - broadening, 21
  - convolution, 21
  - model, 21
  - oversampling, 21
  - precision, 21
  - pw, 22
  - xw, 22
  - yw, 22
- draw\_diffraction\_cone
  - pearl-anglescan-process.ipf, 36
- draw\_hemi\_axes
  - pearl-anglescan-process.ipf, 37
- ds1\_bg
  - pearl-fitfuncs.ipf, 106
- ds2\_bg
  - pearl-fitfuncs.ipf, 106
- ds4\_bg
  - pearl-fitfuncs.ipf, 106
- ds6\_bg
  - pearl-fitfuncs.ipf, 106
- duplicate\_hemi\_scan
  - pearl-anglescan-process.ipf, 37
- EMPTY\_RESULTFILE
  - errorCode, 23
- edit\_offsets
  - pearl-anglescan-tracker.ipf, 55
- edit\_reduction\_params
  - pearl-anglescan-tracker.ipf, 55
- elog\_add\_attachment
  - pearl-elog.ipf, 96

- elog\_config
  - pearl-elog.ipf, 97
- elog\_create\_entry
  - pearl-elog.ipf, 97
- elog\_create\_logbook
  - pearl-elog.ipf, 97
- elog\_init\_pearl\_templates
  - pearl-elog.ipf, 97
- elog\_login
  - pearl-elog.ipf, 98
- elog\_logout
  - pearl-elog.ipf, 98
- elog\_panel\_hook
  - pearl-elog.ipf, 98
- elog\_parse\_id
  - pearl-elog.ipf, 104
- elog\_prompt\_logbook
  - pearl-elog.ipf, 98
- elog\_prompt\_login
  - pearl-elog.ipf, 98
- elog\_success\_msg
  - pearl-elog.ipf, 104
- elog\_validate\_attributes
  - pearl-elog.ipf, 99
- epics\_connect
  - pearl-anglescan-tracker.ipf, 55
- epics\_disconnect
  - pearl-anglescan-tracker.ipf, 55
- epics\_disconnect\_chid
  - pearl-anglescan-tracker.ipf, 55
- errorCode, 22
  - ALREADY\_FILE\_OPEN, 23
  - EMPTY\_RESULTFILE, 23
  - FILE\_NOT\_READABLE, 23
  - INTERNAL\_ERROR\_CONVERTING\_DATA, 23
  - INVALID\_RANGE, 23
  - NO\_FILE\_OPEN, 23
  - NO\_NEW\_BRICKLETS, 23
  - SUCCESS, 23
  - UNKNOWN\_ERROR, 23
  - WAVE\_EXIST, 23
  - WRONG\_PARAMETER, 23
- export\_tracker\_data
  - pearl-anglescan-tracker.ipf, 55
- extend\_data
  - pearl-anglescan-tracker.ipf, 55
- extract\_preview\_image
  - pearl-data-explorer.ipf, 86
- FILE\_NOT\_READABLE
  - errorCode, 23
- fermi-edge-analysis.ipf, 25
  - analyse\_curved\_edge, 26
  - analyser\_energy\_resolution, 26
  - FermiFuncLinDOS2D\_corr, 26
  - hemi\_radius\_mm, 27
  - integrate\_curved\_edge, 26
  - mcp\_radius\_epass, 27
  - mcp\_radius\_mm, 27
  - mcp\_radius\_pix, 27
  - record\_results, 26
  - show\_shift, 26
  - slit\_correction, 27
  - slit\_shift, 27
- FermiFuncLinDOS2D\_corr
  - fermi-edge-analysis.ipf, 26
- FermiGaussConv
  - pearl-fitfuncs.ipf, 106
- find\_attr\_folder
  - pearl-pshell-import.ipf, 121
- find\_hemi\_data
  - pearl-anglescan-process.ipf, 38
- find\_scale\_wave
  - pearl-pshell-import.ipf, 121
- find\_scan\_folder
  - pearl-pshell-import.ipf, 121
- Fit\_DoniachSunjicBroad
  - pearl-fitfuncs.ipf, 107
- fit\_scienta\_ang\_transm
  - pearl-scienta-preprocess.ipf, 138
- fit\_scienta\_poly\_bg
  - pearl-scienta-preprocess.ipf, 138
- format\_url
  - pearl-elog.ipf, 99
- gather\_batch
  - pearl-otf-import.ipf, 116
- gauss2\_reduction
  - pearl-scienta-preprocess.ipf, 138
- get\_default\_panel\_name
  - pearl-elog.ipf, 99
- get\_elog\_df
  - pearl-elog.ipf, 99
- get\_hemi\_nickname
  - pearl-anglescan-process.ipf, 38
- get\_hemi\_prefix
  - pearl-anglescan-process.ipf, 39
- get\_log\_path
  - pearl-elog.ipf, 99
- get\_panel\_attributes
  - pearl-elog.ipf, 100
- get\_panel\_graphs
  - pearl-elog.ipf, 100
- get\_panel\_message
  - pearl-elog.ipf, 100
- get\_source\_image
  - pearl-area-display.ipf, 68
- get\_timestamp
  - pearl-elog.ipf, 100
- get\_view\_folder
  - pearl-area-display.ipf, 68
- get\_view\_image
  - pearl-area-display.ipf, 68
- GetAttrDataFolderDFR
  - pearl-area-import.ipf, 77
- graphname\_from\_dfref
  - pearl-area-display.ipf, 68

- hemi\_add\_anglescan
  - pearl-anglescan-process.ipf, [39](#)
- hemi\_add\_aziscan
  - pearl-anglescan-process.ipf, [39](#)
- hemi\_azi\_cut
  - pearl-anglescan-process.ipf, [39](#)
- hemi\_polar\_cut
  - pearl-anglescan-process.ipf, [40](#)
- hemi\_radius\_mm
  - fermi-edge-analysis.ipf, [27](#)
- INTERNAL\_ERROR\_CONVERTING\_DATA
  - errorCode, [23](#)
- INVALID\_RANGE
  - errorCode, [23](#)
- IgorBeforeNewHook
  - pearl-elog.ipf, [100](#)
- IgorQuitHook
  - pearl-anglescan-tracker.ipf, [55](#)
  - pearl-elog.ipf, [101](#)
- import\_tpi\_scan
  - pearl-anglescan-process.ipf, [40](#)
- import\_tracker\_data
  - pearl-anglescan-tracker.ipf, [56](#)
- init\_package
  - pearl-anglescan-tracker.ipf, [56](#)
  - pearl-data-explorer.ipf, [86](#)
  - pearl-elog.ipf, [101](#)
  - pearl-matrix-import.ipf, [110](#)
- init\_volatile\_vars
  - pearl-elog.ipf, [101](#)
- initStruct
  - pearl-matrix-import.ipf, [110](#)
- int\_linbg\_reduction
  - pearl-scienta-preprocess.ipf, [139](#)
- int\_quadbg\_reduction
  - pearl-scienta-preprocess.ipf, [139](#)
- integrate\_curved\_edge
  - fermi-edge-analysis.ipf, [26](#)
- interpolate\_hemi\_scan
  - pearl-anglescan-process.ipf, [41](#)
- IterateDataFolders
  - pearl-tools.ipf, [141](#)
- IterateWaves
  - pearl-tools.ipf, [142](#)
- itx\_suggest\_foldername
  - pearl-data-explorer.ipf, [86](#)
- kAngleDimLabel
  - pearl-pshell-import.ipf, [136](#)
- kAttachColName
  - pearl-elog.ipf, [104](#)
- kAttachColSel
  - pearl-elog.ipf, [104](#)
- kAttachColTitle
  - pearl-elog.ipf, [104](#)
- kDataDimLabel
  - pearl-pshell-import.ipf, [136](#)
- kDetectorSensitivity
  - pearl-pshell-import.ipf, [136](#)
- kEnergyDimLabel
  - pearl-pshell-import.ipf, [136](#)
- kPreviewDatasets
  - pearl-pshell-import.ipf, [136](#)
- kProjArea
  - pearl-anglescan-process.ipf, [49](#)
- kProjDist
  - pearl-anglescan-process.ipf, [49](#)
- kProjGnom
  - pearl-anglescan-process.ipf, [49](#)
- kProjOrtho
  - pearl-anglescan-process.ipf, [49](#)
- kProjScaleArea
  - pearl-anglescan-process.ipf, [49](#)
- kProjScaleDist
  - pearl-anglescan-process.ipf, [49](#)
- kProjScaleGnom
  - pearl-anglescan-process.ipf, [49](#)
- kProjScaleOrtho
  - pearl-anglescan-process.ipf, [50](#)
- kProjScaleStereo
  - pearl-anglescan-process.ipf, [50](#)
- kProjStereo
  - pearl-anglescan-process.ipf, [50](#)
- kScanDimLabel
  - pearl-pshell-import.ipf, [136](#)
- kScientaScalingDatasets
  - pearl-pshell-import.ipf, [136](#)
- kTransposedDatasets
  - pearl-pshell-import.ipf, [136](#)
- kdfPersistent
  - pearl-elog.ipf, [104](#)
- kdfRoot
  - pearl-elog.ipf, [104](#)
- kdfTemplates
  - pearl-elog.ipf, [105](#)
- kdfVolatile
  - pearl-elog.ipf, [105](#)
- kill\_progress\_panel
  - pearl-gui-tools.ipf, [107](#)
- ks\_filematch\_adh5
  - pearl-data-explorer.ipf, [90](#)
- ks\_filematch\_itx
  - pearl-data-explorer.ipf, [90](#)
- ks\_filematch\_mtrx
  - pearl-data-explorer.ipf, [90](#)
  - pearl-matrix-import.ipf, [115](#)
- ks\_filematch\_pshell
  - pearl-data-explorer.ipf, [90](#)
- lbp\_datasets
  - pearl-data-explorer.ipf, [86](#)
- lbp\_filelist
  - pearl-data-explorer.ipf, [86](#)
- line\_average
  - pearl-anglescan-process.ipf, [41](#)
- list\_logbooks
  - pearl-elog.ipf, [101](#)

- load\_file
  - pearl-data-explorer.ipf, [86](#)
- load\_hdf\_file
  - pearl-data-explorer.ipf, [86](#)
- load\_hemi\_scan
  - pearl-anglescan-process.ipf, [41](#)
- load\_itx\_file
  - pearl-data-explorer.ipf, [87](#)
- load\_mtrx\_file
  - pearl-data-explorer.ipf, [87](#)
- load\_prefs
  - pearl-anglescan-tracker.ipf, [56](#)
  - pearl-data-explorer.ipf, [87](#)
  - pearl-elog.ipf, [101](#)
- load\_pshell\_file
  - pearl-data-explorer.ipf, [87](#)
- load\_selected\_files
  - pearl-data-explorer.ipf, [87](#)
- load\_tracker\_data
  - pearl-anglescan-tracker.ipf, [56](#)
- LoadPearlArpes
  - pearl-menu.ipf, [116](#)
- LoadPearlOptics
  - pearl-menu.ipf, [116](#)
- LoadPearlPreparation
  - pearl-menu.ipf, [116](#)
- mainpage.dox, [27](#)
- make\_hemi\_grid
  - pearl-anglescan-process.ipf, [41](#)
- make\_view\_folder
  - pearl-area-display.ipf, [69](#)
- matrix\_format\_elog\_message
  - pearl-matrix-import.ipf, [110](#)
- matrix\_preview\_2d
  - pearl-matrix-import.ipf, [110](#)
- mcp\_radius\_epass
  - fermi-edge-analysis.ipf, [27](#)
- mcp\_radius\_mm
  - fermi-edge-analysis.ipf, [27](#)
- mcp\_radius\_pix
  - fermi-edge-analysis.ipf, [27](#)
- model
  - DoniachSunjicStruct, [21](#)
- move\_attach\_item
  - pearl-elog.ipf, [102](#)
- mtrx\_create\_folder
  - pearl-matrix-import.ipf, [110](#)
- mtrx\_file\_brickletID
  - pearl-matrix-import.ipf, [110](#)
- mtrx\_get\_cycle\_folder
  - pearl-matrix-import.ipf, [111](#)
- mtrx\_load\_all
  - pearl-matrix-import.ipf, [111](#)
- mtrx\_load\_file
  - pearl-matrix-import.ipf, [111](#)
- mtrx\_load\_info
  - pearl-matrix-import.ipf, [112](#)
- mtrx\_load\_preview
  - pearl-matrix-import.ipf, [112](#)
- mtrx\_open\_file
  - pearl-matrix-import.ipf, [113](#)
- mtrx\_parse\_filename
  - pearl-matrix-import.ipf, [113](#)
- mtrx\_scale\_dataset
  - pearl-matrix-import.ipf, [114](#)
- mtrx\_split\_filename
  - pearl-matrix-import.ipf, [114](#)
- NO\_FILE\_OPEN
  - errorCode, [23](#)
- NO\_NEW\_BRICKLETS
  - errorCode, [23](#)
- normalize\_strip\_2d
  - pearl-anglescan-process.ipf, [42](#)
- normalize\_strip\_phi
  - pearl-anglescan-process.ipf, [42](#)
- normalize\_strip\_theta
  - pearl-anglescan-process.ipf, [43](#)
- normalize\_strip\_x
  - pearl-anglescan-process.ipf, [43](#)
- notebook\_add\_attributes
  - pearl-data-explorer.ipf, [87](#)
- otf\_gather\_batch
  - pearl-otf-import.ipf, [116](#)
- otf\_gather\_iterator
  - pearl-otf-import.ipf, [117](#)
- otf\_interp
  - pearl-otf-import.ipf, [117](#)
- otf\_load\_itx
  - pearl-otf-import.ipf, [117](#)
- otf\_load\_itx\_all
  - pearl-otf-import.ipf, [117](#)
- otf\_load\_itx\_match
  - pearl-otf-import.ipf, [117](#)
- otf\_rename\_folders
  - pearl-otf-import.ipf, [117](#)
- otf\_rename\_folders\_iterator
  - pearl-otf-import.ipf, [117](#)
- otf\_smo\_int
  - pearl-otf-import.ipf, [117](#)
- oversampling
  - DoniachSunjicStruct, [21](#)
- package\_name
  - pearl-data-explorer.ipf, [90](#)
  - pearl-elog.ipf, [105](#)
  - pearl-matrix-import.ipf, [115](#)
- package\_path
  - pearl-anglescan-tracker.ipf, [58](#)
  - pearl-data-explorer.ipf, [90](#)
  - pearl-elog.ipf, [105](#)
  - pearl-matrix-import.ipf, [115](#)
- parse\_result
  - pearl-elog.ipf, [102](#)
- pearl-anglescan-process.ipf, [27](#)
- Calc\_The\_step, [33](#)

- calc\_graph\_azi, 31
- calc\_graph\_polar, 31
- calc\_graph\_radius, 32
- calc\_nth, 32
- calc\_phi\_step, 33
- CalcN\_Theta, 33
- check\_contrast, 33
- clear\_hemi\_grid, 33
- convert\_angles\_ttpa2polar, 33
- convert\_angles\_ttpd2polar, 34
- crop\_strip, 34
- display\_hemi\_scan, 34
- display\_polar\_graph, 35
- display\_scanlines, 36
- draw\_diffraction\_cone, 36
- draw\_hemi\_axes, 37
- duplicate\_hemi\_scan, 37
- find\_hemi\_data, 38
- get\_hemi\_nickname, 38
- get\_hemi\_prefix, 39
- hemi\_add\_anglescan, 39
- hemi\_add\_aziscan, 39
- hemi\_azi\_cut, 39
- hemi\_polar\_cut, 40
- import\_tpi\_scan, 40
- interpolate\_hemi\_scan, 41
- kProjArea, 49
- kProjDist, 49
- kProjGnom, 49
- kProjOrtho, 49
- kProjScaleArea, 49
- kProjScaleDist, 49
- kProjScaleGnom, 49
- kProjScaleOrtho, 50
- kProjScaleStereo, 50
- kProjStereo, 50
- line\_average, 41
- load\_hemi\_scan, 41
- make\_hemi\_grid, 41
- normalize\_strip\_2d, 42
- normalize\_strip\_phi, 42
- normalize\_strip\_theta, 43
- normalize\_strip\_x, 43
- pizza\_service, 44
- pizza\_service\_2, 45
- polar\_graph\_hook, 46
- quick\_pizza\_image, 46
- rotate\_hemi\_scan, 46
- save\_hemi\_scan, 47
- set\_contrast, 47
- set\_polar\_graph\_cursor, 47
- show\_analyser\_line, 47
- strip\_delete\_frames, 48
- trim\_hemi\_scan, 48
- update\_polar\_info, 49
- pearl-anglescan-tracker.ipf, 50
  - add\_image\_data, 52
  - AfterCompiledHook, 52
  - ast\_add\_image, 52
  - ast\_callback\_data, 52
  - ast\_callback\_detector, 52
  - ast\_callback\_manip, 53
  - ast\_close, 53
  - ast\_export, 53
  - ast\_import, 53
  - ast\_prepare, 53
  - ast\_set\_processing, 54
  - ast\_setup, 54
  - ast\_update\_detector, 54
  - ast\_window\_hook, 54
  - bp\_capture, 54
  - edit\_offsets, 55
  - edit\_reduction\_params, 55
  - epics\_connect, 55
  - epics\_disconnect, 55
  - epics\_disconnect\_chid, 55
  - export\_tracker\_data, 55
  - extend\_data, 55
  - IgorQuitHook, 55
  - import\_tracker\_data, 56
  - init\_package, 56
  - load\_prefs, 56
  - load\_tracker\_data, 56
  - package\_path, 58
  - pmp\_data, 56
  - pmp\_data\_mouseup, 56
  - pmp\_parameters, 56
  - pmp\_parameters\_mouseup, 56
  - prefs\_objects, 58
  - process\_image\_data, 56
  - save\_prefs, 56
  - save\_tracker\_data, 57
  - setup\_data, 57
  - setup\_detector, 57
  - setup\_graph, 57
  - toggle\_capture, 57
  - update\_capture, 57
  - update\_data\_graph, 57
  - update\_detector, 57
  - update\_detector\_graph, 58
  - version, 58
- pearl-area-display.ipf, 58
  - ad\_box\_filter, 60
  - ad\_brick\_slicer, 61
  - ad\_calc\_cursor\_profiles, 61
  - ad\_calc\_histogram, 61
  - ad\_calc\_profiles, 61
  - ad\_default\_image\_filter, 62
  - ad\_display, 62
  - ad\_display\_brick, 62
  - ad\_display\_histogram, 63
  - ad\_display\_profiles, 63
  - ad\_display\_slice, 63
  - ad\_export\_profile, 64
  - ad\_gizmo\_set\_plane, 64
  - ad\_profiles\_crosshairs, 65

- ad\_profiles\_cursor\_mode, 65
- ad\_profiles\_hook, 65
- ad\_profiles\_set\_cursor, 65
- ad\_profiles\_set\_slice, 66
- ad\_slicer\_init\_bg, 66
- ad\_slicer\_move\_bg, 66
- ad\_slicer\_start\_bg, 66
- ad\_slicer\_stop\_bg, 67
- ad\_transpose\_filter, 67
- ad\_update\_profiles, 67
- bp\_extract\_slice, 67
- bp\_move\_slice, 67
- bp\_move\_slice\_center, 68
- bp\_reset\_cursors, 68
- get\_source\_image, 68
- get\_view\_folder, 68
- get\_view\_image, 68
- graphname\_from\_dfref, 68
- make\_view\_folder, 69
- pmp\_export, 69
- set\_trace\_colors, 69
- slp\_slice\_position, 69
- svp\_slice\_position, 69
- svp\_smoothing, 69
- update\_slice\_info, 69
- pearl-area-import.ipf, 69
  - ad\_load\_dialog, 71
  - ad\_suggest\_foldername, 71
  - adh5\_default\_reduction, 71
  - adh5\_list\_reduction\_funcs, 72
  - adh5\_load\_complete, 72
  - adh5\_load\_detector, 72
  - adh5\_load\_detector\_image, 73
  - adh5\_load\_detector\_slabs, 73
  - adh5\_load\_info, 73
  - adh5\_load\_preview, 74
  - adh5\_load\_reduced, 74
  - adh5\_load\_reduced\_detector, 75
  - adh5\_loadattr\_all, 75
  - adh5\_redim, 76
  - adh5\_scale, 76
  - adh5\_scale\_scan, 76
  - adh5\_scale\_scienta, 76
  - adh5\_setup\_profile, 76
  - adh5\_test\_reduction\_func, 77
  - BeforeFileOpenHook, 77
  - GetAttrDataFolderDFR, 77
  - read\_attribute\_info, 77
  - reduce\_slab\_image, 77
  - reduce\_slab\_worker, 78
- pearl-area-profiles.ipf, 78
  - ad\_collect\_multiscan\_y, 79
  - ad\_extract\_rod, 79
  - ad\_extract\_rod\_x, 80
  - ad\_extract\_rod\_y, 80
  - ad\_extract\_rod\_z, 80
  - ad\_extract\_slab, 80
  - ad\_extract\_slab\_x, 80
  - ad\_extract\_slab\_y, 81
  - ad\_extract\_slab\_z, 81
  - ad\_profile\_x, 81
  - ad\_profile\_x\_w, 81
  - ad\_profile\_y, 81
  - ad\_profile\_y\_w, 81
  - calc\_y\_profile\_mins, 81
- pearl-arpes.ipf, 81
  - AfterCompiledHook, 82
  - UnloadPearlArpesPackage, 82
- pearl-data-explorer.ipf, 83
  - attributes\_notebook, 85
  - bp\_attr\_notebook, 85
  - bp\_browse\_filepath, 85
  - bp\_dataset\_display, 85
  - bp\_dataset\_folder, 85
  - bp\_dataset\_next, 85
  - bp\_dataset\_prev, 85
  - bp\_file\_next, 85
  - bp\_file\_prev, 85
  - bp\_load\_files, 85
  - bp\_load\_files\_opt, 85
  - bp\_load\_prefs, 85
  - bp\_save\_prefs, 85
  - bp\_update\_datasets, 86
  - bp\_update\_filelist, 86
  - display\_dataset, 86
  - display\_preview\_trace, 86
  - extract\_preview\_image, 86
  - init\_package, 86
  - itx\_suggest\_foldername, 86
  - ks\_filematch\_adh5, 90
  - ks\_filematch\_itx, 90
  - ks\_filematch\_mtrx, 90
  - ks\_filematch\_pshell, 90
  - lbp\_datasets, 86
  - lbp\_filelist, 86
  - load\_file, 86
  - load\_hdf\_file, 86
  - load\_itx\_file, 87
  - load\_mtrx\_file, 87
  - load\_prefs, 87
  - load\_pshell\_file, 87
  - load\_selected\_files, 87
  - notebook\_add\_attributes, 87
  - package\_name, 90
  - package\_path, 90
  - pearl\_data\_explorer, 87
  - pearl\_file\_type, 87
  - PearlDataExplorer, 87
  - preview\_attributes, 88
  - preview\_datafolder, 88
  - preview\_dataset, 88
  - preview\_file, 88
  - preview\_hdf\_file, 88
  - preview\_itx\_file, 88
  - preview\_mtrx\_file, 88
  - preview\_pshell\_file, 89

- preview\_setscale\_x, 89
- prompt\_default\_process, 89
- prompt\_func\_params, 89
- prompt\_hdf\_options, 89
- save\_prefs, 89
- show\_preview\_graph, 90
- test\_attributes\_notebook, 90
- update\_datasets, 90
- update\_filelist, 90
- pearl-elog.ipf, 91
  - AfterFileOpenHook, 94
  - bp\_attach, 94
  - bp\_attach\_allnone, 94
  - bp\_attach\_top, 94
  - bp\_attach\_updown, 94
  - bp\_clear, 94
  - bp\_login, 95
  - bp\_logout, 95
  - bp\_save\_graphs, 95
  - bp\_submit, 95
  - cleanup\_temp\_files, 95
  - create\_cmd\_file, 95
  - create\_graph\_file, 95
  - create\_message\_file, 96
  - elog\_add\_attachment, 96
  - elog\_config, 97
  - elog\_create\_entry, 97
  - elog\_create\_logbook, 97
  - elog\_init\_pearl\_templates, 97
  - elog\_login, 98
  - elog\_logout, 98
  - elog\_panel\_hook, 98
  - elog\_parse\_id, 104
  - elog\_prompt\_logbook, 98
  - elog\_prompt\_login, 98
  - elog\_success\_msg, 104
  - elog\_validate\_attributes, 99
  - format\_url, 99
  - get\_default\_panel\_name, 99
  - get\_elog\_df, 99
  - get\_log\_path, 99
  - get\_panel\_attributes, 100
  - get\_panel\_graphs, 100
  - get\_panel\_message, 100
  - get\_timestamp, 100
  - IgorBeforeNewHook, 100
  - IgorQuitHook, 101
  - init\_package, 101
  - init\_volatile\_vars, 101
  - kAttachColName, 104
  - kAttachColSel, 104
  - kAttachColTitle, 104
  - kdfPersistent, 104
  - kdfRoot, 104
  - kdfTemplates, 105
  - kdfVolatile, 105
  - list\_logbooks, 101
  - load\_prefs, 101
  - move\_attach\_item, 102
  - package\_name, 105
  - package\_path, 105
  - parse\_result, 102
  - pearl\_elog, 102
  - PearlElogPanel, 102
  - prepare\_command\_line, 102
  - prepare\_graph\_attachments, 103
  - save\_prefs, 103
  - set\_panel\_attributes, 103
  - set\_panel\_graphs, 103
  - set\_panel\_message, 104
  - update\_attach\_items, 104
- pearl-fitfuncs.ipf, 105
  - Au4f, 106
  - Au4f\_2p2, 106
  - Au4f\_2p3, 106
  - Calc\_DoniachSunjicBroad, 106
  - DoniachSunjicBroad, 106
  - DoniachSunjicBroadS, 106
  - ds1\_bg, 106
  - ds2\_bg, 106
  - ds4\_bg, 106
  - ds6\_bg, 106
  - FermiGaussConv, 106
  - Fit\_DoniachSunjicBroad, 107
  - ShowComponents\_Au4f\_2p2, 107
  - ShowComponents\_Au4f\_2p3, 107
  - version, 107
- pearl-gui-tools.ipf, 107
  - display\_progress\_panel, 107
  - kill\_progress\_panel, 107
  - update\_progress\_panel, 107
- pearl-matrix-import.ipf, 108
  - AfterFileOpenHook, 109
  - BeforeFileOpenHook, 109
  - check\_package\_folder, 109
  - init\_package, 110
  - initStruct, 110
  - ks\_filematch\_mtrx, 115
  - matrix\_format\_elog\_message, 110
  - matrix\_preview\_2d, 110
  - mtrx\_create\_folder, 110
  - mtrx\_file\_brickletID, 110
  - mtrx\_get\_cycle\_folder, 111
  - mtrx\_load\_all, 111
  - mtrx\_load\_file, 111
  - mtrx\_load\_info, 112
  - mtrx\_load\_preview, 112
  - mtrx\_open\_file, 113
  - mtrx\_parse\_filename, 113
  - mtrx\_scale\_dataset, 114
  - mtrx\_split\_filename, 114
  - package\_name, 115
  - package\_path, 115
  - preview\_matrix\_file, 114
  - subtract\_line\_bg, 115
- pearl-menu.ipf, 115

- Display2dProfiles, 115
- Display3dSlicer, 115
- DisplayGizmoSlicer, 115
- LoadPearlArpes, 116
- LoadPearlOptics, 116
- LoadPearlPreparation, 116
- PearlAnglescanTracker, 116
- PearlLiveDisplay, 116
- PearlMenuEnableFunc, 116
- pearl-otf-import.ipf, 116
  - gather\_batch, 116
  - otf\_gather\_batch, 116
  - otf\_gather\_iterator, 117
  - otf\_interp, 117
  - otf\_load\_itx, 117
  - otf\_load\_itx\_all, 117
  - otf\_load\_itx\_match, 117
  - otf\_rename\_folders, 117
  - otf\_rename\_folders\_iterator, 117
  - otf\_smo\_int, 117
- pearl-polar-coordinates.ipf, 117
  - cart2polar, 118
  - cart2polar\_wave, 118
  - polar2cart, 118
  - polar2cart\_wave, 118
  - polar\_distance, 118
- pearl-pshell-import.ipf, 118
  - find\_attr\_folder, 121
  - find\_scale\_wave, 121
  - find\_scan\_folder, 121
  - kAngleDimLabel, 136
  - kDataDimLabel, 136
  - kDetectorSensitivity, 136
  - kEnergyDimLabel, 136
  - kPreviewDatasets, 136
  - kScanDimLabel, 136
  - kScientaScalingDatasets, 136
  - kTransposedDatasets, 136
  - ps\_detect\_scale, 121
  - ps\_scale\_dataset, 122
  - ps\_scale\_dataset\_2, 122
  - ps\_scale\_datasets, 123
  - ps\_set\_dimlabels, 123
  - ps\_set\_dimlabels2, 124
  - psh5\_close\_file, 124
  - psh5\_list\_scan\_datasets, 124
  - psh5\_list\_scan\_regions, 125
  - psh5\_list\_scans, 125
  - psh5\_load\_complete, 125
  - psh5\_load\_dataset, 126
  - psh5\_load\_dataset\_meta, 127
  - psh5\_load\_dataset\_reduced, 127
  - psh5\_load\_dataset\_slab, 128
  - psh5\_load\_dataset\_slabs, 128
  - psh5\_load\_info, 129
  - psh5\_load\_preview, 129
  - psh5\_load\_reduced, 130
  - psh5\_load\_scan\_attrs, 131
  - psh5\_load\_scan\_complete, 131
  - psh5\_load\_scan\_data, 132
  - psh5\_load\_scan\_info, 132
  - psh5\_load\_scan\_meta, 133
  - psh5\_load\_scan\_preview, 133
  - psh5\_load\_scan\_section, 134
  - psh5\_open\_file, 134
  - reduce\_slab\_image, 135
  - reduce\_slab\_worker, 135
  - select\_dataset, 135
  - twave2list, 135
  - wave2list, 135
- pearl-scienta-preprocess.ipf, 137
  - capture\_int\_linbg\_cursors, 138
  - csr\_int\_linbg\_reduction, 138
  - fit\_scienta\_ang\_transm, 138
  - fit\_scienta\_poly\_bg, 138
  - gauss2\_reduction, 138
  - int\_linbg\_reduction, 139
  - int\_quadbg\_reduction, 139
  - prompt\_Shockley\_anglefit, 139
  - prompt\_gauss2\_reduction, 139
  - prompt\_int\_linbg\_reduction, 139
  - prompt\_int\_quadbg\_reduction, 139
  - prompt\_redim\_linbg\_reduction, 139
  - redim\_linbg\_reduction, 139
  - scienta\_ang\_transm, 140
  - scienta\_norm, 140
  - scienta\_poly\_bg, 140
  - Shockley\_anglefit, 140
  - test\_gauss2\_reduction, 140
  - test\_int\_linbg, 141
  - test\_int\_quadbg, 141
  - test\_shockley\_anglefit, 141
- pearl-tools.ipf, 141
  - AppendToGraphIterator, 141
  - DefaultFolderIterator, 141
  - DefaultWaveIterator, 141
  - IterateDataFolders, 141
  - IterateWaves, 142
  - SumWavesIterator, 142
- pearl-vector-operations.ipf, 142
  - create\_rotation\_matrix\_free, 142
  - rotate2d\_x, 142
  - rotate2d\_y, 142
  - rotate\_x\_wave, 142
  - rotate\_y\_wave, 142
  - rotate\_z\_wave, 142
  - set\_rotation\_x, 143
  - set\_rotation\_y, 143
  - set\_rotation\_z, 143
- pearl\_data\_explorer
  - pearl-data-explorer.ipf, 87
- pearl\_eolog
  - pearl-eolog.ipf, 102
- pearl\_file\_type
  - pearl-data-explorer.ipf, 87
- PearlAnglescanProcess, 17

- PearlAnglescanTracker
  - pearl-menu.ipf, [116](#)
- PearlAreaDisplay, [17](#)
- PearlAreaImport, [18](#)
- PearlAreaProfiles, [18](#)
- PearlArpes, [18](#)
- PearlDataExplorer, [18](#)
  - pearl-data-explorer.ipf, [87](#)
- PearlElog, [19](#)
- PearlElogPanel
  - pearl-elog.ipf, [102](#)
- PearlLiveDisplay
  - pearl-menu.ipf, [116](#)
- PearlMatrixImport, [19](#)
- PearlMenuEnableFunc
  - pearl-menu.ipf, [116](#)
- PearlPShellImport, [19](#)
- PearlScientaPreprocess, [19](#)
- pizza\_service
  - pearl-anglescan-process.ipf, [44](#)
- pizza\_service\_2
  - pearl-anglescan-process.ipf, [45](#)
- pmp\_data
  - pearl-anglescan-tracker.ipf, [56](#)
- pmp\_data\_mouseup
  - pearl-anglescan-tracker.ipf, [56](#)
- pmp\_export
  - pearl-area-display.ipf, [69](#)
- pmp\_parameters
  - pearl-anglescan-tracker.ipf, [56](#)
- pmp\_parameters\_mouseup
  - pearl-anglescan-tracker.ipf, [56](#)
- polar2cart
  - pearl-polar-coordinates.ipf, [118](#)
- polar2cart\_wave
  - pearl-polar-coordinates.ipf, [118](#)
- polar\_distance
  - pearl-polar-coordinates.ipf, [118](#)
- polar\_graph\_hook
  - pearl-anglescan-process.ipf, [46](#)
- precision
  - DoniachSunjicStruct, [21](#)
- prefs\_objects
  - pearl-anglescan-tracker.ipf, [58](#)
- prepare\_command\_line
  - pearl-elog.ipf, [102](#)
- prepare\_graph\_attachments
  - pearl-elog.ipf, [103](#)
- preview\_attributes
  - pearl-data-explorer.ipf, [88](#)
- preview\_datafolder
  - pearl-data-explorer.ipf, [88](#)
- preview\_dataset
  - pearl-data-explorer.ipf, [88](#)
- preview\_file
  - pearl-data-explorer.ipf, [88](#)
- preview\_hdf\_file
  - pearl-data-explorer.ipf, [88](#)
- preview\_itx\_file
  - pearl-data-explorer.ipf, [88](#)
- preview\_matrix\_file
  - pearl-matrix-import.ipf, [114](#)
- preview\_mtrx\_file
  - pearl-data-explorer.ipf, [88](#)
- preview\_pshell\_file
  - pearl-data-explorer.ipf, [89](#)
- preview\_setscale\_x
  - pearl-data-explorer.ipf, [89](#)
- process\_image\_data
  - pearl-anglescan-tracker.ipf, [56](#)
- prompt\_Shockley\_anglefit
  - pearl-scienta-preprocess.ipf, [139](#)
- prompt\_default\_process
  - pearl-data-explorer.ipf, [89](#)
- prompt\_func\_params
  - pearl-data-explorer.ipf, [89](#)
- prompt\_gauss2\_reduction
  - pearl-scienta-preprocess.ipf, [139](#)
- prompt\_hdf\_options
  - pearl-data-explorer.ipf, [89](#)
- prompt\_int\_linbg\_reduction
  - pearl-scienta-preprocess.ipf, [139](#)
- prompt\_int\_quadbg\_reduction
  - pearl-scienta-preprocess.ipf, [139](#)
- prompt\_redim\_linbg\_reduction
  - pearl-scienta-preprocess.ipf, [139](#)
- ps\_detect\_scale
  - pearl-pshell-import.ipf, [121](#)
- ps\_scale\_dataset
  - pearl-pshell-import.ipf, [122](#)
- ps\_scale\_dataset\_2
  - pearl-pshell-import.ipf, [122](#)
- ps\_scale\_datasets
  - pearl-pshell-import.ipf, [123](#)
- ps\_set\_dimlabels
  - pearl-pshell-import.ipf, [123](#)
- ps\_set\_dimlabels2
  - pearl-pshell-import.ipf, [124](#)
- psh5\_close\_file
  - pearl-pshell-import.ipf, [124](#)
- psh5\_list\_scan\_datasets
  - pearl-pshell-import.ipf, [124](#)
- psh5\_list\_scan\_regions
  - pearl-pshell-import.ipf, [125](#)
- psh5\_list\_scans
  - pearl-pshell-import.ipf, [125](#)
- psh5\_load\_complete
  - pearl-pshell-import.ipf, [125](#)
- psh5\_load\_dataset
  - pearl-pshell-import.ipf, [126](#)
- psh5\_load\_dataset\_meta
  - pearl-pshell-import.ipf, [127](#)
- psh5\_load\_dataset\_reduced
  - pearl-pshell-import.ipf, [127](#)
- psh5\_load\_dataset\_slab
  - pearl-pshell-import.ipf, [128](#)

- psh5\_load\_dataset\_slabs
  - pearl-pshell-import.ipf, [128](#)
- psh5\_load\_info
  - pearl-pshell-import.ipf, [129](#)
- psh5\_load\_preview
  - pearl-pshell-import.ipf, [129](#)
- psh5\_load\_reduced
  - pearl-pshell-import.ipf, [130](#)
- psh5\_load\_scan\_attrs
  - pearl-pshell-import.ipf, [131](#)
- psh5\_load\_scan\_complete
  - pearl-pshell-import.ipf, [131](#)
- psh5\_load\_scan\_data
  - pearl-pshell-import.ipf, [132](#)
- psh5\_load\_scan\_info
  - pearl-pshell-import.ipf, [132](#)
- psh5\_load\_scan\_meta
  - pearl-pshell-import.ipf, [133](#)
- psh5\_load\_scan\_preview
  - pearl-pshell-import.ipf, [133](#)
- psh5\_load\_scan\_section
  - pearl-pshell-import.ipf, [134](#)
- psh5\_open\_file
  - pearl-pshell-import.ipf, [134](#)
- pw
  - DoniachSunjicStruct, [22](#)
- quick\_pizza\_image
  - pearl-anglescan-process.ipf, [46](#)
- read\_attribute\_info
  - pearl-area-import.ipf, [77](#)
- record\_results
  - fermi-edge-analysis.ipf, [26](#)
- redim\_linbg\_reduction
  - pearl-scienta-preprocess.ipf, [139](#)
- reduce\_slab\_image
  - pearl-area-import.ipf, [77](#)
  - pearl-pshell-import.ipf, [135](#)
- reduce\_slab\_worker
  - pearl-area-import.ipf, [78](#)
  - pearl-pshell-import.ipf, [135](#)
- rotate2d\_x
  - pearl-vector-operations.ipf, [142](#)
- rotate2d\_y
  - pearl-vector-operations.ipf, [142](#)
- rotate\_hemi\_scan
  - pearl-anglescan-process.ipf, [46](#)
- rotate\_x\_wave
  - pearl-vector-operations.ipf, [142](#)
- rotate\_y\_wave
  - pearl-vector-operations.ipf, [142](#)
- rotate\_z\_wave
  - pearl-vector-operations.ipf, [142](#)
- SUCCESS
  - errorCode, [23](#)
- save\_hemi\_scan
  - pearl-anglescan-process.ipf, [47](#)
- save\_prefs
  - pearl-anglescan-tracker.ipf, [56](#)
  - pearl-data-explorer.ipf, [89](#)
  - pearl-elog.ipf, [103](#)
- save\_tracker\_data
  - pearl-anglescan-tracker.ipf, [57](#)
- scienta\_ang\_transm
  - pearl-scienta-preprocess.ipf, [140](#)
- scienta\_norm
  - pearl-scienta-preprocess.ipf, [140](#)
- scienta\_poly\_bg
  - pearl-scienta-preprocess.ipf, [140](#)
- select\_dataset
  - pearl-pshell-import.ipf, [135](#)
- set\_contrast
  - pearl-anglescan-process.ipf, [47](#)
- set\_panel\_attributes
  - pearl-elog.ipf, [103](#)
- set\_panel\_graphs
  - pearl-elog.ipf, [103](#)
- set\_panel\_message
  - pearl-elog.ipf, [104](#)
- set\_polar\_graph\_cursor
  - pearl-anglescan-process.ipf, [47](#)
- set\_rotation\_x
  - pearl-vector-operations.ipf, [143](#)
- set\_rotation\_y
  - pearl-vector-operations.ipf, [143](#)
- set\_rotation\_z
  - pearl-vector-operations.ipf, [143](#)
- set\_trace\_colors
  - pearl-area-display.ipf, [69](#)
- setup\_data
  - pearl-anglescan-tracker.ipf, [57](#)
- setup\_detector
  - pearl-anglescan-tracker.ipf, [57](#)
- setup\_graph
  - pearl-anglescan-tracker.ipf, [57](#)
- Shockley\_anglefit
  - pearl-scienta-preprocess.ipf, [140](#)
- show\_analyser\_line
  - pearl-anglescan-process.ipf, [47](#)
- show\_preview\_graph
  - pearl-data-explorer.ipf, [90](#)
- show\_shift
  - fermi-edge-analysis.ipf, [26](#)
- ShowComponents\_Au4f\_2p2
  - pearl-fitfuncs.ipf, [107](#)
- ShowComponents\_Au4f\_2p3
  - pearl-fitfuncs.ipf, [107](#)
- slit\_correction
  - fermi-edge-analysis.ipf, [27](#)
- slit\_shift
  - fermi-edge-analysis.ipf, [27](#)
- slp\_slice\_position
  - pearl-area-display.ipf, [69](#)
- strip\_delete\_frames
  - pearl-anglescan-process.ipf, [48](#)

- subtract\_line\_bg
  - pearl-matrix-import.ipf, [115](#)
- SumWavesIterator
  - pearl-tools.ipf, [142](#)
- svp\_slice\_position
  - pearl-area-display.ipf, [69](#)
- svp\_smoothing
  - pearl-area-display.ipf, [69](#)
- test\_attributes\_notebook
  - pearl-data-explorer.ipf, [90](#)
- test\_gauss2\_reduction
  - pearl-scienta-preprocess.ipf, [140](#)
- test\_int\_linbg
  - pearl-scienta-preprocess.ipf, [141](#)
- test\_int\_quadbg
  - pearl-scienta-preprocess.ipf, [141](#)
- test\_shockley\_anglefit
  - pearl-scienta-preprocess.ipf, [141](#)
- toggle\_capture
  - pearl-anglescan-tracker.ipf, [57](#)
- trim\_hemi\_scan
  - pearl-anglescan-process.ipf, [48](#)
- twave2list
  - pearl-pshell-import.ipf, [135](#)
- UNKNOWN\_ERROR
  - errorCode, [23](#)
- UnloadPearlArpesPackage
  - pearl-arpes.ipf, [82](#)
- update\_attach\_items
  - pearl-elog.ipf, [104](#)
- update\_capture
  - pearl-anglescan-tracker.ipf, [57](#)
- update\_data\_graph
  - pearl-anglescan-tracker.ipf, [57](#)
- update\_datasets
  - pearl-data-explorer.ipf, [90](#)
- update\_detector
  - pearl-anglescan-tracker.ipf, [57](#)
- update\_detector\_graph
  - pearl-anglescan-tracker.ipf, [58](#)
- update\_filelist
  - pearl-data-explorer.ipf, [90](#)
- update\_polar\_info
  - pearl-anglescan-process.ipf, [49](#)
- update\_progress\_panel
  - pearl-gui-tools.ipf, [107](#)
- update\_slice\_info
  - pearl-area-display.ipf, [69](#)
- version
  - pearl-anglescan-tracker.ipf, [58](#)
  - pearl-fitfuncs.ipf, [107](#)
- WAVE\_EXIST
  - errorCode, [23](#)
- WRONG\_PARAMETER
  - errorCode, [23](#)
- wave2list
  - pearl-pshell-import.ipf, [135](#)
- xdw
  - DoniachSunjicStruct, [22](#)
- xw
  - DoniachSunjicStruct, [22](#)
- yw
  - DoniachSunjicStruct, [22](#)